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RepeatOne[®]

User Manual

1. General	4
1.1. Design Principle of RepeatOne [®]	4
2. Volume Selection and Choice of TipOne [®] Repeat	5
2.1. TipOne [®] Repeat Dispenser Tips.	6
2.2. 25 ml and 50 ml TipOne [®] Repeat Dispenser Tip	6
3. Operation	7
3.1. Inserting a TipOne [®] Repeat Dispenser Tip	7
3.2. Filling the TipOne [®] Repeat Dispenser Tip	7
3.3. Dispensing.	8
3.4. Removing the TipOne [®] Repeat Dispenser Tip	9
4. Safety Precautions and Operational Limitations	9
5. Troubleshooting	10
6. Maintenance, Service and Warranty.	11
7. Ordering Information.	12
8. Technical Data.	13
9. Quick Reference Guide	14
10. Declaration of Conformity	63

1. General

RepeatOne® is a hand-held device for repetitive dispensing. Up to 48 dispenses of the same sample can be made from a single filling, resulting in a considerable reduction in time and expense.

RepeatOne® allows the selection of five separate volumes from each TipOne® Repeat Dispenser Tip. Seven different TipOne® Repeat Dispenser Tip sizes provide the most important volumes from 10 to 5000 µl.

Use only original STARLAB TipOne® Repeat Dispenser Tips to guarantee the pipette's performance.

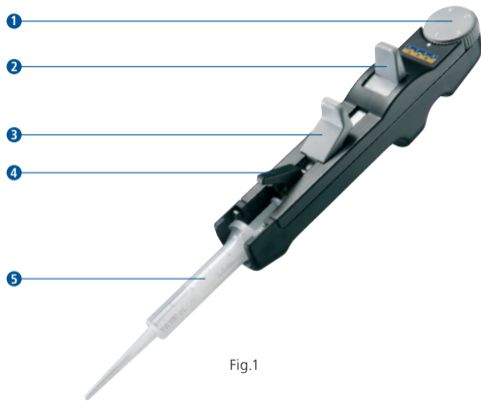
Fig.1

1.1. Design Principle of RepeatOne® (Fig.1)

- 1 Volume selection dial**
To determine the dispensing volume, the dial setting (1–5) is multiplied by the minimum pipetting volume of the TipOne® Repeat Dispenser Tip.
- 2 Dispensing lever**
The volume is dispensed by pressing the dispensing lever down until it stops.
- 3 Filling lever**
The TipOne® Repeat Dispenser Tip is filled by sliding this lever upward.
- 4 Locking clamp**
The locking clamp secures the TipOne® Repeat Dispenser Tip firmly into RepeatOne®.
- 5 TipOne® Repeat Dispenser Tip**

2. Volume Selection and Choice of TipOne® Repeat

For repetitive dispenses of a particular volume:

- Look for the desired volume in the upper horizontal line.
- In the column below, you will find the number of dispensing steps together with the appropriate setting of the volume selection dial (shaded).
- Select the corresponding TipOne® Repeat Dispenser Tip from the column on the right.

Pipetting Volume TipOne® Repeat																							TipOne® Repeat				
10 µl	20 µl	25 µl	30 µl	40 µl	50 µl	75 µl	100 µl	125 µl	150 µl	200 µl	250 µl	300 µl	400 µl	500 µl	750 µl	1 ml	1.25 ml	1.5 ml	2 ml	2.5 ml	3 ml	4 ml	5 ml				
48	23		15	11	8																				0.5 ml		
1	2		3	4	5																						
		48			23	15	11	8																		1.25 ml	
		1			2	3	4	5																			
					48		23		15	11	8															2.5 ml	
					1		2		3	4	5																
							48		23		15	11	8													5.0 ml	
							1		2		3	4	5														
									48		23	15	11	8												12.5 ml	
									1		2	3	4	5													
											48		23		15	11	8									25 ml*	
											1		2		3	4	5										
															48		23		15	11	8				50 ml*		
															1		2		3	4	5						

Number of dispensing steps

23
2

Volume selection dial

*Adapter necessary for TipOne® Repeat 25 ml and 50 ml.

2.1. TipOne® Repeat Dispenser Tips

The minimum dispensing volume ① and the maximum filling capacity ② is stated on each TipOne® Repeat Dispenser Tip (Fig. 2). The scale facilitates drawing off partial quantities and helps you to estimate the amount of liquid still available for operation.

The volume to be dispensed is obtained by multiplying the set number located on the volume selection dial by the minimum dispensing volume of the TipOne® Repeat Dispenser Tip.

Eg. 1.25 ml TipOne® Repeat, Minimum volume: 1=25 µl at setting 4: 4 x 25 µl = dispensed volume: 100 µl

Better precision is achieved by using a TipOne® Repeat Dispenser Tip of small capacity with a high dial setting.

Fig. 2

Number of Dispensing Steps	Position of Volume Selection Dial
48	1
23	2
15	3
11	4
8	5

The correlation between the position of the volume selection dial and the number of dispensing steps is as shown in the table above.

2.2. 25 ml and 50 ml TipOne® Repeat Dispenser Tips (Fig.3)

Note: An adapter, included with each RepeatOne®, is necessary for use with the 25 ml and 50 ml TipOne® Repeat Dispenser Tip.

Attach the adapter to the 25 ml or 50 ml TipOne® Repeat Dispenser Tip by turning it clockwise; remove by turning it counter clockwise.

Fig. 3

3. Operation

3.1. Inserting a TipOne® Repeat Dispenser Tip

- Slide the filling lever ① down until it stops.
- Raise the locking clamp ② upward.
- Insert the TipOne® Repeat Dispenser Tip until it clicks into position. Ensure the TipOne® Repeat's plunger is fully inserted into the barrel before attaching it to the RepeatOne® (Fig. 4).

Fig. 4

Fig. 5

- Ensure the filling lever is completely down, then lower the locking clamp to secure the TipOne® Repeat Dispenser Tip in place (Fig. 5).

3.2. Filling the TipOne® Repeat Dispenser Tip

The 0.5 ml, 1.25 ml and 2.5 ml TipOne® Repeat Dispenser Tip can be used with a standard 200 µl tip. Attaching a tip facilitates filling from narrow-necked vessels, piercing sealing foil

and makes dispensing of small volumes more precise.

- Immerse the tip into the liquid.
- Fill by slowly sliding the filling lever upward (Fig. 6).
- Discard the first dispensing step.

RepeatOne® is now ready for operation.

Fig. 6

Additional Information

Sliding the filling lever too quickly can cause excessive vacuum. This can cause tiny air bubbles to accumulate in the liquid which may lead to dispensing inaccuracies.

Important: The first dispensing step must be discarded in order to release any residual pressure from the dispensing system after filling and to prepare the system for precise dispensing. Small bubbles in the TipOne® Repeat Dispenser Tip beneath the piston will not affect dispensing accuracy since the residual stroke lock prevents the dispensing of any remaining sample after the last dispense.

It is not necessary to completely fill the TipOne® Repeat Dispenser Tip. Partially filling will not cause dispensing inaccuracies.

3.3. Dispensing (Fig. 7)

- Check the volume selection dial.
- Wipe the tip of the TipOne® Repeat Dispenser Tip with lint-free tissue.
- Hold the dispensing tip so it touches the inner wall of the vessel.
- Dispense the liquid by completely depressing the dispensing lever.

Fig. 7

Additional Information

Dispensing with a 5.0 ml or 12.5 ml TipOne® Repeat Dispenser Tip is normally carried out without attaching a pipette tip.

Never attach a pipette tip when using the 25 ml or 50 ml TipOne® Repeat Dispenser Tips.

A new, TipOne® Repeat Dispenser Tip must be used each time to avoid carry-over and contamination.

3.4. Removing the TipOne® Repeat Dispenser Tip

(Fig. 8)

- Hold the RepeatOne® over a beaker.
- Slowly depress the filling lever down to the stop. This will completely empty the TipOne® Repeat Dispenser Tip.
- Raise the locking clamp upward.
- Remove the TipOne® Repeat Dispenser Tip.

Note: When using the 25 ml or 50 ml TipOne® Repeat Dispenser Tips, the reusable adapter is removed by turning it counter clockwise.

Fig. 8

4. Safety Precautions and Operational Limitations

When removing TipOne® Repeat Dispenser Tips, please note the following:

- Do not touch the tip end, especially when using hazardous reagents!
- Empty the dispenser tips into a proper discard receptacle to avoid splashing and aerosols.
- Treat any retained fluid as a bio-hazard. Do not allow the fluid to come into contact with laboratory personnel.
- If dispensing a liquid with a surface tension significantly different from water, inaccuracies or leaking can occur.
- The RepeatOne® with an attached, partially filled TipOne® Repeat Dispenser Tip should be stored in a stand to prevent air entering the dispenser tip.

5. Troubleshooting

Error	Cause	Solution
TipOne® Repeat cannot be inserted.	Filling lever not fully in the down position; therefore, locking clamp cannot be completely raised.	Depress the filling lever downward and raise the locking clamp as far as it will go.
	The dispenser tip piston is not completely inserted in the barrel.	Slide the piston into the proper position.
Unable to fully lift the filling lever.	Locking clamp incompletely depressed.	Completely depress the locking clamp.
Excessive air bubbles when filling the TipOne® Repeat Dispenser Tip.	Filling lever lifted slightly before immersing into the liquid.	Draw up some solution and then discard it (using filling lever). Then refill the TipOne® Repeat Dispenser Tip.
	Leaking TipOne® Repeat Dispenser Tip.	Replace the dispenser tip with a new one.
Inconsistent dispensing.	Volume selection dial has not clicked into position.	Click volume selection dial into position.
Dripping TipOne® Repeat Dispenser Tip.	Leaking TipOne® Repeat Dispenser Tip.	Replace with a new dispenser tip.
	Attached TipOne® Repeat Dispenser Tip is loose.	Use more force to attach the TipOne® Repeat Dispenser Tip.
RepeatOne® jams or does not function.		Return the RepeatOne® for repair or replacement.

6. Maintenance, Service and Warranty

Maintenance

External surfaces of RepeatOne® can be cleaned with warm water or isopropyl alcohol.

Apart from the need to clean externally, the RepeatOne® is completely maintenance free.

Important: under no circumstances should the RepeatOne® be dismantled!

RepeatOne® and TipOne® Repeat are not autoclavable!

Service

If a problem cannot be solved with the aid of the Troubleshooting Guide, return RepeatOne® to your regional STARLAB office or distributor for repair or replacement.

Important: Please decontaminate the dispenser before return. Proper decontamination has to be confirmed in a separate “declaration of decontamination”.

Please contact STARLAB for more detailed information.

Warranty

Please check that the unit has been received complete and undamaged. Save all packaging and documents until you have thoroughly inspected your shipment; if you find that your order is incorrect or damaged please call STARLAB for instructions for its return.

This warranty is valid for 24 months, but only if the product has been used and cared for according to this User Manual. No liability is accepted for loss or damage arising from incorrect use.

The liability is limited to the repair or replacement of the unit, or a refund of the purchase price, at STARLAB's option. STARLAB is not liable for any consequential damages. We refer to our general terms of sale. STARLAB reserves the right to alter the specifications without prior notice.

7. Ordering Information

RepeatOne® Dispenser and Adapter		
Cat-No.	Description	Pack Size
S4026-1500	Dispenser RepeatOne®, incl. Adapter for 25 ml and 50 ml Dispenser Tips	1
S4026-1525	Adapter for 25 and 50 ml TipOne® Repeat Dispenser Tips*	1

* Sterile adapters are available on special request

Important Note:

Please only use the accessories recommended by STARLAB. Using non-recommended disposables can reduce the precision and accuracy, and life, of RepeatOne®. We do not honour any warranty or accept any responsibility for damage resulting from such action.

TipOne® Repeat Dispenser Tips		
Cat-No.	Description	Pack Size
S4770-0050	TipOne® Repeat, 0.5 ml Dispenser Tips	100
S4772-0125	TipOne® Repeat, 1.25 ml Dispenser Tips	100
S4774-0250	TipOne® Repeat, 2.5 ml Dispenser Tips	100
S4776-0500	TipOne® Repeat, 5.0 ml Dispenser Tips	100
S4778-1250	TipOne® Repeat, 12.5 ml Dispenser Tips	100
S4780-2500	TipOne® Repeat, 25 ml Dispenser Tips	50
S4782-5000	TipOne® Repeat, 50 ml Dispenser Tips	25

TipOne® Repeat Dispenser Tips (Sterile)		
Cat-No.	Description	Pack Size
S4784-0050	TipOne® Repeat, 0.5 ml Dispenser Tips (Sterile)	100
S4786-0125	TipOne® Repeat, 1.25 ml Dispenser Tips (Sterile)	100
S4788-0250	TipOne® Repeat, 2.5 ml Dispenser Tips (Sterile)	100
S4790-0500	TipOne® Repeat, 5.0 ml Dispenser Tips (Sterile)	100
S4792-1250	TipOne® Repeat, 12.5 ml Dispenser Tips (Sterile)	100
S4794-2500	TipOne® Repeat, 25 ml Dispenser Tips (Sterile)	50
S4796-5000	TipOne® Repeat, 50 ml Dispenser Tips (Sterile)	25

8. Technical Data

TipOne® Repeat Dispenser Tip Size	Dispensing Volume			Systematic Error (Inaccuracy)			Random Error (Imprecision; CV)		
	Min.		Max.	Min.		Max.	Min.		Max.
0.5 ml	10 µl	30 µl	50 µl	±2.0 %	±1.0 %	±1.0 %	±2.0 %	±1.0 %	±0.8 %
1.25 ml	25 µl	75 µl	125 µl	±0.9 %	±0.9 %	±0.9 %	±0.9 %	±0.7 %	±0.6 %
2.5 ml	50 µl	150 µl	250 µl	±0.7 %	±0.7 %	±0.7 %	±0.9 %	±0.7 %	±0.6 %
5 ml	100 µl	300 µl	500 µl	±0.6 %	±0.6 %	±0.6 %	±0.6 %	±0.5 %	±0.3 %
12.5 ml	250 µl	750 µl	1250 µl	±0.5 %	±0.5 %	±0.5 %	±0.6 %	±0.5 %	±0.3 %
25 ml	500 µl	1500 µl	2500 µl	±0.4 %	±0.4 %	±0.4 %	±0.6 %	±0.4 %	±0.25 %
50 ml	1000 µl	3000 µl	5000 µl	±0.3 %	±0.3 %	±0.3 %	±0.5 %	±0.4 %	±0.25 %

Test conditions in accordance with EN ISO 8655

Liquid: Water bi-distilled, degassed

Test Temperature: 20–25°C, ±0.5 % constant

(Data for TipOne® Repeat Dispenser Tips 0.5 ml, 1.25 ml and 2.5 ml are determined with attached 200 µl pipette tip)

Operating Temperature: +4 °C to +40 °C

Measurement: Performed with original TipOne® Repeat Dispenser Tips in accordance with EN ISO 8655

Technical specifications are subject to change without notice!

9. Quick Reference Guide

Important: The Quick Reference Guide should only be used if you are thoroughly acquainted with the detailed instructions, otherwise errors may occur.

- ✦ Look for desired volume in the volume selection chart (See 2.).
- ✦ Select appropriate TipOne® Repeat Dispenser Tip by number of dispensing steps.
- ✦ Set volume selection dial to corresponding position.
- ✦ Slide filling lever down.
- ✦ Raise locking clamp upward.
- ✦ Insert TipOne® Repeat Dispenser Tip into position.
- ✦ Lower locking clamp and press closed.
- ✦ Immerse the tip of the TipOne® Repeat Dispenser Tip into liquid.
- ✦ Slide filling lever upward.
- ✦ Discard first dispensing step.
- ✦ Dispense.
- ✦ Empty TipOne® Repeat Dispenser Tip completely.
- ✦ Raise locking clamp upward.
- ✦ Remove TipOne® Repeat Dispenser Tip.

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Declaration of Conformity

Product Description
Catalogue Number

Manual Dispenser RepeatOne®
S4026-1500

STARLAB herewith certifies that the product has been manufactured and inspected in accordance with the following International and European standards:

EN ISO 8655
ASTM E1154-89 (2003)
98/79/EEC

Piston-Operated Volumetric Apparatus
Standard Specification for Piston or Plunger Operated Volumetric Apparatus
In Vitro Diagnostic Medical Devices

*As an ISO 9001 and ISO 13485 certified company, our working procedures and processes are checked and audited regularly.
Additionally, the quality of our products is continuously monitored.*

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**STARLAB Sientific
India Private Limited**

First Floor, Old No. 322, New No. 24
Thambu Chetty Street
Chennai-600001
India
T: +91(0)44 42130702/03
F: +91(0)44 42130701
info@starlab.co.in
www.starlab.co.in

www.starlabgroup.com

Tolerance Check · Meßprotokoll · Protocole de mesure
Protocollo di miscura · Protocolo de medición

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STARLAB Srl.