Plate Centrifuge

Instruction Manual

N2631-0008

Plate Centrifuge

www.starlabgroup.com

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N2631-0008 Plate Centrifuge

Important Information

Latest and up-to-date version of the declaration of conformity is available at:

www.starlab.click/doc-platecentrifuge

Plate Centrifuge

Cat. No. N2631-0008

Safety Precautions

Please read the entire instruction manual before operating the Plate Centrifuge for the first time.

Please observe the following notes:

Never put hands in the rotor area unless the rotor is completely stationary.

Never move the centrifuge while the rotor is spinning.

Never use solvents or flammables near this or other electrical equipment.

Never centrifuge flammable, explosive or corrosive materials.

Never centrifuge hazardous materials outside of a hood or proper containment facility.

Never continue to operate a centrifuge that is vibrating excessively.

Never use the centrifuge in any manner not specified in these instructions.

Always load the rotor symmetrically. Each plate should be counter balanced by another plate of the same brand, type and volume.

Always use plates that have been sealed in some manner to avoid spillage.

Always load plates so the sealed top faces the centre of the rotor and the bottom of the plate faces out.

Always be sure the plates are pushed all the way to the bottom of the rotor slot before operating the centrifuge.

Always locate the centrifuge within easy access to an electrical outlet.

Product Description

The Plate Centrifuge is designed for the centrifugation of small samples in PCR plates with 96 or 384 Wells. All standard PCR plates are suitable: skirted, non-skirted or semi-skirted.

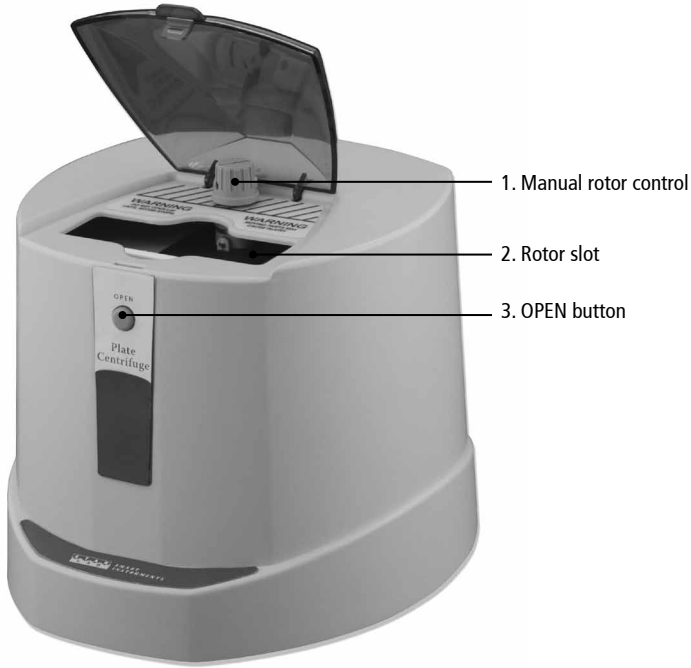
Figure 1: Control elements

First Steps

Packaging: Take the centrifuge from the package. Keep the packaging for shipping or storage. Check the contents.

Contents: Plate centrifuge x 1. Power cord x 1. Instruction manual x 1.

Setting Up the Plate Centrifuge

Choose a location with an easily accessible power outlet.

1. Choose a clean, clutter-free surface and make sure that the feet of the centrifuge are free from obstructions.
2. Place the centrifuge on a flat, stable bench.
3. Make sure that the ventilation of the centrifuge is not impeded and keep sufficient distance to other units.

Note: Both the connection for the power cord and the power switch can be found at the back of the centrifuge. An indication of the requirements for the power supply and frequency (which must be matched) can be found on the label on the underside of the unit.

4. Plug the power cord into the back of the unit.
5. Connect the power cord to the power socket.
6. Turn the power switch at the back of the unit to the 'ON' position.

Operation

Loading: Plates must to be sealed carefully with a heat-sealing or self-adhesive plate seal and loaded in the rotor in a vertical position. See Fig. 2.

Important: If you have only one PCR plate to centrifuge you **MUST** load a 'dummy' plate of the same brand, style and sample volume (water can be used) in the other rotor to balance the centrifuge.

Essential! The rotor must be stationary before loading a plate.

1. Press the OPEN button (3) to release the lid of the centrifuge.
2. Open the lid completely
3. Position the rotor by spinning the manual rotor control (1) until a complete rotor slot is visible.

Note: The manual rotor control (1) should be easy to rotate. If you find it difficult to turn, close the lid and run the unit for a few seconds to resolve the friction brake. Stop the unit by pressing the OPEN button (3) and wait until the rotor has come to a complete standstill before repeating steps 2 to 3.

During insertion of the plates, make sure that the sealed top of the plate faces the centre of the rotor and the bottom of the plate faces out. Be sure that the plate is pushed all the way down into the slot.

4. Place the first plate into the rotor slot (2).
5. Reposition the rotor by spinning the manual rotor control (1), until the second rotor slot is visible.
6. Insert second plate.

Figure 2: Correct loading of a PCR plate

To Start Centrifuging

1. Close the lid of the centrifuge. Centrifugation starts automatically.

A run of 15 seconds is sufficient to spin samples to the bottom of the wells.

Note: A slight vibration may be felt at the centrifuge's full speed of 2500 rpm. Excessive vibration indicates that the rotor is not balanced. In this case stop the run immediately.

DO NOT run the centrifuge unbalanced. Always use a balance plate of the same brand, type and sample volumes.

DO NOT centrifuge with a damaged rotor.

To Stop Centrifuging

1. Press the OPEN button. The lid is released, the friction brake is activated and the rotor slows to a stop.
2. DO NOT open the lid until the rotor is completely still.
3. Reposition the rotor by spinning the manual rotor control (1) until a rotor slot is accessible.

Note: The manual rotor control (1) should be easy to rotate. If you find it difficult to turn, close the lid and run the unit for a few seconds to resolve the friction brake. Stop the unit by pressing the OPEN button and wait until the rotor has come to a complete standstill before repeating steps 2 and 3.

4. Remove the first plate from the rotor slot.
5. Reposition the rotor by spinning the manual rotor control (1) until the second rotor slot is visible.
6. Remove the second plate from the rotor slot.

Cleaning

Note: Use only pH-neutral detergents (pH-value 6–8) for cleaning.

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DO NOT immerse the centrifuge in liquid.

Always keep the rotor and rotor slots clean.

1. Wipe the outside of the centrifuge with a clean, damp cloth.
2. The rotor slots may be cleaned with a sponge.

The rotor can be dismantled from the centrifuge when a more thorough cleaning is necessary, see opposite for instructions.

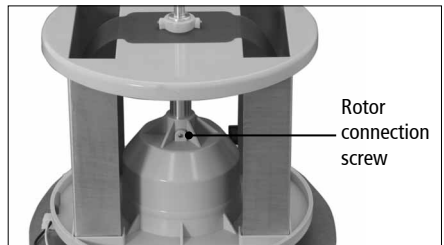
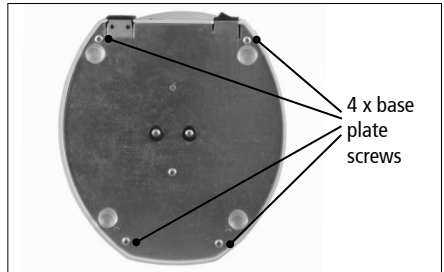
Rotor Disinfection

Note: Remove the rotor from the centrifuge before you start disinfecting. Instructions on how to remove the rotor is described below.

Decontamination of the centrifuge is required if infectious material is passing into the interior of the rotor. Use neutral disinfectants only. A disinfectant spray is ideal to reach inaccessible places.


Dismounting the Rotor for Cleaning

1. Disconnect the unit from the power supply.
2. Turn the unit upside down and lay on a soft cloth to avoid scratches.
3. Remove all four outer screws on the base plate.
4. Turn the centrifuge right way up
5. Loosen the rotor knob set screw on the manual rotor control (1) with a 1/16" Allen key.
6. Pull off the manual rotor control (1).
7. Lift off housing.
8. Loosen the rotor connection screw with a .050" Allen key.
9. Remove the rotor by lifting upward.

Reinstalling the Rotor

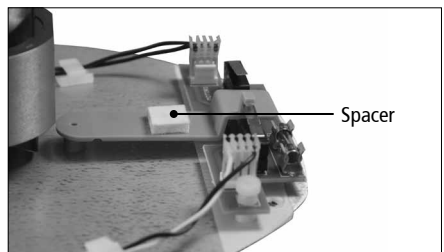Note: Hold 0.5" distance from the bottom plate of the centrifuge. Use a suitable spacer.

Maintenance and Service

Disconnect the centrifuge from the mains before you carry out any maintenance or service work. The Plate Centrifuge with its brushless motor does not require regular maintenance or service work. Repairs should be performed only by qualified service personnel or the STARLAB service. The warranty is void if repairs were previously performed by untrained personnel.

Warranty

The Plate Centrifuge and rotor are covered by a two-year warranty which covers the centrifuge and rotor. This warranty is void in the event of incorrect operation, use of non-standard spare parts or accessories, and unauthorised modification of the rotor or centrifuge. The rotor warranty does not cover normal wear and tear, nor does it cover damage and/or weakening caused by radioactive contamination, cleaning agents, or over tightening of the Allen screw. Fuses are not covered by the warranty.

Should you not be satisfied with your product, or defects/damage to the unit are found, please contact your local STARLAB office or distributor (see www.starlabgroup.com).

Technical Specifications

Model	Plate Centrifuge
Maximum speed	2,500 rpm
Maximum RCF	500 x g
Maximum rotor capacity	2 PCR plates
Power supply	230 V / 50~60 Hz; 0.4 A (Fuse 0.5 AT)
Dimensions (WxDxH)	190 x 210 x 183 mm
Weight	1.5 kg

Features and specifications subject to change without notice

Ordering Information

N2631-0008	Plate Centrifuge
N2631-0018*	Adapter for Plate Centrifuge

* Adapter plate available for use with 0.2ml PCR tubes and PCR strips and 48-well PCR Plates.

Please see our website www.starlabgroup.com for information on our wide range of PCR Consumables, comprising plates, tubes, strips and plate seals.

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STARLAB GmbH
STARLAB (UK), Ltd
STARLAB S.r.l.