This is your **Vortex** user manual.

USER MANUAL · BEDIENUNGSANLEITUNG
MODE D'EMPLOI ISTRUZIONI PER L'USO · INSTRUCCIONES DE MANEJO

Passionate for science.

EC Declaration of Conformity

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Wir / We / Nous:

Starlab International GmbH

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EN 61010-1 **Safety Requirements for Electrical Equipment for Measurement
Control, and Laboratory Use – Part 1: General Requirements**

EN 61326-1 **Electrical equipment for measurement, control and laboratory
use - EMC requirements - Part 1: General requirements**

EN 61000-3-2 **Electromagnetic compatibility (EMC) - Part 3-2: Limits - Limits for
harmonic current emissions**

EN 61000-3-3 **Electromagnetic compatibility (EMC) - Part 3-3: Limits - Limita-
tion of voltage changes, voltage fluctuations and flicker**

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Authorised Managing Director
Klaus Ambos

Registry Court
Amtsgericht [first instance court]
Hamburg
Register number: HRB 122020
VAT ID no. pursuant to
§ 27 a German
VAT Act DE 214 891 717

ARTICLE NO. N2400-6110

1. Safety instructions

Before operation

- Read the Instruction Manual carefully before use.
- Ensure that only trained staff use the instrument. The instrument may only be opened by a qualified and trained technician.
- Make sure the socket is earthed before use.
- Set the instrument on a stable, clean, dry, non-slip and fireproof surface.
- Do not operate in hazardous areas.
- Safe operation is only guaranteed with the original accessories from STARLAB. You will find the complete list in chapter 8.
- Accessories must be securely attached.
- Check instrument and accessories for damage before each use. Do not use damaged components.
- The operation voltage mentioned on the label must match.
- Always disconnect the plug before fitting accessories.
- When working, wear personal protection to avoid the risk from splashing of liquids and possible glass breakage which can result from mechanical vibrations.

During operation

- Hold the tube as vertically as possible.
- Place a tube at the centre of the mixing head. A sharp-edged tube can wear the mixing head.
- Set the “speed-control knob“ to the lowest position then gradually increase the speed.
- Reduce the speed if the medium splashes out of the vessel or the instrument is unstable on the bench.

2. Product description

The instrument is designed for mixing liquids.

Normal use:

The instrument is designed for use in professional laboratories. Do not use it in residential areas.

3. Inspection

Unpack the equipment carefully and check for any damages. If the instrument is faulty you should contact Starlab or your Starlab distributor promptly.

The packing includes:

- 1 Vortex incl. universal attachment N2400-6011
- 1 EU-power supply for N2400-6010 or 1 UK-power supply for N2400-6110
- 1 Instruction manual

Please check all items are present.

4. Ambient conditions

- Power supply: 100 – 240 V 50/60Hz
- Ambient temperature: 4 – 40 °C
- Relative humidity: < 80 %

Place the instrument in safe and stable surface.

5. Operation

5.1 Control Elements

5.2 Operating

Permanent operation

1. Place the Vortex on the table, turn the speed control knob on the lowest position and plug-in the power supply.
2. Push the two-direction switch leftward to ON. The power LED lights and the instrument is in the “standby” mode.
3. Turn on the Vortex by pushing the speed control knob clockwise. Set a rated speed. The mixing head begins to vibrate and you can press your test tube vertically on the mixing head.
4. If you wish to finish, push the two-direction-switch rightwards to OFF, LED light goes off and the instrument is turned off.

Touch operation

1. Place the Vortex on the table and plug-in the power supply
2. Push the two-direction switch rightward to the touch mode. The power LED lights and the instrument is in the “standby” mode.
3. Turn on the Vortex by turning the speed control knob clockwise. Set a rated speed. Press the test tube vertically on to the mixing head, and the instrument begins to work.
4. If you wish to finish, push the two-direction-switch leftwards to OFF. The power LED is now off and instrument turns off.

6. Maintenance and cleaning

- Proper maintenance and cleaning will keep the instrument in good working order and will lengthen its life time.
- Be careful not spray any cleaning fluid into the instrument. Unplug the power cord when cleaning.
- Please use only Isopropyl alcohol and water containing tenside. Before using other methods or cleaning agents, the user must ascertain with the manufacturer that this will not destroy the instrument.
- Wear protective gloves during cleaning of the instrument.
- When not in use for long periods, switch off the instrument and store it in a dry, clean and stable place at room temperature.

Warranty

This instrument is under warranty to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty claim expires if the instrument is opened by untrained or unqualified staff. Our “General terms and conditions” apply.

Service

If help is needed, please contact Starlab or your Starlab distributor. See www.starlabgroup.com for your local contact.

Please provide the customer care representative with the following information:

- Serial Number (on the rear panel)
- Description of problem (i.e., hardware or software)
- Methods and procedures adopted to resolve the problems
- Your contact information

7. Technical specifications

Parameter	Data
Voltage	220 – 230/110 – 120 V
Frequency	50/60 Hz
Motor power	60 W
Motor rating input	58 W
Motor rating output	10 W
Shaking movement	Orbital
Orbital diameter	4 mm
Motor type	Shaded pole motor
Speed range	0 – 2,500 rpm
Dimension	127 x 130 x 160 mm
Net weight	3.5 kg
Permitted ambient temperature	4 – 40 °C
Permitted relative humidity	< 80 %
IP-protection	IP21

8. ordering information

Article No.	Description
N2400-6010	Vortex with EU-plug
N2400-6110	Vortex with UK-plug

8.1 Accessories

Article No.	Description
N2400-6011	Replacement universal cup Universal top plate for vessels up to 30 mm Ø
N2400-6012*	Tube adapter for 15 x 1.5/2.0 ml microcentrifuge tubes (up to 10 mm Ø). Use in combination with N2400-6015
N2400-6013*	Tube adapter for 8 x 15 ml centrifuge tubes (up to 16 mm Ø). Use in combination with N2400-6015
N2400-6014	Adapter for 96-well microtitre plates
N2400-6015	Vortex Adapter Plate for use with N2400-6012 or N2400-6013

* N2400-6012 and N2400-6013 can also be used as an attachment on Starlab's Microplate Mixer in combination with its adapter N2400-8043. See website for ordering details

Please visit our website www.starlabgroup.com, for details on our comprehensive range of plastic consumables for mixing.

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Starlab (UK), Ltd
Starlab S.r.l.