This is your user manual for Thermoblocks.

USER MANUAL · BEDIENUNGSANLEITUNG
MODE D'EMPLOI ISTRUZIONI PER L'USO · INSTRUCCIONES DE MANEJO

Passionate for science.

User Manual	3
Bedienungsanleitung	9
Manuel d'utilisation	15
Istruzioni per l'uso	21
Ordering information	30-31

1 USER INSTRUCTIONS

1.1 Using this manual

These instructions for the use of exchangeable thermoblocks are a supplement to the operating manual for the Mixer HC. They are not intended as a replacement for that operating manual.

Therefore, please also read the operating manual of the Mixer HC before using the exchangeable thermoblock for the first time.

2 PRODUCT DESCRIPTION

2.1 Features

The exchangeable thermoblocks for the Mixer HC are used for convenient temperature control and mixing of all standard lab tubes, Multiwell Plates and Deepwell Plates. The following exchangeable thermoblocks can be set on the Mixer HC:

Description	Speed Range in rpm
Thermoblock for 24 x 0.5 ml microcentrifuge tubes	300 to 1500
Thermoblock for 24 x 1.5 ml microcentrifuge tubes	300 to 1400
Thermoblock for 24 x 2.0 ml microcentrifuge tubes	300 to 1400
Thermoblock for 24 x microcentrifuge tubes, 12 mm	300 to 1400
Thermoblock for 8 x 5.0 ml preparation tubes	300 to 750
Thermoblock for 8 x 15 ml conical centrifuge tubes	300 to 750
Thermoblock for 4 x 50 ml conical centrifuge tubes	300 to 750
Thermoblock for 24 x 1.5/2.0 ml cryo tubes	300 to 1400
Thermoblock for Multiwell Plates / Deepwell Plates with lid	300 to 1400
Adapter plate 96 x 0.2 ml for PCR plates and tubes for MTP / DWP thermoblock	300 to 800 when loading >200 g

3 INTENDED USE

The exchangeable thermoblocks are intended exclusively for indoor use to hold tubes and plates during sample preparation.

The exchangeable thermoblocks must only be operated by appropriately qualified and instructed personnel. Carefully read the operating manual before working with the exchangeable thermoblocks and familiarize yourself with their use.

4 INSTALLATION AND OPERATION

4.1 Mounting exchangeable thermoblocks

Warning! Risk of burns from hot surfaces.

The exchangeable thermoblock and the thermal base can be very hot after heating and cause burns.

- › Allow heated exchangeable thermoblocks and the thermal base to cool down completely before removing the exchangeable thermoblock.
-

Warning! Injury or property damage from improperly attached exchangeable thermoblocks.

If the exchangeable thermoblocks are not attached properly, they can become detached from the device or break.

- › Tighten the screws of the exchangeable thermoblocks so they are hand-tight.
-

Warning! Contamination due to opening tube lids.


In the following cases, the lids of tubes can spring open and the sample material can escape:

- high vapor pressure of the content,
 - improperly sealed lid,
 - damaged sealing lip.
- › Always check that all tubes have been sealed tightly before use.
-

To insert the exchangeable thermoblocks, use the accessories supplied (eg. screw driver, screws, flat washers and lock washers).

4.1.1 Exchangeable thermoblocks for micro test tubes (microcentrifuge / cryogenic / microtubes)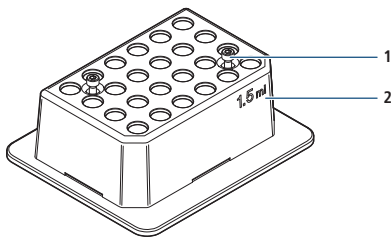

1. Place the exchangeable thermoblock (2) on the thermosocket with both flat edges and the writing facing the front.
2. Tighten the captive screws (1).


4.1.2 Exchangeable thermoblock for Multiwell Plates / Deepwell Plates

Caution! Microplates melt at extremely high temperatures.

- Only heat microplates made of polystyrene to a maximum 70 °C.

-
1. Insert the short screws in the MTP exchangeable thermoblock with the flat washers.
 2. Secure the screws from below using the lock washers.
The lock washers prevent the screws from falling out of the bores if the exchangeable thermoblock is not fixed on the unit.
 3. Place the MTP thermoblock on the device in such a way that the spring plate is to the front left.
 4. Tighten the screws.

4.1.3 Adapter plate for 0.2 ml tubes or 96-well PCR plates

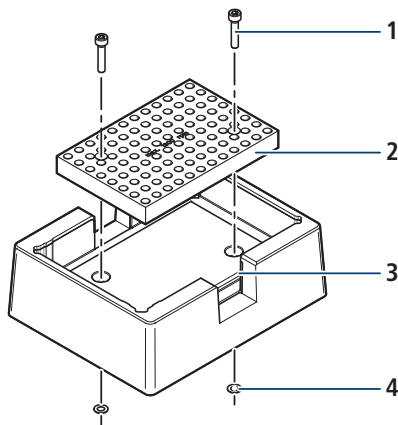1. Remove the lock washers (4) from the short screws in the exchangeable thermoblock.
2. Remove the short screws.
3. Connect the adapter plate (2) to the MTP thermoblock (3) using the long screws (1).
4. Secure the screws from below using the lock washers.
The lock washers prevent the screws from falling out of the bores if the exchangeable thermoblock is not fixed on the unit.
5. Place the adapter plate on the device with the MTP thermoblock.
6. Tighten the screws.

4.2 Inserting tubes and plates

Caution! Microplates melt at extremely high temperatures.

- › Only heat microplates made of polystyrene to a maximum 70 °C.
-

1. Place the tubes or plate on the thermoblock by pressing gently down.
For optimum temperature control, transfer the micro test tubes to the exchangeable thermoblocks using Iso-Racks. The Iso-Rack reduces the heat exchange of the thermoblock with the environment. If necessary close the Iso-Rack lid.
3. If using the microplate thermoblock: put on the lid.

5 MAINTENANCE

5.1 Cleaning

Danger! Electric shock as a result of penetration of liquid.

- › Switch off the Mixer HC and disconnect it from the power supply before starting cleaning or disinfecting the thermoblocks.
 - › Do not allow any liquids to enter the inside of the housing.
 - › Do not use spray disinfection.
 - › Only reconnect the unit to the power supply once it is completely dry.
-

Clean the exchangeable thermoblocks immediately if sample fluid enters the bore holes or the surface.

1. Clean with a mild, soap-based solution, using a lint-free cloth.
2. Wipe off the soap solution with distilled water.
3. Dry all cleaned parts.

5.2 Disinfection / Decontamination

1. Disinfect the unit.
2. Select a disinfection method which complies with the legal requirements and regulations applicable to your range of application.
3. Wipe the surfaces with the lint-free cloth and disinfectant.

If you are shipping the unit to the authorised Technical Service for repairs or to your local distributor for disposal, please note the following:

Warning! Risk to health from contaminated unit

- Follow the instructions for the decontamination supplied by your local distributor.
 - Decontaminate all the parts you want to dispatch.
 - Enclose the fully-completed decontamination certificate for returned goods (incl. the serial number of the unit) with the dispatch.
-

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Starlab (UK), Ltd
Starlab S.r.l.