# This is your hand controller 8-channel adapter for pipette tips user manual.

USER MANUAL · BEDIENUNGSANLEITUNG

Passionate for science.

 starlab

<b>1</b>	<b>Safety information</b>	<b>4</b>
1.1	Important information	4
1.2	General information	5
1.3	Intended use	5
1.4	Setting up	5
1.5	Ambient conditions	6
1.6	Operating conditions	6
1.7	Safety during operation	7
1.8	Maintenance and repair	8
<b>2</b>	<b>Technical data</b>	<b>8</b>
2.1	Materials of the 8-channel adapter	9
2.2	System parts	9
<b>3</b>	<b>Use and operation</b>	<b>10</b>
3.1	Assembling the adapter	10
3.2	Assembling the bench stand	11
3.3	Prior to aspiration	12
3.4	Aspiration of liquids	12
3.5	Ejecting the tips	12
3.6	After aspiration	13
<b>4</b>	<b>Cleaning and decontamination</b>	<b>14</b>
<b>5</b>	<b>Accessories and spare parts</b>	<b>16</b>
<b>6</b>	<b>Troubleshooting</b>	<b>17</b>
<b>7</b>	<b>Repair - Maintenance - Return</b>	<b>18</b>
<b>8</b>	<b>Return Form</b>	<b>19</b>

# 1 Safety information

## 1.1 Important information

### WARNING

- > Keep this manual complete and accessible to personnel at all times!
- > Read this manual carefully before installing or operating the equipment. Observe the instructions contained in this manual.
- > Do not modify the equipment without authorization.

### NOTICE

**This manual is an integral part of the equipment described therein. It describes the safe and proper use of the equipment. Make operating personnel aware of dangers arising from the device and the used substances. Starlab cannot accept any liability for inappropriate use of the equipment or for damage from failure to follow instructions contained in this manual.**

This manual is only to be used and distributed in its complete and original form. It is strictly the users' responsibility to check carefully the applicability of the guidance in this manual with respect to his product.

01/02/2023

**The following signal word panels and safety symbols are used throughout this manual:**

This is the safety alert symbol. It is used to alert you to potential personal injury hazards. Obey all safety messages that follow this symbol to avoid possible injury and death.

### DANGER

- > DANGER indicates a hazardous situation which, if not avoided, will result in death or serious injury.

### WARNING

- > WARNING indicates a hazardous situation which, if not avoided, could result in death or serious injury.

### CAUTION

- > CAUTION indicates a hazardous situation which, if not avoided, could result in minor or moderate injury.

### NOTICE

NOTICE is used to address practices not related to personal injury.

## 1.2 General information

### WARNING

- › Read and comply with this manual before installing or operating the equipment.
- › Before operating the equipment read and comply with the section "Cleaning and decontamination"!

### NOTICE

Remove all packing material, remove the product from its packing-box and inspect the equipment.

If the equipment is damaged, notify the supplier within three days; state the item number of the product together with your order number and the supplier's invoice number. Retain all packing material for inspection.

**Do not use the equipment if it is damaged.**

## 1.3 Intended use

### WARNING

- › The device may only be used by well trained and skilled personnel.
- › The device may not be used for in vivo applications (e. g., not humans or animals).

### NOTICE

Ensure that the equipment and all components are suitable for the intended application. Use the system only for **aspiration of liquids**.

## 1.4 Setting up

### NOTICE

The 8-channel tip adapter with ejector is intended to be connected to a Starlab hand controller. The hand controller is intended to be connected to a suitable suction system, e. g., Starlab Aspirator systems.

Comply with all applicable and relevant safety requirements (regulations and guidelines), implement the required actions and adopt suitable safety measures.

## 1.5 Ambient conditions

- > The systems are designed for **ambient temperatures** during operation between +10 °C and +40 °C. Check the maximum temperatures.

## 1.6 Operating conditions

- > The devices are **not suitable** to aspirate:
  - flammable mixtures
  - unstable substances
  - substances which react explosively under impact (mechanical stress) and/or when being exposed to elevated temperatures without air,
  - self inflammable substances,
  - substances which are inflammable without air and
  - explosive substance
- > The devices have **no approval** for operation below ground.

- > The devices are not suitable for aspiration of dust.

- > **Take into consideration interactions and chemical reactions of the aspirated media.**
- > Ensure that the materials of the wetted parts are compatible with the aspirated substances, see section "Technical data".

### NOTICE

If aspirating **different substances**, it is recommended that the aspiration hose is purged prior to changing the media in order to rinse out residues and to avoid reactions of the substances either with each other or with the materials.

## 1.7 Safety during operation

### DANGER

- › Ensure that while operating the device no persons can be endangered.
- › Wear personal protection equipment.
- › Comply with requirements when using infectious liquids, pathogenic germs, toxic, radioactive or aggressive substances.
- › Prior to operation check all relevant safety and health requirements and determine the applicability of regulatory limitations prior to use if necessary.
- › Comply with material safety data sheets and notes on safe use of from the manufacturer.
- › Comply with the biosafety level of the laboratory.
- › Use appropriate safety work materials and methods if necessary, e. g., steam sterilization, sterilization indicators or germicides. For notes concerning the sterilization of wetted parts, see the "Technical Data" section and "Cleaning and decontaminating". It is the user's responsibility to ensure effective sterilization.
- › Never operate a defective or damaged device.

### WARNING

- › Comply with applicable regulations when disposing of chemicals. Take into consideration that chemicals may be polluted.
- › Use only **OEM spare parts and accessories**; otherwise safety and performance of the equipment as well as the electromagnetic compatibility of the equipment might be reduced.
- › In case of leaks, aspirated substances might be released into the environment. Comply especially with notes on operation and use.

### NOTICE

Ensure that, in case of failure, the system will always return to a safe status. Provide appropriate protective measures (i.e., precautions which allow for the requirements of the respective application) even for the case of failure and **malfunction**.

## 1.8 Maintenance and repair

- > **Attention:** The device might be contaminated with noxious or otherwise dangerous substances that have been aspirated during operation. Ensure that the device is decontaminated before maintenance.

- > Take adequate precautions to protect people from the effects of dangerous substances if contamination has occurred. Wear appropriate safety clothing when you come into contact with contaminated components.

### NOTICE

Ensure that **maintenance** is carried out by suitably trained technicians. Ensure that the maintenance technician is familiar with the safety procedures which relate to the products processed by the system.

In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments returned to the manufacturer can be repaired only when certain procedures (see section "**Repair - Maintenance - Return**") are followed.

## 2 Technical data

### Type

### 8-channel adapter for pipette tips for hand controller

Permissible ambient temperature storage / operation	-10 °C to +60 °C/ +10 °C to +40 °C (14 °F to 140 °F/ 50 °F to 104 °F)
Autoclaving	121°C and 2 bar absolute (1 bar overpressure)
Adapter for	pipette tips 2 - 200 µl and 5 - 300 µl
Distance between nose cones	9 mm
Dimensions W x D x H approx.	80 mm x 42 mm x 118 mm
Weight approx.	80 g

## 2.1 Materials of the 8-channel adapter

### Wetted parts


Stainless steel, PP with and without glass fibre, PPS, PVDF, FKM, EPDM, silicone

### Not wetted parts

Stainless steel, PTFE with carbon, PP with and without glass fibre, PPS with glass fibre

We reserve the right to change technical specifications without prior notice.

## 2.2 System parts

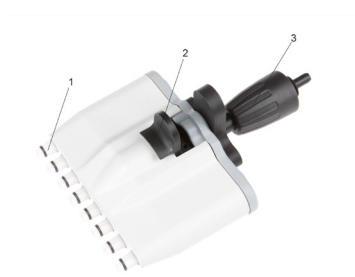Position	Description
1	nose cone with O-ring
2	ejector
3	screw cap

### 3 Use and operation

Unpack the equipment.


#### Scope of delivery

- 8-channel adapter (1)
- bench stand (2)
- mounting tool for nose cones (3)
- reagent reservoir / bowl for disinfectant (4)

#### NOTICE

The 8-channel tip adapter with ejector is in-tended to be connected to a Starlab hand controller.

#### 3.1 Assembling the adapter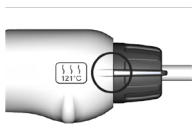

Ensure that there is no liquid in the tube to avoid risk of contamination.

Aspirate until the aspiration tube is empty and decontaminate if necessary.


Select continuous aspiration.

We recommend to switch off vacuum supply or to close connection to the vacuum supply.

Remove pipette if necessary.


- > **Attention:** Liquid in the tube may leak. Adopt suitable measures to prevent the release of dangerous, toxic, explosive, corrosive, noxious or polluting fluids, vapors and gases. In such cases, use an appropriate collecting and disposal system and take protective action for personnel, equipment and environment.

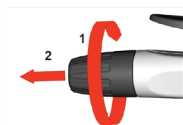Turn cap (1) and remove (2).

Pull existing adapter with hose out of the hand controller.  
Remove adapter from hose.

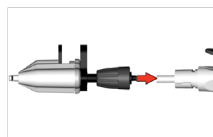
Attach 8-channel adapter to the hose.  
Pull the 8-channel adapter with hose back in the hand controller.

Align mark at the cap when reassembling the cap.


Screw cap hand tight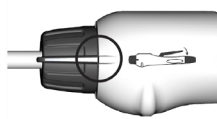

Set the hand controller to “Aspiration” if necessary.

### 3.2 Assembling the bench stand

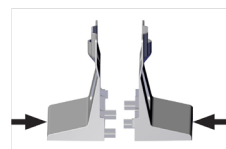Push together both halves of the stand correctly.

#### NOTICE

**Attention:** Once assembled, it is not possible to disassemble the bench stand without damage.

### 3.3 Prior to aspiration

Attach pipette tips.  
Ensure that the tips are securely seated. Align the adapter.  
The adapter can turn freely in both directions.

**Attention:** Pipette tips are disposables items. Never attach pipette tips that contain liquid due to risk of contamination and infection. Wear protective gloves!

### 3.4 Aspiration of liquids

Use only appropriate vessels. Dip pipette tip into the liquid.

**Setting aspiration** at the hand controller: Press the lever until the liquid has been aspirated. Release the lever to stop aspiration.

**Setting continuous aspiration** at the hand controller: In setting continuous aspiration it is not necessary to press the lever. Remove pipette from liquid to stop aspiration. If necessary set the adjusting knob to "Aspiration" to stop continuous aspiration.

### 3.5 Ejecting the tips

Hold the device over a suitable disposal container and press the ejector to the stop.

Never lay the instrument horizontal when the tips contain liquid. Liquid may enter and contaminate the instrument or the environment.

### 3.6 After aspiration

At the end of aspiration, continue to aspirate until there is no more liquid in the aspiration tubing and purge with cleaning or disinfectant agent (suck through) if necessary. Use reagent reservoir / bowl supplied for disinfectant if necessary.

Eject pipette tips.


**Attention:** Risk of contamination and infection. Wear protective gloves and eye protection! Dispose of pipette tips according to appropriate regulations.

We recommend storing the device without tips, placed upright in the bench stand supplied.

## 4 Cleaning and decontamination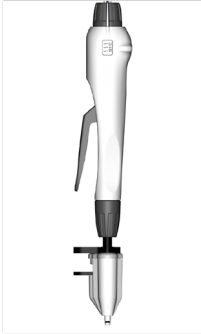

### Autoclaving


The hand controller with adapters and aspiration hose is designated for steam sterilization at 121 °C and 2 bar absolute (1 bar overpressure). Time of exposure according to DIN 58946  $t_e = 20$  minutes.

DO NOT exceed a temperature of 121 °C.

Comply with instructions for use of the manufacturer of the autoclave.

### NOTICE

It is the user's responsibility to ensure effective autoclaving.

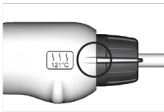
### NOTICE

Prior to autoclaving set the hand controller to "autoclaving".

Prior to autoclaving loosen the cap.

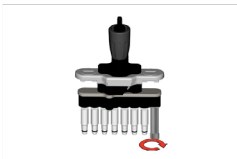### NOTICE

After autoclaving tighten the nose cones if necessary with the mounting tool supplied.

**NOTICE**


It is not possible to disassemble the 8-channel adapter completely.

It is possible to replace the nose cones, see section "Replacing the nose cones".

**NOTICE**

It is not possible to disassemble the hand controller completely.

It is possible to disassemble and replace the adapters, the screw cap and the aspiration hose, see instructions for use of the hand controller. If the housing opens slightly by mistake, e. g., if the handle is dropped, press the gap together by hand. Perform a function test.

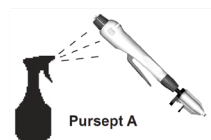**Chemical disinfection**

Wiping and spray disinfection is possible by using the germicide Pursept A.

**NOTICE**

Over time, discoloration and material changes (e.g., resiliency, elasticity/tightness, cracking) due to repeated steam sterilizations may occur.

## Important notes on use of disinfectants

Aggressive disinfectants which release chlorine or oxygen radicals, e.g., sodium hypochlorite (chlorine bleach) or peroxide compounds can corrode the materials.

- > Therefore, prolonged use of these disinfectants is strongly discouraged. After brief disinfecting operations, rinse thoroughly to avoid leaving residues of disinfectant.

The use of the disinfectant **Sekusept® Plus** (Manufacturer: Ecolab GmbH & Co OHG, Düsseldorf, Germany) in extensive inhouse testing did not cause any damage. Hence the disinfectant Sekusept® Plus is preferred. Comply with use and safety instructions of the manufacturer!

- > Using corrosive disinfectants may result in damage, malfunction and/or failure of the equipment.
- > Chlorine may damage the materials of the equipment or the vacuum supply.
- > Leaking of liquid from a damaged aspiration system may lead to exposure of personnel and material or to damage/destruction of wetted equipment or laboratory furniture.
- > For information about the compatibility with the materials, ask the manufacturer of the disinfectant.
- > The materials of the system are listed in the section "Technical data".

## 5 Accessories and spare parts

**Hand controller for Aspirator** **N2400-9001**

**Tubing for hand controller, 2.5 m** **N2400-9005**

## 5 Troubleshooting

Fault	Possible cause	Remedy
No aspiration at the hand controller. Pipette drips.	<ul style="list-style-type: none"> <li>› Adapter or aspiration hose defective?</li> <li>› Aspiration hose kinked?</li> <li>› Aspiration hose clogged or agglutinated</li> <li>› Poorly fitting pipette tips?</li> <li>› Leak in the system?</li> <li>› No vacuum?</li> </ul>	<ul style="list-style-type: none"> <li>☑ Replace adapter or aspiration hose.</li> <li>☑ Correct hose guide.</li> <li>☑ Remove clogged or agglutinated hose section or replace hose.</li> <li>☑ Attach tips securely or choose appropriate tips.</li> <li>☑ Check the system for leaks</li> <li>☑ Check the vacuum system</li> </ul>
Pipette loose in the adapter.	<ul style="list-style-type: none"> <li>› Adapter cone wet or greasy?</li> <li>› Tips with improper diameter selected?</li> <li>› Seal at the nose cone defective?</li> </ul>	<ul style="list-style-type: none"> <li>☑ Clean or dry nose cones.</li> <li>☑ Choose appropriate pipette or change tips.</li> <li>☑ Replace nose cone.</li> </ul>
Liquid aspirates slowly or not at all.	<ul style="list-style-type: none"> <li>› Pipette tip or nose cone clogged or damaged?</li> </ul>	<ul style="list-style-type: none"> <li>☑ Replace pipette tips or clean nose cones.</li> </ul>

## 7 Repair - Maintenance - Return

### IMPORTANT

In case of service requests contact your local Starlab sales representative or Starlab service technician. Contact details can be found "Starlab Service Hotline" or the back page.

# Return Form

Return code

Kindly enclose this form with every return!

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## Reason for return (Please tick)

Wrong delivery  Product Defect  Wrong item ordered

Delivery note N°: \_\_\_\_\_

## Details :

Returned Items			internal, filled in by the recipient	
Item N°	Lot N° / Serial N°	Quantity	Complete Quantity	Saleable Units
1.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
2.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
3.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
4.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
5.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
6.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
7.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
8.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:

## Return handling

Exchange item  Credit Note

## Number of boxes

Total number of returned boxes: \_\_\_\_\_

Was a decontamination of the returned goods necessary?  Yes  No

If yes, which was the nature of the contamination?

.....

Method of decontamination: .....

I confirm that the information on this form is accurate and the items detailed are safe to handle.

Date

Signature

**Starlab International GmbH**  
Neuer Hötligbaum 38  
22143 Hamburg, Germany  
E-Mail: [info@starlab.de](mailto:info@starlab.de)  
[www.starlabgroup.com](http://www.starlabgroup.com)  
Tel.: +49 (0)40 675 99 39 0  
Fax: +49 (0)40 675 9939 20

Authorised Managing Director  
**Klaus Ambos**

Registry Court  
Amtsgericht (first instance court)  
Hamburg  
Register number: HRB 122020  
VAT ID no. pursuant to  
§ 27 a German  
VAT Act DE 214 891 717

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**Starlab (UK), Ltd**
**Starlab S.r.l.**