This is your **hand controller** user manual.

USER MANUAL · BEDIENUNGSANLEITUNG · MODE D'EMPLOI

Passionate for science.

User Manual	3–26
Bedienungsanleitung	27–50
Mode d'emploi	51–74

1	Safety information	4
1.1	Important information	4
1.2	General information	5
1.3	Intended use	5
1.4	Setting up	5
1.5	Ambient conditions	6
1.6	Operating conditions	6
1.7	Safety during operation	7
1.8	Maintenance and repair	8
2	Technical data	9
2.1	Wetted parts	9
2.2	System parts	10
3	Use and operation	10
3.1	Assembling the wall holder	11
3.2	Adapters	13
3.3	Changing the adapter	14
3.4	Settings	15
3.5	Handling	16
3.6	Storage following use	18
4	Cleaning and decontaminating	18
5	Accessories and spare parts	20
6	Troubleshooting	21
7	Replacing the aspiration hose	22
8	Repair - Maintenance - Return	24
9	Return Form	25

1 Safety information

1.1 Important information

WARNING

- > **Keep this manual complete and accessible to personnel at all times!**
- > **Read this manual carefully before installing or operating the equipment.**
Observe the instructions contained in this manual.
- > **Do not modify the equipment without authorization.**

NOTICE

This manual is an integral part of the equipment described therein. It describes the safe and proper use of the equipment. Make operating personnel aware of dangers arising from the equipment and the used substances. Starlab cannot accept any liability for inappropriate use of the equipment or for damage from failure to follow instructions contained in this manual.

This manual is only to be used and distributed in its complete and original form. It is strictly the users' responsibility to check carefully the applicability of the guidance in this manual with respect to this product.

01/02/2023

The following signal word panels and safety symbols are used throughout this manual:

This is the **safety alert** symbol. It is used to alert you to potential personal injury hazards. Obey all safety messages that follow this symbol to avoid possible injury and death.

DANGER

- > **DANGER** indicates a hazardous situation which, if not avoided, will result in death or serious injury.

WARNING

- > **WARNING** indicates a hazardous situation which, if not avoided, could result in death or serious injury.

CAUTION

- > **CAUTION** indicates a hazardous situation which, if not avoided, could result in minor or moderate injury.

NOTICE

NOTICE is used to address practices not related to personal injury

1.2 General information

- › **Read and comply with this manual before installing or operating the equipment.**
- › **Before operating the equipment read and comply with the section "Cleaning and decontamination"!**

NOTICE

Remove all packing material, remove the product from its packing box, remove the protective covers from the inlet and outlet ports and keep. Inspect the equipment. If the equipment is damaged, notify the supplier within three days; state the item number of the product together with your order number and the supplier's invoice number. Retain all packing material for inspection.

Do not use the equipment if it is damaged.

1.3 Intended use

- › The device may only be used by well trained and skilled personnel.
- › The device may not be used for in vivo applications (e. g., not humans or animals).

NOTICE

Ensure that the equipment and all components are suitable for the intended application. Use the system only for **aspiration of liquids**.

1.4 Setting up

NOTICE

The Hand Controller is intended to be connected to a suitable suction system, e.g., Starlab Aspirator.

Comply with all **applicable and relevant safety requirements** (regulations and guidelines), **implement the required actions and adopt suitable safety measures**.

1.5 Ambient conditions

- > The systems are designed for **ambient temperatures** during operation between +10 °C and +40 °C.
Check the maximum temperatures.

1.6 Operating conditions

- > The devices are not suitable to aspirate:
 - **unstable substances**
 - substances which react explosively under impact (mechanical stress) and/or when being exposed to **elevated temperatures** without air,
 - **self inflammable** substances,
 - substances which are **inflammable without air** and
 - **explosive substance**
- > The devices have **no approval** for operation below ground.

- > The equipment is **not suitable** for aspiration of dust.

- > **Take into consideration interactions and chemical reactions of the aspirated media.**
- > Ensure that the materials of the wetted parts are compatible with the aspirated substances, see section "Technical data".

If aspirating **different substances**, it is recommended that the aspiration hose be purged prior to changing the media in order to rinse out residues and to avoid reactions of the substances with each other and with the materials.

1.7 Safety during operation

DANGER

- › Ensure that while operating the device no persons can be endangered.
- › Wear personal protection equipment.
- › Comply with requirements when using infectious liquids, pathogenic germs, toxic, radioactive or aggressive substances.
- › Prior to operation check all relevant safety and health requirements and determine the applicability of regulatory limitations prior to use if necessary.
- › Comply with material safety data sheets and notes on safe use of from the manufacturer.
- › Comply with the biosafety level of the laboratory.
- › Use appropriate safety work materials and methods if necessary, e.g., steam sterilization, sterilization indicators or germicides. For notes concerning the sterilization of wetted parts, see the "Technical Data" section and "Cleaning and decontaminating". It is the user's responsibility to ensure effective sterilization.
- › Never operate a defective or damaged device.

WARNING

- › Comply with applicable regulations when disposing of chemicals. Take into consideration that chemicals may be polluted.
- › Use only **OEM spare parts and accessories**; otherwise safety and performance of the equipment as well as the electromagnetic compatibility of the equipment might be reduced.
- › In case of leaks in the aspiration hose, aspirated substances might be released into the environment.
Comply especially with notes on operation and use.

NOTICE

Ensure that, in case of failure, the system will always return to a safe status. Provide appropriate protective measures (i.e., precautions which allow for the requirements of the respective application) even for the case of failure and **malfunction**.

1.8 Maintenance and repair

- > **Attention:** The device might be contaminated with noxious or otherwise dangerous chemicals that have been aspirated during operation. Ensure that the device is decontaminated before maintenance.

- > Take adequate precautions to protect people from the effects of dangerous substances if contamination has occurred. Wear appropriate safety-clothing when you come in contact with contaminated components.

Ensure that maintenance is carried out only by suitably trained technicians. Ensure that the maintenance technician is familiar with the safety procedures which relate to the products processed by the system.

In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments returned to the manufacturer can be repaired only when certain procedures (see section "**Repair - Maintenance - Return**") are followed.

2 Technical data


Type	Hand controller
Permissible ambient temperature storage / operation	-10 to +60 / +10 to +40 °C (14 to 140 / 50 to 104 °F)
Permissible relative atmospheric moisture during operation (no condensation)	30 to 85 %
Adapters for	Pipettes with outer diameter 5.5 - 8 mm, Pipette tips 2 - 200 µl and 50 - 1000 µl
Length of aspiration hose	2.5 m
Dimensions W x D x H approx. (without aspiration hose)	185 x 30 x 38 mm
Weight (with aspiration hose) approx.	125 g

2.1 Wetted parts

Components	Wetted parts
Aspiration hose	Silicone rubber
Adapter for pipettes	TPE / PP
Adapter for pipette tips	PP
Screw cap, adjusting knob, operating lever	PPS, glass fiber reinforced
Housing	PP, glass fiber reinforced


We reserve the right for technical modifications without prior notice.

2.2 System Parts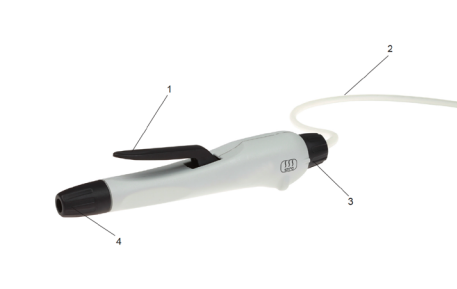


Position	Designation
1	operating lever
2	aspiration hose
3	adjusting knob
4	screw cap

3 Use and operation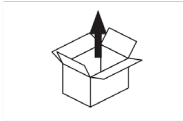


Unpack the equipment.

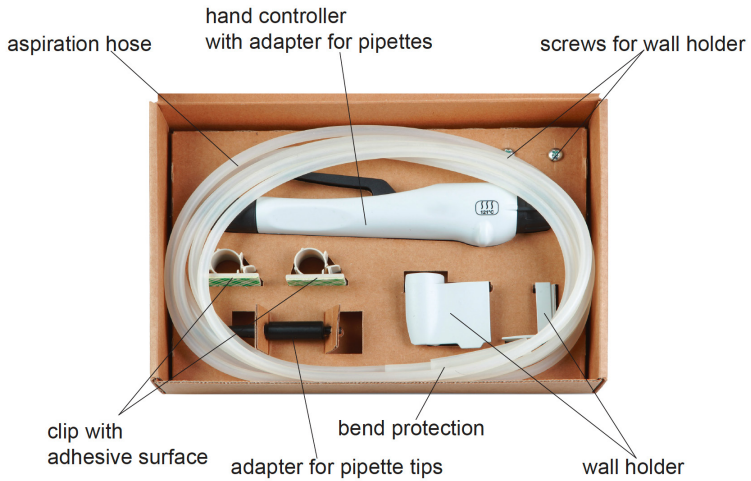
Attach the aspiration hose of the hand controller to a suitable aspiration system, e. g., Starlab Aspirator.

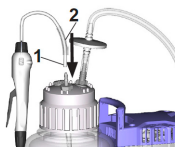
Attach the aspiration hose (1) first, then slide on the bend protection (2).

Enclosed are two clips with an adhesive surface. Affix these clips to the biosafety cabinet to guide the aspiration tubing.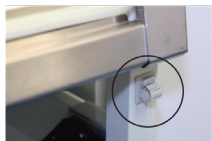

3.1 Assembling the wall holder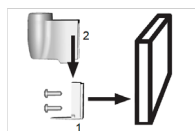

The surface of the wall has to be smooth, dry and free of grease.


Remove the protective foil from the adhesive tape at the wall holder part 1. Align part 1 with cover directed downward and press firmly on the wall for 1 minute.

The full load is achieved after 72 h at 20 °C.

Alternatively screw part 1 of the wall holder to the wall using the two screws enclosed in the delivery.

Assemble part 2 of the wall holder to part 1 and click into place.

Disassembling

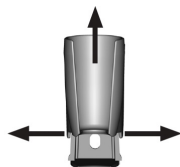Push both wings of the wall holder (part 2) apart and slide the part upwards.

NOTICE

When replacing the adhesive tape to fix the wall holder we recommend to use the 3M™ VHB™ high performance connection system 4932P. 3M™ and VHB™ are trade marks or registered trade marks of 3M.

Positioning the hand controller in the wall holder.


Alternatively, keep the hand controller in the table stand (N2400-9004) - see section "Accessories and spare parts".

3.2 Adapters

Two adapters are supplied with the hand controller: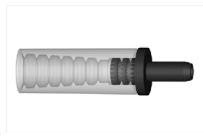

Adapter for pipettes with outer diameter of 5.5 - 8.0 mm

- Pasteur pipette
- Aspiration pipette
- Graduated pipettes

www.starlab.click/aspirationpipettes

Adapter for pipette tips

- Pipette tips 2 - 200 μ l
- Pipette tips 50 - 1000 μ l


TipOne®

www.starlab.click/tipone

A single-channel adapter for pipette tips 2 - 200 μ l with tip ejector, and an 8-channel tip adapter with ejector for pipette tips 2 - 200 μ l and 50 - 1000 μ l are also available to purchase separately as accessories. See section "Accessories and spare parts".

3.3 Changing the adapter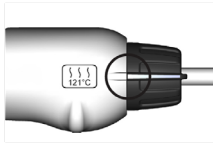

Ensure that there is no liquid in the tube to avoid risk of contamination.

Aspirate until the aspiration tube is empty and decontaminate if necessary.


Select continuous aspiration.

We recommend to switch off vacuum supply or to close connection to the vacuum supply.


Remove pipette if necessary.


- Attention: Liquid in the tube may leak. Adopt suitable measures to prevent the release of dangerous, toxic, explosive, corrosive, noxious or polluting fluids, vapors and gases. In such cases, use an appropriate collecting and disposal system and take protective action for personnel, equipment and environment.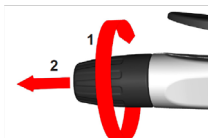

Turn cap (1) and remove (2).

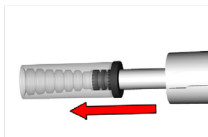Pull adapter with hose out of the hand controller.

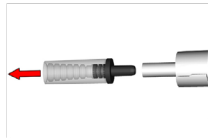
Remove adapter from hose.


Attach new adapter to the hose.

Pull the adapter with hose back into the hand controller.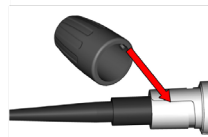


Align mark at the cap when reassembling the cap.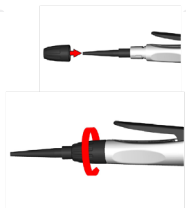


Attach cap over the adapter to the hand controller.

Screw cap.


3.4 Settings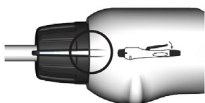

Setting "Aspiration".

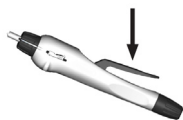
When the hand controller is set to "Aspiration" press the lever to aspirate.

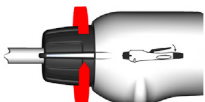The adjustment knob can be turned through 360°, but the knob is not removable.

Setting "Continuous aspiration" and "Autoclaving".

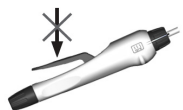Continuous aspiration / Autoclaving

The setting "continuous aspiration / autoclaving" is designated to aspirate larger amounts of liquid. The hand controller aspirates without pressing the lever.

Select this setting also for autoclaving, see section "Cleaning and decontamination".

3.5 Handling

Hold the hand controller so it is comfortable for working, see examples below:

Prior to aspiration

Choose the appropriate adapter and carefully insert the pipette into the adapter.

To ease the assembly, loosen the screw cap prior to inserting the pipette and then retighten cap if necessary.

Attention: Never insert a filled pipette due to risk of contamination and infection. Wear protective gloves!

Aspiration of liquid

Use only appropriate vessels. Dip pipette tip into the liquid.

Setting aspiration: Press the lever until the liquid has been aspirated. Release the lever to stop aspiration

Setting continuous aspiration: When setting continuous aspiration it is not necessary to press the lever. Remove pipette from liquid to stop aspiration. If necessary, set the adjusting knob to “Aspiration” to stop continuous aspiration.

After aspiration

At the end of aspiration, continue to aspirate until there is no more liquid in the aspiration tubing and purge (suck through) with cleaning or disinfectant agent if necessary.

Carefully pull pipette out of the adapter. To ease the disassembly, loosen the screw cap prior to removing the pipette and then retighten cap if necessary.

Attention: Risk of contamination and infection. Wear protective gloves and eye protection! Dispose of pipettes according to appropriate regulations.

- > **Attention:** When removing the hand controller, liquid in the tube may leak - especially when using systems without couplings! At the end of aspiration, continue to aspirate until there is no more liquid in the aspiration tubing and purge with cleaning or disinfectant agent if necessary.


Adopt suitable measures to prevent the release of dangerous, toxic, explosive, corrosive, noxious or polluting fluids, vapors and gases. In such cases, use an appropriate collecting and disposal system and take protective action for personnel, equipment and environment.

3.6 Storage following use

NOTICE

At the end of aspiration, continue to aspirate until there is no more liquid in the aspiration tubing and purge with cleaning or disinfectant agent if necessary.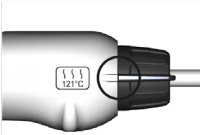

Set the hand controller to continuous aspiration to avoid a possible agglutination of the aspiration hose.

To avoid a permanent load on the vacuum supply, we recommend to switch off vacuum supply or to close connection to the vacuum supply.

Store the device in dry conditions.

4 Cleaning and decontaminating

Autoclaving

The hand controller with adapters and aspiration hose is designated for steam sterilization at 121 °C and 2 bar absolute (1 bar overpressure).

Time of exposure according to DIN 58946 $t_e = 20$ minutes.



DO NOT exceed a temperature of 121 °C. Comply with instructions for use of the manufacturer of the autoclave.

NOTICE

It is the user's responsibility to ensure effective autoclaving.

NOTICE

Prior to autoclaving set the hand controller to "autoclaving".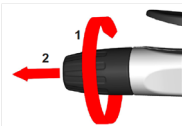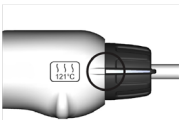

Prior to autoclaving turn cap (1) and remove (2)


NOTICE

It is not possible to disassemble the hand controller completely.

It is possible to disassemble and replace the adapters, the screw cap and the aspiration hose, see “Accessories and spare parts”.

If the housing opens slightly by mistake, e. g., if the handle is dropped, press the gap together by hand. Perform a function test.

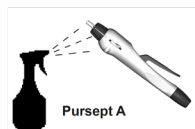**Chemical disinfection**

Wiping or spray disinfection of the hand controller is possible by using the germicide Pursept A.

NOTICE

Over time, discoloration and material changes (e.g., resiliency, elasticity/tightness, cracking) due to repeated steam sterilizations may occur.

Important notes on use of disinfectants

Aggressive disinfectants which release chlorine or oxygen radicals, e.g., sodium hypochlorite (chlorine bleach) or peroxide compounds can corrode the materials. The use of the disinfectant **Sekusept® Plus** (Manufacturer: Ecolab GmbH & Co OHG, Düsseldorf, Germany) in extensive inhouse testing did not cause any damage. Hence the disinfectant Sekusept® Plus is preferred. Comply with use and safety instructions of the manufacturer!

CAUTION

- > Using corrosive disinfectants may result in damage, malfunction and/or failure of the equipment.
- > Chlorine may damage the materials of the equipment or the vacuum supply.
- > Leaking of liquid from a damaged aspiration hose may lead to exposure of personnel and material or to damage/destruction of wetted equipment or laboratory furniture.
- > For information about the compatibility with the materials, ask the manufacturer of the disinfectant.
- > The wetted materials of the system are listed in the section “Technical data”.

5 Accessories and spare parts

Hand controller for Aspirator

N2400-9001

Tubing for hand controller, 2.5 m

N2400-9005

Connection Kit for 2nd an additional hand controller

N2400-9006

(to be assembled at a Aspirator, does not include hand controller or quick coupling)

Quick coupling hand controller to bottle

N2400-9008

(suitable as extension kit for a 2nd an additional hand controller)

Quick coupling made of PVDF, with adapter to connect a hand controller to a collection bottle. When disconnected, the collection bottle closes vacuum tight.

Attention: The quick coupling accessory set is not suitable for use with sodium hypochlorite (chlorine bleach).

Stand for hand controller

N2400-9004

Single channel adapter for pipette tips for hand controller

N2400-9002

for pipette tips 2 - 200 μ l

8-channel adapter for pipette tips

N2400-9003


for pipette tips 2 - 200 μ l and 5 - 300 μ l

6 Troubleshooting

Fault	Possible cause	Remedy
No aspiration at the hand controller, pipette drips.	<ul style="list-style-type: none"> › Adapter or aspiration tube defective? › Aspiration hose clogged or agglutinated? › Poorly fitting pipette or tips? › Leak in the system? › No vacuum? 	<ul style="list-style-type: none"> ☑ Replace adapter with aspiration hose. ☑ Remove clogged or agglutinated hose section or replace hose. ☑ Connect pipette securely or choose an appropriate adapter. ☑ Check the system for leaks. ☑ Check the vacuum system.
Pipette loose in the adapter.	<ul style="list-style-type: none"> › Adapter wet or greasy? › Pipette with improper diameter selected? 	<ul style="list-style-type: none"> ☑ Clean or replace adapter. ☑ Choose appropriate pipette or change adapter.
Broken pipette in the adapter.	<ul style="list-style-type: none"> › The pipette was damaged or forced into the adapter? 	<ul style="list-style-type: none"> ☑ Remove the glass splinters by means of forceps, replace adapter and aspiration hose if necessary.

7 Replacing the aspiration hose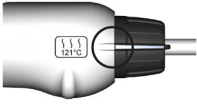

Ensure that there is no liquid in the tube to avoid risk of contamination.

Aspirate until the aspiration tube is empty and decontaminate if necessary.


Select continuous aspiration.

Switch off vacuum supply or close connection to the vacuum supply.


Remove pipette if necessary.

- > **Attention:** Liquid in the tube may leak. Adopt suitable measures to prevent the release of dangerous, toxic, explosive, corrosive, noxious or polluting fluids, vapors and gases. In such cases, use an appropriate collecting and disposal system and take protective action for personnel, equipment and environment.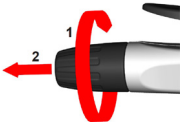

Turn cap (1) and remove (2).

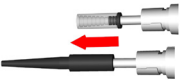Pull adapter with hose out of the hand controller.

Remove adapter from hose.

Pull aspiration hose out of the hand controller.


Insert new aspiration hose into the hand controller and push through.


Attach adapter to the hose.

Pull the adapter with hose back into the hand controller.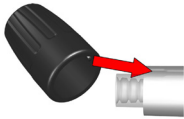

Align mark at the cap when reassembling the cap.

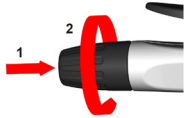Attach cap over the adapter to the hand controller.
Screw cap.

8 Repair - Maintenance - Return

IMPORTANT

In case of service requests contact your local Starlab sales representative or Starlab service technician. Contact details can be found "Starlab Service Hotline" or the back page.

Return Form

Return code

Kindly enclose this form with every return!

Recipient	
STARLAB INTERNATIONAL GmbH Neuer Höltigbaum 38 D-22143 Hamburg	
<input type="checkbox"/> Warehouse	<input type="checkbox"/> Technical Service

Sender	
Company:	
Name:	
Address:	
Customer N°:	
Tel. / E-Mail:	

Reason for return (Please tick)	
<input type="checkbox"/> Wrong delivery	<input type="checkbox"/> Product Defect
<input type="checkbox"/> Wrong item ordered	

Delivery note N°:

Details :

Returned Items			internal, filled in by the recipient	
Item N°	Lot N° / Serial N°	Quantity	Complete Quantity	Saleable Units
1.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
2.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
3.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
4.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
5.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
6.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
7.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:
8.			<input type="checkbox"/> yes <input type="checkbox"/> no	Units:

Return handling	
Exchange item	<input type="checkbox"/>
Credit Note	<input type="checkbox"/>

Number of boxes
Total number of returned boxes:

Was a decontamination of the returned goods necessary? Yes No

If yes, which was the nature of the contamination?
.....

Method of decontamination:

I confirm that the information on this form is accurate and the items detailed are safe to handle.

Date _____ Signature _____

Starlab International GmbH
Neuer Höltigbaum 38
22143 Hamburg, Germany
E-Mail: info@starlab.de
www.starlabgroup.com
Tel.: +49 (0)40 675 99 39 0
Fax: +49 (0)40 675 9939 20

Authorised Managing Director
Klaus Ambos

Registry Court
Amtsgericht (first instance court)
Hamburg
Register number: HRB 122020
VAT ID no. pursuant to
§ 27 a German
VAT Act DE 214 891 717

User Manual	3–26
Bedienungsanleitung	27–50
Mode d'emploi	51–74

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Starlab (UK), Ltd
Starlab S.r.l.