Strength and Durability
for reliable protection

StarGuard® Protect

In Compliance with:

Personal Protective Equipment EU Regulation 2016/425, PPE Cat. III, EN ISO 21420:2020, EN ISO 374-1:2016 +A1:2018, EN 16523-1:2015, EN 374-2:2019 Level 3 (AQL 0.65), EN 374-4:2019, EN ISO 374-5:2016, EN 455 parts 1, 2, 3 and 4, Protection against Bacteria and Fungi: PASS, Protection against Viruses: PASS

The following is a link to view the Declaration of Conformity: <https://www.starlab.click/doc-starguard-protect>

Notified Body 2777: SATRA Technology Europe Limited. Bracetown Business Park. Clonee. D15YN2P. Republic of Ireland. This NB carried out all EU Type Examination and ongoing conformity.

AQL
0.65

Available sizes:

100 per box

100 per box

100 per box

100 per box

100 per box

CAUTION:

Protection is only for the hand. Before donning inspect the integrity of the glove to assure there are no tears, cracks, holes, bumps color or other visual defects and dispose of any gloves presenting such defects. Upon request a list of substances known to cause allergies will be provided to users. Requests may be submitted to info@starlab.de

WARNING:

EN 374-4:2013 Degradation levels indicate the change in puncture resistance of the gloves after exposure to the challenge chemical. This information does not reflect the actual duration of protection in the workplace and the differentiation between mixtures and pure chemicals.

The chemical resistance has been assessed under laboratory conditions from samples taken from the palm only (except in cases where the glove is equal to or over 400mm - where the cuff is tested also) and relates only to the chemical tested. It can be different if the chemical is used in a mixture.

It is recommended to check that the gloves are suitable for the intended use because the conditions at the workplace may differ from the type test depending on temperature, abrasion and degradation.

When used, protective gloves may provide less resistance to the dangerous chemical due to the changes in the physical properties. Movements, snagging, rubbing, degradation caused by the chemical contact etc may reduce the actual use time significantly. For corrosive chemicals, degradation can be the most important factor to consider in selection of chemical resistant gloves.

The penetration resistance has been assessed under laboratory conditions and relates only to the intended specimen.

Chemical	Performance Level	Degradation % change
	(EN ISO 374-1: 2016)	(EN 374-4: 2013)
(J) n-Heptan	3	38.3 %
(K) Sodium Hydroxide 40%	6	-8.4 %
(P) Hydrogen Peroxide 30%(R)	4	31.6 %
(T) Formaldehyde 37%	6	6.2 %

Performance Levels	Breakthrough Times (Minutes)
Level 1	> 10 minutes
Level 2	> 30 minutes
Level 3	> 60 minutes
Level 4	> 120 minutes
Level 5	> 240 minutes
Level 6	> 480 minutes

Proper donning and doffing of gloves are essential in protecting the user and safe guarding the integrity of the gloves being worn. The following are suggested steps to properly wear and remove ambidextrous gloves:

Donning

Step 1. After inspection of the glove, Insert all five fingers into the cuff of the glove, and pull the cuff over your wrist until the glove is properly in place. Assure the cuff is pulled up completely to avoid any substances entering the glove.

Step 2. Check that the glove's fit is secure around the fingers and palm. Check the cuff, which should have a secure fit around your wrist.

Doffing

Step 1. When removing gloves, Using a gloved hand, take hold of the outside edge of your glove near your wrist with two fingers, careful to not touch your skin. Peel the glove away from your hand being sure to turn it inside out as you do. Grasp the removed glove in your hand that is still gloved.

Step 2. Slide your ungloved finger under the wrist of the remaining glove. Using your finger that's inside the glove, peel the remaining glove away, continuing to grasp the removed glove with your gloved hand. Once glove is removed properly dispose of the gloves.

Strength and Durability
for reliable protection

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GB **Caution:** These nitrile gloves are not intended to provide a barrier against all chemicals and are not suitable for contact with aggressive chemicals requiring a specialised heavy-duty protection glove. Quoted breakthrough times for chemicals based on laboratory tests may differ from actual times at the workplace due to factors such as abrasion and temperature degradation. Compounds used in the manufacture of this product may cause allergic reactions in some users. In case of an allergic reaction, discontinue use and seek medical advice. **Instructions for use:** For single and transient use only. Do not reuse. Immediately discontinue use following damage. Dispose of according to your local regulations. **Transport and storage:** This product to be stored in carton during transportation. Store product in a cool (<40°C), dry and well-ventilated location. Avoid exposure to sunlight, fluorescent lighting and x-rays.

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Starlab UK Ltd., 5 Tanners Drive,
Starlab International GmbH