



This is your **3D Rocking Shaker** pro locking hose replacement instruction.

1

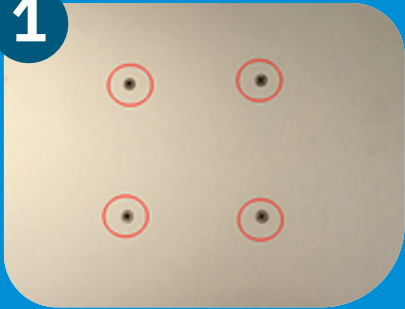


Remove the 4 screws of the shaking plate marked by red circles and keep them well

2

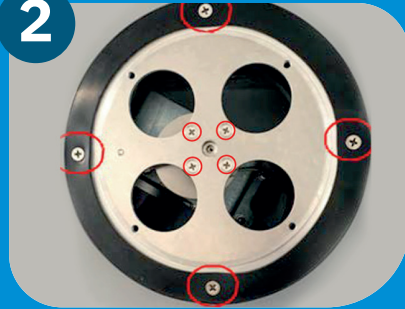


Remove the 8 screws marked by red circles and keep them well

3

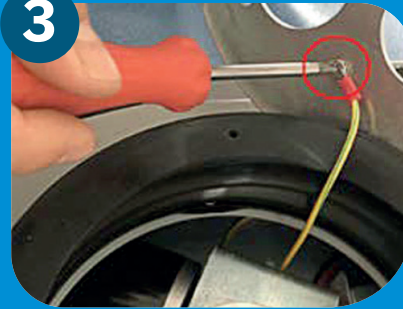


Remove the screw marked by red circle, separate ground wire and metal plate

4

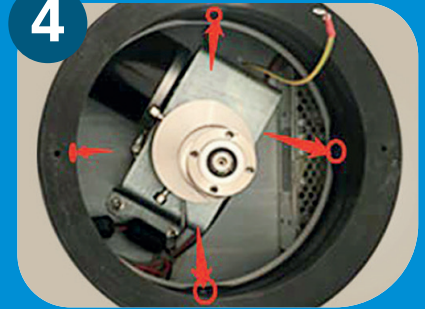


Remove 4 screws which marked with red circle that inside of the rubber locking hose

5

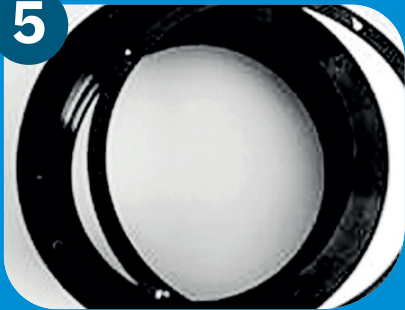


Remove the metal ring from the old cuff and set this into the new one. Fix the 4 screws as in step 4

6

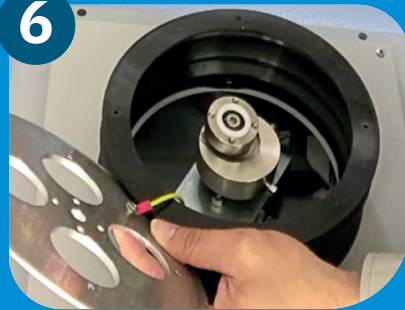

Connect the ground wire with metal flange as in step 3

7

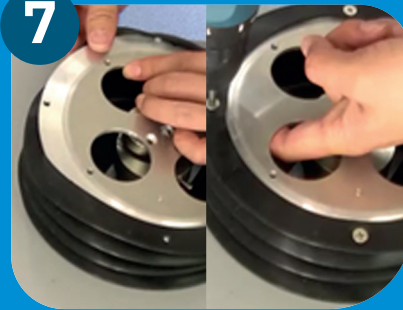


Put the flange into the rubber locking hose and fix it with screws (outer screws as marked in step 2)

8



Fix the middle 4 screws

9

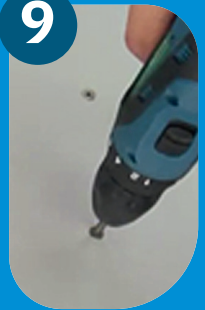

Install the shaking plate and fix it with screws as in step 1