

The Starlab Vortex Mixer.

Good vibrations. **Great results.**



Love at first vibe.

INNOVATIVE DESIGN

The Starlab Vortex Mixer combines stylish design with an innovative, easy to use touchscreen for quick speed and time settings.



A screen that will touch you.



EASY TO USE GLOVE COMPATIBLE TOUCHSCREEN

Precise speed settings. Quick select presets for speed and time. Speed and time custom settings. One-touch start / stop in continuous mode.

Great features. Good vibrations.

THE STARLAB VORTEX MIXER

With its compact and stylish design, the Starlab Vortex Mixer is the perfect bench partner for your lab. The unique touchscreen provides precise speed settings for the gentle mixing of samples right up to vigorous vortex mixing and the resuspending of cells or chemical pellets.

- Glove compatible touch screen
- Precise vortexing speeds, from 300 to 3,500 rpm
- One-touch pre-set options for speed / time
- Touch or continuous operation
- End of run audible alarm
- Power-saving sleep mode
- Optional accessories



WATCH VIDEO

Starlab Vortex Mixer.

ANTI-SLIP FEET
provides stability on the
bench during mixing

TOUCH MODE

Minimal pressure required
to maintain vibration.
Mix single or multiple
tubes up to 3,500 rpm.

SPEED RANGE
300 – 3,500 RPM

**GLOVE
COMPATIBLE
TOUCHSCREEN**
for precise speed and
time settings

**POWER-SAVING
SLEEP MODE**

**SMOOTH, QUIET
OPERATION**



OPTIONAL ACCESSORIES



Adapter for microplates

Mix one plate with a SBS/SLAS footprint with speeds up to 1,500 rpm




Adapter for 1.5/2.0 ml tubes

Mix up to 15 tubes at the same time with speeds up to 1,500 rpm

THE EASY TO USE TOUCHSCREEN

Tap speed or time to access the preset or custom settings. Once set, speed and duration is visible on the home screen. Speed and time settings can be changed during operation.

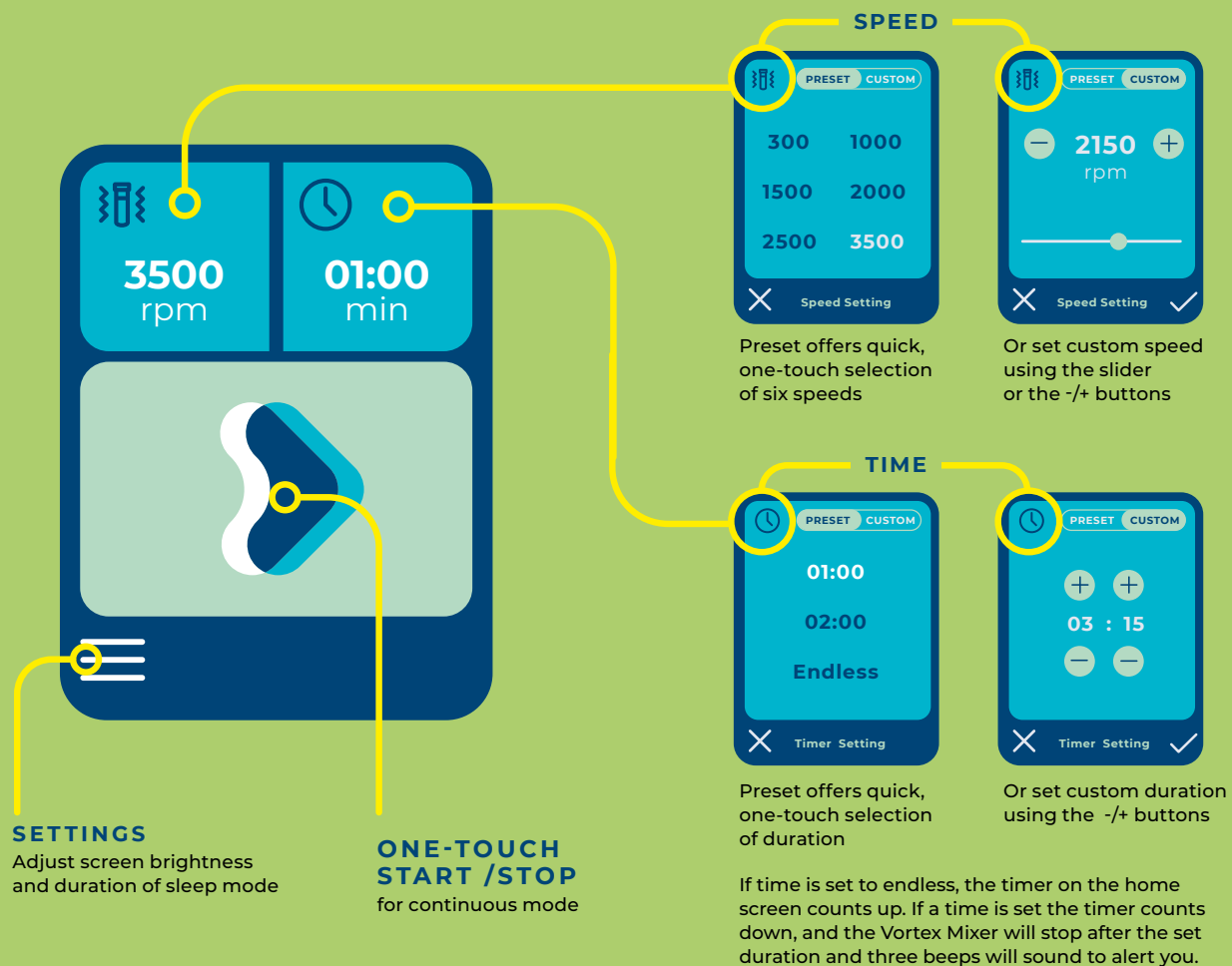

Screens shown are with the Standard Attachment installed. With the optional adapters attached, the maximum speed defaults to 1,500 rpm.
NOTE: Touch mode is not available when using the optional adapters. Microplate adapter is not suitable for deepwell plates.

A close-up photograph of a Starlab Vortex Mixer. The device is a dark, rectangular unit with a grid of small, raised circular buttons on its top surface. A central circular button is slightly larger and more prominent. The device is resting on a light-colored surface, and a portion of a blue and white striped tube is visible in the foreground. The background is a solid light blue color with a decorative white wavy line at the top.

Bringing your lab good vibes only.

TOUCH MODE

Minimal pressure required to maintain vibration.
Mix single or multiple tubes up to 3,500 rpm.
The large surface area allows different shaped tubes and containers to be vortexed, whilst the recessed centre provides the benefits associated with traditional, cup-style attachments.

Subscribe to the vibe.

ORDER INFO & TECHNICAL SPECIFICATIONS



Description	Article No.
 Vortex Mixer ¹⁾	S8010-0000
Accessories	
 Replacement Standard Attachment	S8010-0010
 Adapter for 1.5 / 2.0 ml tubes	S8010-0011
 Adapter for microplates	S8010-0012
 Replacement power adapter with four exchangeable plugs, 24 V, 1 A for EU, UK, US & AUS	S8010-5012

¹⁾ Vortex Mixer is supplied with one standard attachment and a power adapter with four exchangeable plugs.
2-year warranty.

Technical specifications	
Speed range	Standard attachment: 300 to 3,500 rpm Adapter for 1.5/2.0 ml tubes: 300 to 1,500 rpm Adapter for microplates: 300 to 1,500 rpm
Speed increments	50 rpm
Timing capacity	1 second to 99:59 minutes
Type of movement	Orbital
Operation modes	Two: touch and continuous
Noise level	≤ 55 dBA
USB output	5 VDC 1 A
Permissible ambient temperature	2 °C – 40 °C
Permissible relative humidity	≤ 80 %
Power supply	100 – 240 V, 50 Hz – 60 Hz
Unit input	24 V DC 12 W
Weight	≤ 2.8 kg
Dimensions (W x D x H)	200 x 130 x 103 mm



 starlab.click/shop-vortex



T: +49 (0)40 675 99 39 0
info@starlab.de
www.starlab.de



T: +44 (0)1908 283800
info@starlab.co.uk
www.starlab.co.uk



T: +33 (0)1 60 13 71 70
info@starlab.fr
www.starlab.fr



For our international
dealers, please go to
starlab.click/**distributors**



T: +39 (0)2 7020 1040
info@starlab.it
www.starlab.it



T: +41 56 6645 980
info@starlab.ch
www.starlab.ch



THE STARLAB VORTEX MIXER. GOOD VIBRATIONS. GREAT RESULTS.

With its compact and stylish design, the Starlab Vortex Mixer is the perfect bench partner for your lab.

The unique touchscreen provides precise speed settings for the gentle mixing of samples right up to vigorous vortex mixing and the resuspending of cells or chemical pellets.

 starlab.click/shop-vortex