
This is your battery safety information.

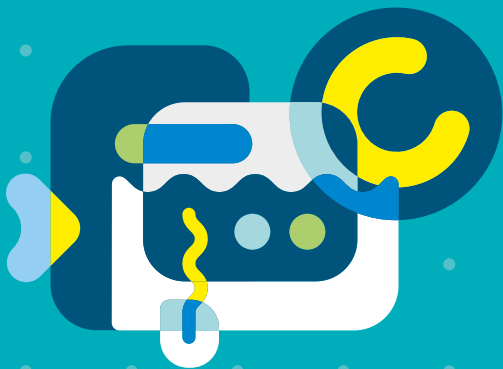Passionate for science.

Safety instructions	3–7
Sicherheitsbestimmungen	9–13
Règles de sécurité	15–19
Norme di sicurezza	21–25
Normas de seguridad	27–31
Normas de segurança	33–37
1 Battery and accumulator safety	4
2 Markings	5
3 Batteries	5
4 Accumulators	6

1. Battery and accumulator safety

When used correctly, batteries and accumulators are a safe and reliable source of energy. However, improper use may result in spills, vapors or, in extreme cases, explosions and/or fires. Therefore, observe the warnings in this document.

The warnings apply to:

- › ErgoOne® E
- › ErgoOne® Fast
- › GuardOne®

Starlab reacts to Regulation (EU) 2023/1542 (Battery Regulation) and thus helps to ensure the safe use of the batteries and rechargeable batteries by users. The user uses the batteries and rechargeable batteries at his laboratory workstation. The batteries and rechargeable batteries are replaced or charged there and the following warnings apply in this context.

According to the regulation, this document is valid until 02/18/2027.

2. Markings

	With this mark, we confirm that the product complies with the requirements set out in the EC Directives and has been subjected to the specified testing procedures.
	The device or the battery should be disposed of properly.
	This electrical instrument may not be disposed of in household waste.

3. Batteries

Handling

- › Do not insert batteries incorrectly. Observe the polarity marking on the battery and the device.
- › Do not short-circuit batteries.
- › Do not use batteries of different types or manufacturers together.
- › Keep batteries away from fire, heat and direct sunlight.
- › Do not disassemble or damage the batteries.
- › Do not modify the batteries.
- › Remove empty batteries from the laboratory equipment immediately and discard them.

Storage

- › Store unused batteries in their original packaging. If you store unpackaged batteries together, a short circuit and heat build-up may occur.
- › Store and store batteries in their original packaging in a well-ventilated, dry and cool place.
- › Protect batteries from direct sunlight and moisture.
- › Batteries are normally stored at temperatures between -20 °C and + 40 °C and relative humidity of 5% to 95%.

4. Accumulators

Handling

- › To ensure that the battery is used correctly, read the specifications carefully.
- › Keep the battery away from fire, heat and direct sunlight.
- › Do not connect the battery to a charger or device with reversed positive and negative terminals.
- › Do not short-circuit the battery.
- › Do not expose the battery to physical shock or vibration.
- › Do not disassemble the battery.
- › Do not change the battery.
- › Charge the battery in a suitable charger.
- › Do not use modified, damaged or battery-unsuitable chargers.
- › Do not discharge the battery if the ambient temperature is outside the temperature limits specified in the instructions for use.
- › Protect the battery from moisture or moisture.

- › Never use the battery if it is damaged. This could be, for example: Decreased performance, damaged and deformed „inflated“ housings, fluid spilled, and battery management system error messages for lithium-ion batteries on the device.
- › Dispose of the battery in accordance with local regulations.

Storage of NiMH batteries

- › Store NiMH batteries at temperatures between +10°C and +40 °C.

Storage of lithium-ion batteries

- › The lithium-ion battery should be stored in a cool, dry and well-ventilated place. Protect the battery from fire, heat and direct sunlight.
- › Store lithium-ion batteries at temperatures between -0°C and +35 °C.

Safety instructions	3–7
Sicherheitsbestimmungen	9–13
Règles de sécurité	15–19
Norme di sicurezza	21–25
Normas de seguridad	27–31
Normas de segurança	33–37

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Starlab (UK), Ltd
Starlab S.r.l.