# This is your **GuardOne®** user manual.

USER MANUAL · BEDIENUNGSANLEITUNG  
MODE D'EMPLOI · ISTRUZIONI PER L'USO

Passionate for science.

# EC Declaration of Conformity

The CE marking affixed to the equipment indicates that the  
**STARLAB GuardOne® Workstation for process protection**  
**S8040-0000:** LFG032SL (Laminar Flow GuardOne 32 inch)  
**S8040-0001:** LFG048SL (Laminar Flow GuardOne 48 inch)  
**S8040-1000:** DAG032SL (Dead Air GuardOne 32 inch)  
**S8040-1001:** DAG048SL (Dead Air GuardOne 48 inch)  
meets the requirements of the following directives:

## 2014/35/EU Low Voltage Directive (LVD)

EN 61010-1:2010 (Safety requirements for electrical equipment for measurement, control, and laboratory use - General requirements)

## 2014/30/EU EMC Directive

EN 55011:2009+A1:2010 (ISM equipment - Radio-frequency disturbance characteristics - Limits and methods of measurement)

EN 61326-1:2013 (Electrical equipment for measurement, control and laboratory use - EMC requirements - General requirements)

## Product-specific Requirements

IEC 62471:2006 (Photobiological safety of lamps and lamp systems)

ISO 14644-1:2015 Class 5 (Cleanrooms and associated controlled environments - Classification of air cleanliness by particle concentration)

## Additional Legal Requirements

2011/65/EU (RoHS – Directive on the restriction of the use of certain hazardous substances in electrical and electronic equipment)

EN 50581: 2012 (Technical documentation for the assessment of electrical and electronic products with respect to the restriction of hazardous substances)

2012/19/EU (WEEE – Directive on waste electrical and electronic equipment)

EN 50419:2006 (Marking of electrical and electronic equipment in accordance with Directive WEEE)

EC 1907/2006 (Directive on REACH)

94/62/EC (Directive on packaging and packaging waste)

## Additional Requirements

EN ISO 14971:2012 (Medical devices – Application of risk management to medical devices)

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### Product-specific Requirements

IEC 62471:2006 (Photobiological safety of lamps and lamp systems)  
ISO 14644-1:2015 Class 5 (Cleanrooms and associated controlled environments - Classification of air cleanliness by particle concentration)

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### 2014/35/EU Low Voltage Directive

UK Electrical Equipment (Safety) Regulations 2016

### 2014/30/EU Electromagnetic Compatibility Directive

UK Electromagnetic Compatibility Regulations 2016

### 2012/19/EU WEEE

### 2011/65/EU RoHS Directive

UK Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

### EC 1907/2006 Directive on REACH

### 94/62/EC Packaging

### Additional requirements:

EN 50581: 2012  
EN 50419:2006  
EN 61010-1:2010  
EN 55011:2009+A1:2010  
EN 61326-1:2013

Certified by:

Klaus Ambos  
CEO STARLAB International GmbH

Hamburg, 28.01.2022

Passionate for science.

Starlab International GmbH  
Neuer Höltingbaum 38  
22143 Hamburg, Germany  
E-Mail: [info@starlab.de](mailto:info@starlab.de)  
[www.starlabgroup.com](http://www.starlabgroup.com)  
Tel.: +49 (0)40 675 99 39 0  
Fax: +49 (0)40 675 9939 20

Authorised Managing Director  
Klaus Ambos

Registry Court  
Amtsgericht [first instance court]  
Hamburg  
Register number: HRB 122020  
VAT ID no. pursuant to  
§ 27 a German  
VAT Act DE 214 891 717

GuardOne® · Manual · 3

1 Overview of the GuardOne® Workstation

1.1 Setup and Function

Setup

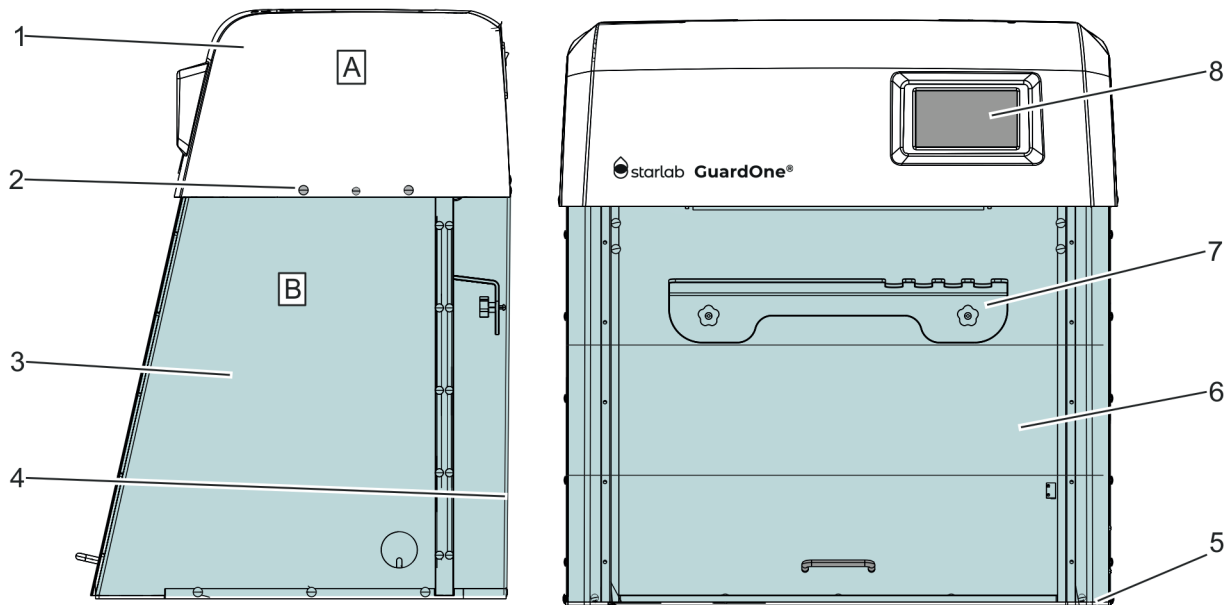
Fig. 1: GuardOne® Workstation

A	Top housing unit	4	Back panel
B	PC (Polycarbonate) glass chamber	5	Base plate
1	Top housing with pre-installed pre-filter and HEPA filter	6	Pre-installed front door with two folding segments
2	Plastic rivets	7	Multifunctional shelf
3	Side panel	8	Touchscreen display with graphical user interface

Models

There are two different kinds of models of the GuardOne® Workstation, a Dead Air and a Laminar Flow model. Each of them you can order in two sizes.

The two sizes of the GuardOne® Workstation are:

- > 32": In this size, the GuardOne® Workstation has one UV lamp and, assuming it is the Laminar Flow model, one pre- and one HEPA filter.
- > 48": In this size, the GuardOne® Workstation has two UV lamps and, assuming it is the Laminar Flow model, two blower motors, two pre- and two HEPA filters.

The dimensions can be found Chapter 10 "Specifications".

Dead Air

- > This type of the GuardOne® Workstation lacks the airflow function to create a clean air environment. Nonetheless, you can decontaminate the working area using the UV lamps.

Laminar Flow

- > Depending on its size, the Laminar Flow model is delivered with one (in 32" version) or two (in 48" version) blower motors and one or two pre- and one or two HEPA filters. In addition, one or two UV lamps are included.
- > This type of GuardOne® Workstation can create a clean air environment using the airflow function and it decontaminates the working area using the UV lamps.

**i** In this operating manual, all functions with regards to the pre- and HEPA filter as well as the UV lamp are described. If you have a Dead Air model, you can ignore the instructions for operating, replacing, maintaining and troubleshooting the filters.

**i** This operating manual provides the instructions for the Laminar Flow, size 32" with one HEPA and prefilter and one UV lamp. If you have a model Laminar Flow or Dead Air of size 48", follow the instructions for installation and maintenance of HEPA filters, UV lamps and pre-filters twice to install and replace the lamps and filters.

## Function

The GuardOne® Workstation is a containment equipment using air filtration technology and UV irradiation to create a working environment in which the risk of cross-contamination under standard laboratory working conditions is minimized.

To achieve this, the GuardOne® Workstation uses the following components:

- **Create a clean air environment**  
Combination of low-noise blower motor, standard HEPA filter and pre-filter.
- **Decontamination and sterilization of the working area**  
UV light with a fixed wavelength of 254 nm.

## Starlab Service Hotline

For further information, contact your local Starlab service or your local Starlab sales representative on [www.starlabgroup.com](http://www.starlabgroup.com).

1.2 Scope of Delivery

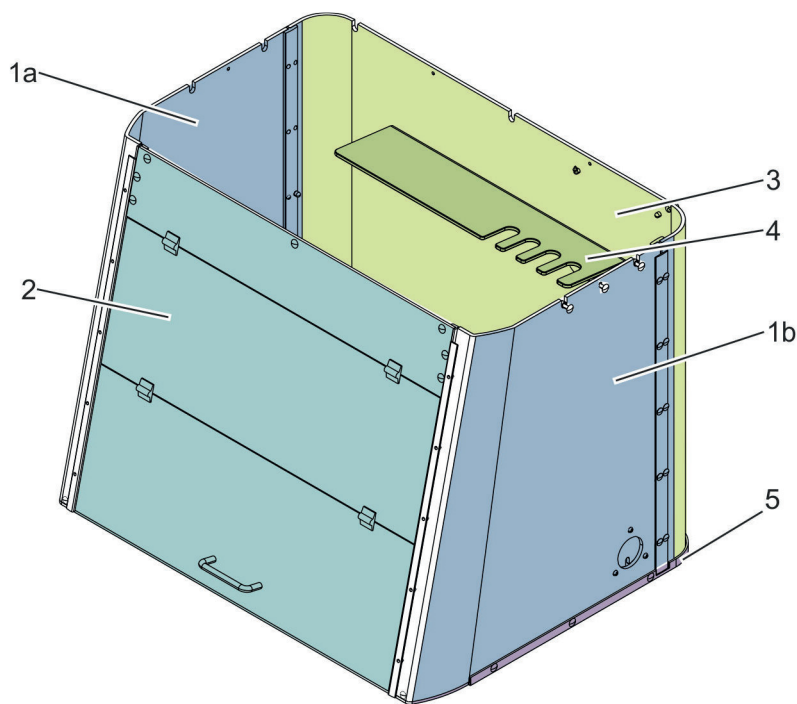
Fig. 2: Installed PC glass chamber of the model Laminar Flow, size 32"

1a + b	Side panels	3	Back panel	5	Base plate
2	Front door	4	Multifunctional shelf		

Plate panels

Just like the front and the back panel, the side panels are made of 6-mm clear polycarbonate (PC) glass, with overlap between the segments to eliminate potential exposure to harmful shortwave UV irradiation during UV 254 decontamination. The right side panel is additionally equipped with a cable duct, through which cables can be routed from the outside to the interior of the PC glass chamber. UV irradiation cannot escape here because it has a separate PC glass cover.

Back plate

The back plate is made from 6-mm clear polycarbonate glass, with overlap between the segments to eliminate potential exposure to harmful shortwave UV irradiation during UV 254 decontamination.

Multifunctional shelf

The multi-functional shelf is mounted on the PC back panel of the glass chamber indoors. It is used to set up samples and to store pipettes or other additional tools. You can also store samples there that need to be decontaminated because the shelf is within the irradiation area of the UV lamp.

Base panel

The base panel is made of a polypropylene blend that reflects UV light, just like the front, side and back panels do.

Front door

The front panel is made from 6-mm clear polycarbonate glass with overlap between the segments to eliminate potential exposure to harmful shortwave UV irradiation during UV 254 decontamination. The front door can be opened in two steps.

- › The opening of the lowest segment facilitates the user’s work process at the GuardOne® Workstation.
  - › Further opening the middle segment makes it easier for the user to clean and maintain the PC glass chamber from the inside.
- It is not recommended to completely open both segments for the work process, as a clean air environment cannot be guaranteed otherwise.

The front door is additionally equipped with a magnetic switch, which registers when the front door is open. In the case of an open front door, the UV light source is automatically switched off. The same applies to the airflow: When closing the front door, the blower automatically stops.

## Top housing unit

The top housing is a preinstalled unit in which the blower motor, pre- and HEPA filter are already installed, provided that it is the Laminar Flow model of the GuardOne® Workstation.

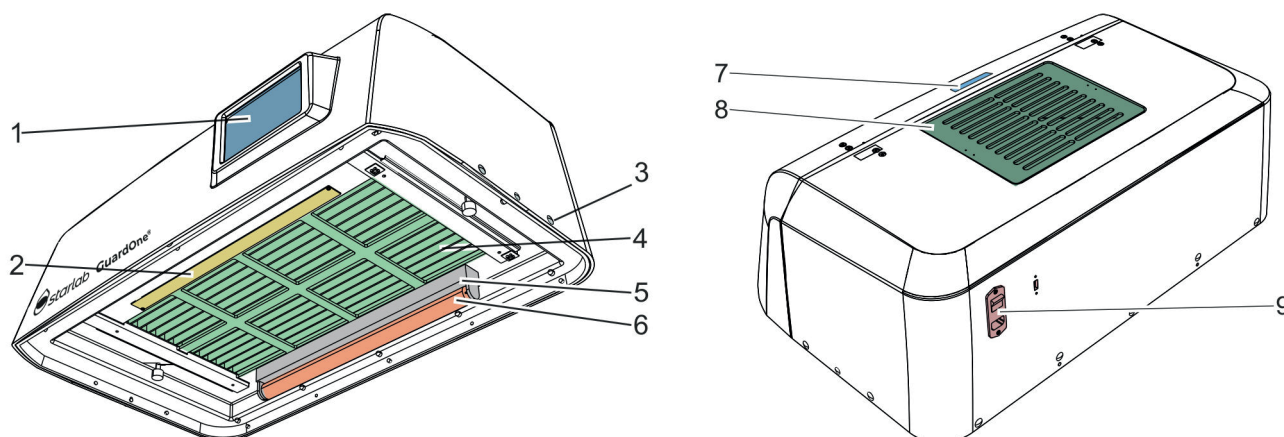
Fig. 3: Pre-installed top housing unit of the model Laminar Flow, size 32"

- |   |   |
|---|---|
| 1 Touchscreen display with graphical user interface | 6 UV lamp   |
| 2 LED light panel                                   | 7 Recess  |
| 3 Plastic rivets                                    | 8 Coarse- and pre-filter                                      |
| 4 Filter grid with HEPA filter inside               | 9 Main power switch, main power connection and USB connection |
| 5 UV bulb holder                                    |   |

### The top housing unit contains following components:

- › Coarse- and pre-filter consisting of a filter cover and filter fleece on the roof
  - On the roof of the top housing unit there is access to the pre-filter.
  - Protected by a removable coarse filter, the pre-filter can easily be reached.
- › Integrated HEPA filter
  - The HEPA filter is located in the interior of the top housing unit.
  - The life time of the HEPA filter is 600 days.
- › Safety contact in form of a magnetic switch in the inside
  - Like the front door, the top housing unit can be opened at its front side.
  - It is equipped with a magnetic switch that automatically switches off the airflow and the touchscreen display when the top housing flap is open.
- › UV light bulb holder in the back underneath the HEPA filter
  - The GuardOne® Workstation uses UV light with a fixed wavelength of 254 nm for decontamination and sterilization of the working area.
  - Depending on the model, one or two UV lamps marked with a serial number are included. The UV lamps must be installed in the bulb holders after the assembly of the glass chamber and the top housing unit.
  - The bulb holder for the UV lamp protrudes a few centimetres from the top housing unit.
  - The life time of the UV lamps is 9000 hours.
- › LED light panel in the front underneath the HEPA filter
  - The LED light panel is pre-installed.
  - Depending on the photosensitivity of the samples, you can switch the LED light on and off and change the settings for LED brightness or the LED colour between normal white and red.
- › Touchscreen display with graphical user interface on the front
  - The operating unit is located on the front side. The GuardOne® Workstation is operated via a touchscreen display with the graphical user interface.
  - The display with the graphical user interface can be used to set the settings for the airflow, the LED and UV light. The display also provides information about the manufacturer and health of the GuardOne® Workstation.
- › Main power switch, main power connection and USB connection at the back
  - The main power switch and connection as well as a USB connection are located on the back. The latter can be used for installing software updates.
- › Recess in the roof to keep the front door open to clean or maintain the PC glass chamber
  - When the front door is folded up, the handle can be placed in the recess to prevent slipping.
- › Plastic rivets, thumbscrews and two cables in a separate bag
  - For assembly, plastic rivets, thumbscrews and two cables are included.

**i** Spare parts must only be ordered from the manufacturer Starlab.  
For further information, contact your local Starlab sales representative.

## 2 Safety

### 2.1 Explanation of Safety Symbols

#### Safety instructions

Safety instructions are indicated by symbols in this manual. The safety instructions are initiated by signal words that indicate the degree of the danger.

#### **DANGER!**

This combination of symbol and signal word indicates an imminently dangerous situation which, if not avoided, will result in death or serious injury.

#### **WARNING!**

This combination of symbol and signal word indicates a potentially dangerous situation which, if not avoided, could result in death or serious injury.

#### **CAUTION!**

This combination of symbol and signal word indicates a potentially dangerous situation which, if not avoided, could result in slight or minor injuries.

#### **NOTICE!**

This combination of symbol and signal word indicates a potentially dangerous situation which, if not avoided, could result in damage to property.

#### **ENVIRONMENT!**

This combination of symbol and signal word indicates potential dangers to the environment.

This symbol highlights useful tips and recommendations as well as information for an efficient and trouble-free operation.

#### Safety instructions in procedures

Safety instructions can refer to specific, individual instructions. Embedded in the procedure, such safety instructions do not interrupt the flow of reading when performing the action. The signal words described above will be used.

#### For example:

1. Loosen the screw.
2.  **CAUTION!**  
Risk of entrapment on the panel!  
Close the panel carefully.
3. Tighten the screw.

#### Safety labelling at the GuardOne® Workstation

#### **WARNING!**

**Danger due to illegible labeling!** Over time, stickers and signs can soil so that dangers cannot be detected and necessary operating instructions cannot be followed. As a result, there is a risk of injury.

- Always keep safety, warning and operating instructions in a legible condition.
- Renew damaged stickers and signs immediately.

#### **UV-C irradiation**

**Danger due to UV-C irradiation!** UV-C irradiation from the biocidal UV lamp is harmful to eyes and skin.

- › Avoid eye and skin exposure to unshielded radiation.
- › Only use in an enclosed environment which shields users from the irradiation like the GuardOne® Workstation interior.
- › Turn off the UV irradiation when you finish the decontamination of the GuardOne® Workstation.

#### **Electrical hazard**

**Danger due to electrical hazard!** Marked component can be energized.

Check the operating manual (Chapter 4.6) or contact your local Starlab service technician for assistance.

#### **Refer to manual for safety information**

This symbol indicates hazardous areas. Read the operating manual before performing the action.

## 2.2 Intended Use

The GuardOne® Workstation is to be used only for the intended use described herein. The GuardOne® Workstation is a containment equipment using air filtration technology and UV irradiation to create a working environment which minimizes the risk of cross-contamination under standard laboratory working conditions. The GuardOne® Workstation exclusively serves for process protection. The compliance with all instructions in this operating manual is part of the intended use. Any use beyond the intended is considered as misuse. All country-specific safety requirements for operating electrical equipment in the laboratory must be observed.

### **WARNING!**

#### **Risk of misuse!**

Misuse of the GuardOne® Workstation can lead to dangerous situations for persons and property.

- Only install Starlab spare parts and wear parts in the GuardOne® Workstation. The installation of spare parts and wear parts not authorized by Starlab results in malfunction of the GuardOne® Workstation, contamination of the work environment, and damage to the sample material.
- Use the GuardOne® Workstation exclusively within the specified conditions of use which are supplied in the specifications (Chapter 10 "Specifications").

### **WARNING!**

#### **Danger due to property protection only!**

The GuardOne® Workstation exclusively serves for process protection. There is no personal protection included.

- Do not use pathogens within the GuardOne® Workstation.

## 2.3 Safety Installations

### **Main power switch**

The main power switch is also designed as an emergency stop switch. It is located at the rear of the top housing unit in the upper right corner. Turning the main switch to position „0“, results in the airflow and UV decontamination stopping immediately.

The GuardOne® Workstation is switched off. Electrical elements, such as the display, are de-energized.

### **Safety contact on the front door**

To ensure that access to the inside is only possible with the UV lamp switched off, the lower segment of the front door is equipped with safety contacts. At the same time, the safety contacts prevent the airflow from blowing when the front door is closed. The UV decontamination of the GuardOne® Workstation can only be started when the front door is closed. At the same time, closing the door means that the airflow stops working.

## 2.4 Responsibilities of the Operator

### **Operator**

The operator is the person who operates the GuardOne® Workstation for commercial, economic or research purposes.

### **Operator obligations**

In addition to the safety instructions in this operating manual, the operator has to observe the safety, occupational safety and environmental regulations. The following applies in particular:

- › The operator must be aware of the applicable health and safety regulations. He has to identify hazards in a risk assessment that are caused by the specific working conditions in the work area of the GuardOne® Workstation.
- › During the entire operating time of the GuardOne® Workstation, the operator has to check if these operating instructions comply with the current state of the regulations. If necessary, the operator has to adjust them.
- › The operator has to regulate and set the responsibilities for installation, operation, troubleshooting, maintenance and cleaning.
- › The operator has to ensure that all persons who handle the GuardOne® Workstation have read and understood this operating manual. Additionally, the operator has to keep the staff trained and informed about dangers on a regular basis.

Furthermore, the operator is responsible for ensuring that the GuardOne® Workstation is always in perfect technical condition.

Therefore:

- › The operator must ensure that the maintenance intervals are observed.

2.5 Personnel Requirements

Personnel

Only persons who can be expected to perform their work reliably are allowed to perform the installation, operation and maintenance of the GuardOne® Workstation. Persons whose reactivity is impaired, for example by drugs, alcohol or medications, are not allowed to install, operate or maintain the GuardOne® Workstation. Observe the age and job-specific regulations that apply at the installation site when selecting the personnel.

There is no distinction between installation, operation and maintenance personnel. Suitable staff is allowed to complete all tasks after a briefing by the operator. Only use staff with a training in laboratory tasks for performing tasks with the GuardOne® Workstation. Usually, this is a laboratory specialist. As such, the personnel is educated in his field of activity and knows the relevant standards and provisions. Due to their professional training and work experience, the personnel can execute work along the biological and chemical process chain in the field of process, production and biotechnology. The personnel can recognize possible dangers independently and avoid them. The personnel considers the regulations in health and environmental protection, occupational safety and quality assurance.

The personnel is responsible for the following tasks:

- > Setting up the GuardOne® Workstation for the first time
- > Assembling the GuardOne® Workstation (Chapter 4 “Assembly”)
- > Establishing and checking the operational readiness of the GuardOne® Workstation
- > Performing a function test
- > Commissioning
- > Changing the factory settings
- > Operating the graphical user interface
- > Setting the GuardOne® Workstation for work
- > Adjusting the airflow, operate LED light and UV decontamination
- > Maintaining the GuardOne® Workstation (replace UV bulb, pre- and HEPA filter)
- > Cleaning and decontaminating the GuardOne® Workstation
- > Performing troubleshooting
- > Disassembling and disposing the GuardOne® Workstation

2.6 Personal Protective Equipment

Personal protective equipment protects personnel from impairment of safety and health at work. The work at the GuardOne® Workstation does not require any personal protective equipment. The responsibility for wearing the necessary personal protective equipment lies with the operator.

CAUTION!  
Health hazard due to inadequate protection!

- Depending on the type of sample being processed in the workstation, a personal protective equipment may be required.
- The responsibility for the sample processing and the provision of the necessary personal protective equipment lies with the operator. As such, he must provide the necessary protective equipment.
  - Always wear suitable protective equipment when working with potentially harmful specimens.

2.7 Hazards when Used as Intended

The following section identifies residual risks that exist during intended use. To reduce the risks of personal injury and property damage and to avoid dangerous situations, observe the safety instructions in the following sections of this operating manual.

## Dangers to persons

### **DANGER!**

#### **Danger to life due to electric current!**

Contact with live parts causes immediate danger to life by electric shock. A damage to the insulation or of individual components can be life-threatening.

- In case of damage to the isolation, turn off the power supply immediately and arrange a repair.
- Before performing any cleaning work, switch the main power to OFF.

### **WARNING!**

#### **Risk of injury due to high-power UV irradiation!**

The UV lamp emits high-power UV irradiation. A long-term exposure to UV light results in damage to health.

- Avoid eye and skin exposure to unshielded radiation.
- Only use in an enclosed environment which shields users from the irradiation like the GuardOne® Workstation interior.
- Turn off the UV irradiation when you finish the decontamination of the GuardOne® Workstation.
- In case of a malfunction of the UV lamp, contact the Starlab service technician immediately.

### **WARNING!**

#### **Risk of injury due to rubbing alcohol (ethanol)!**

Ethanol is an eye irritant. The alcohol vapor can cause light-headedness and drowsiness.

- Before working with ethanol, switch the main power to OFF.
- Ensure proper ventilation when working with ethanol. If you feel tired or light-headed, immediately breathe fresh air.
- Before a longer stay, ventilate the work area.
- In case of persistent drowsiness, seek medical assistance.
- Always keep ethanol storage containers tightly closed.
- Do not eat, drink or smoke in the working area.
- In case of accidental contact with ethanol, immediately initiate first aid measures:
  - Take off contaminated clothing immediately.
- **Inhalation:** Provide for fresh air. In case of complaints, consult a doctor.
- **Skin contact:** Wash off the ethanol with soap and plenty of water. In case of complaints, consult a doctor.
- **Eye contact:** Rinse the eyes immediately with water. If possible, remove inserted contact lenses and continue to rinse with water. Consult a doctor.
- **Ingestion:** Consult a doctor immediately.

### **WARNING!**

#### **Risk of fire due to objects with high heat development!**

Working with open flames or objects with high heat development in the GuardOne® Workstation can cause serious injuries and property damage.

- Do not smoke in the working area of the GuardOne® Workstation. Avoid handling open light, fire and ignition sources of all kinds.
- Take the following measures in case of fire outbreak:
  - Switch off the main power of the GuardOne® Workstation. Unplug the cable connector.
  - If there is no danger to your own health, save persons from the danger zone. If necessary, initiate first aid measures.
  - Alert fire and/or rescue services.
  - If there is no danger to your own health, contain or extinguish the fire using adequate equipment. Continue the fire-fighting until the arrival of the fire services.
  - Inform the person in charge at the location.
  - Keep access routes free for rescue vehicles. Instruct rescue vehicles.

### **CAUTION!**

#### **Risk of injury due to tripping over objects, cables and wires cluttering the space!**

Objects in the workspace of the GuardOne® Workstation are tripping hazards.

- Keep the working area clean.
- Store tools, spare parts, wear parts and detergents away from the GuardOne® Workstation.
- Remove items from the workspace when they are no longer needed.
- Bundle cables and wires.
- Identify unavoidable tripping hazards (for example, cable lines) and mark them.

Property damage

**NOTICE!**  
**Property damage due to operation of the touchscreen with sharp objects!**

The display on the top housing has a touch screen. Due to the pressure sensitivity, there is a risk of scratches or breakage.

- Never operate the touch screen with sharp objects, for examples biros.
- Always operate the touchscreen with your fingers. Apply only light pressure.

**NOTICE!**  
**Property damage to plastic materials during prolonged irradiation!**

Some types of plastic, such as pipette tips, can become brittle when exposed to prolonged irradiation.

- Do not place plastic containers that are not UV stable in the GuardOne® Workstation during prolonged exposure to UV light.

**NOTICE!**  
**Broken screws due to misposition!**

If inserted out of position, screws can break out. This can cause damage to the top housing and the glass chamber of the GuardOne® Workstation. A decontaminated work environment can no longer be guaranteed.

- Carefully insert the screws into the aligned holes during installation.
- Follow the instructions in this operating manual. Install the PC glass chamber and the top housing of the GuardOne® Workstation carefully.

Dangers to the environment

**ENVIRONMENT!**  
**Dangers to the environment due to incorrect handling of environmentally hazardous substances!**

Incorrect handling and disposal of environmentally hazardous substances can cause significant damage to the environment.

- Always observe the instructions for handling environmentally hazardous substances and their disposal below.
- If environmentally hazardous substances escape into the environment accidentally, take appropriate measures immediately. If in doubt, inform the responsible authority about the damage and ask for appropriate action to take.

The following environmentally hazardous substances are used:

**Cleaning liquids**

Solvent-based cleaning agents contain toxic substances. They must not escape into the environment. Disposal must be carried out by a disposal specialist and is the responsibility of the operator of the GuardOne® Workstation.

**UV lamps**

Defective UV lamps are hazardous waste. They contain quicksilver and cannot be recycled. Contact your in-house or local waste management system representative for assistance or further information.

## 2.8 Behaviour in Case of Emergency

### Behaviour in case of accidents

#### Preventive measures

- > Always be prepared for accidents!
- > Keep first aid facilities (first aid kit, blankets etc.) in working order and ready at hand.
- > Make sure that the staff is familiar with accident reporting, first aid and rescue facilities.
- > Keep access routes free for rescue vehicles.

#### Measures in case of accidents

- > Switch the main power of the GuardOne® Workstation off (Chapter 4.6 “Complete Installation”).
- > If there is no danger to your own health, save persons from the danger zone.
- > If necessary, initiate first aid measures.
- > Alert fire and/or rescue services.
- > Inform the person in charge at the location.
- > Keep access routes free for rescue vehicles.
- > Instruct rescue vehicles.

### Behaviour in case of fire outbreak

#### Preventive measures

- > Always be prepared for a fire outbreak!
- > Keep first aid facilities (first aid kit, blankets etc.) and firefighting equipment in working order and ready at hand.
- > Make sure that the staff is familiar with accident reporting, first aid and rescue facilities.
- > Keep access routes free for rescue vehicles.

#### Measures in case of fire outbreak

- > Switch off the main power of the GuardOne® Workstation (Chapter 4.6 “Complete Installation”) and unplug the cable connector.
- > If there is no danger to your own health, save persons from the danger zone.
- > If necessary, initiate first aid measures.
- > Alert fire and/or rescue services.
- > If there is no danger to your own health, contain or extinguish the fire using adequate equipment. Continue the fire-fighting until the arrival of the fire services.
- > Inform the person in charge at the location.
- > Keep access routes free for rescue vehicles.
- > Instruct rescue vehicles.

### Behaviour in case of a power failure

#### Measures after a power failure

1. Turn the main power switch to ON.
2. Continue with your application.

**i** You do not have to reapply previous settings. Previous settings are retained in the event of a power failure.

## 3 Transport, Packing and Storage

### 3.1 Safety during Transport, Packing and Storage

#### Safety indications

##### **WARNING!**

##### **Risk of injury due to heavy loads!**

During lifting operations and when setting up the package, loads can fall off. This can cause serious injury.

- Consider dimensions and weight of the GuardOne® Workstation.
- Components of the GuardOne® Workstation are heavy. Lift only with the help of a second person or suitable equipment.

##### **NOTICE!**

##### **Damage to the UV lamp due to bumping of the package!**

As a result of shocks to the package, the UV lamp can be damaged.

- Transport the package carefully.
- In the event of damage claims, contact your local Starlab sales representative immediately.

##### **NOTICE!**

##### **Property damage due to placing on a hard base!**

The bulb holder for the UV lamp protrudes a few centimeters from the top housing unit. If you place the unit flat on a hard base, the bulb holders may be pressed in and can be damaged.

- Do not store the top housing unit unpacked on a hard base.

**i** In case of incompleteness or damage, contact your local Starlab sales representative immediately. Contact details can be found “Starlab Service Hotline” or the back page.

### 3.2 Inspection of Transport and Delivery

#### **Delivery directly through the manufacturer**

The first delivery of the GuardOne® Workstation takes place via parcel service directly from the manufacturer. Protected by foam insulation and bubble wrap, all included components are delivered in one package.

Below, you find further information about the scope of delivery as well as the correct procedure for dealing with an incomplete or damaged delivery.

#### **The scope of delivery comprises the following components:**

- > 1x top housing unit (with 1 or 2 pre-installed pre- and HEPA filter).
- > 1x base plate
- > 1x front door (pre-assembled)
- > 1x back panel
- > 2x side panels
- > 1x multifunctional shelf
- > 1 or 2x UV lamps (with serial number on package)
- > 1x bag with two power cords (IEC 60320 C13/C14-type with earth contact)
- > 1x bag with plastic rivets

#### **Transport inspection**

Check the delivery upon receipt for completeness and transport damage. In case of externally visible transport damage, proceed as follows:

- > Do not accept the delivery or accept only with reservation.
- > Note the extent of damage on the transport documents.
- > Initiate the complaint.

**i** Claim for any damages as soon as they are recognized. Claims for damages can only be made within 15 working days after purchase of the GuardOne® Workstation.

**i** In case of incompleteness or damage contact your local Starlab sales representative immediately. Contact details can be found “Starlab Service Hotline” or the back page.

## In-house transport

### WARNING!

#### Risk of injury due to heavy loads!

During lifting operations and when setting up the package, loads can falling off. This can cause serious injury.

- Consider dimensions and weight of the GuardOne® Workstation.
- Components of the GuardOne® Workstation are heavy. Lift only with the help of a second person or suitable equipment.

You can move the GuardOne® Workstation to another location with the help of a second person without disassembling it.

## 3.3 Handling of Packaging Material

### Packaging material

Depending on parcel service and/or inhouse transport directly to the final recipient, the manufacturer delivers the GuardOne® Workstation in a secured state. To protect against damage and contamination during transport, the GuardOne® Workstation is packed in a single cardboard package with individual pieces additionally wrapped in protective foil.

### Disposal

For disposal of packaging materials, pay attention to the applicable legal and local regulations.

### ENVIRONMENT!

#### Danger to the environment due to incorrect disposal!

Packaging materials are valuable raw materials. Sensibly processed and recycled, they can be used again in many cases. Dangers for the environment can arise from incorrect disposal of packaging materials.

- Dispose of packaging materials in an environmentally friendly manner.
- Observe the locally applicable disposal regulations. If necessary, commission a specialist company with the disposal.
- Dispose of the pallet in an environmentally friendly way or store dry until reuse.

## 3.4 Storage

### Requirements for the storage location

Choose the storage location considering the following aspects:

- > Do not store the GuardOne® Workstation outdoors.
- > Store dry and dust-free.
- > Do not expose to aggressive media.
- > Protect from sunlight.
- > Permissible storage temperature: -5 – +55 °C
- > Permissible humidity: 10 – 95 %

### Requirements for the storage

Store the GuardOne® Workstation under the following conditions:

- > The GuardOne® Workstation is out of service.
- > Do not put any objects on the GuardOne® Workstation.
- > When storing for more than 3 months, regularly check the general condition of all parts.
- > Before putting into service again, check the pre-filter visually and replace it if heavily soiled.

## 4 Assembly

### NOTICE!

#### Property damage due to incorrect installation!

- The installation must only be carried out by experienced staff.

### NOTICE!

#### Property damage at the pre-filter due to obstruction!

An obstructed pre-filter will have a negative impact on air filtration efficiency and can cause blocked filter alarms.

- Make sure that the pre-filter is not obstructed in any way at any time.

### Choosing an installation site

Before starting the assembly, make sure that the requirements at the installation site specified below are met.

- > The GuardOne® Workstation must be placed on a firm and level work surface.
- > A safe standing of the GuardOne® Workstation must be ensured.
- > The surface of the installation site has, at minimum, the same dimensions as the GuardOne® Workstation itself.
- > A minimum distance of 10 cm to adjacent objects and walls must be kept to both sides and the back of the GuardOne® Workstation.
- > The GuardOne® Workstation must be easily accessible from all sides. Especially at the back of the GuardOne® Workstation, there must be enough space to operate the main power switch on the right rear side of the top housing.
- > Above the GuardOne® Workstation, there must be a minimum of at least 25 cm free space to ensure free and unimpaired airflow.
- > The laboratory or workroom must be of sufficient size to provide ample space around the GuardOne® Workstation.
- > Adequate lighting must be available.
- > Adequate ventilation must be available.
- > A connection option to the power supply must be available.
- > Escape routes and rescue equipment are freely accessible.
- > All local safety requirements are fulfilled.
- > The installation site is dry and frost-free.
  - A minimum relative humidity of 40% is guaranteed.
  - A minimum room temperature of + 15 °C and a maximum room temperature of + 40 °C is guaranteed.
- > The GuardOne® Workstation is not exposed to direct sunlight.
- > The GuardOne® Workstation must be set up in a place that can easily be kept clean and dust free.

### Preparing the installation

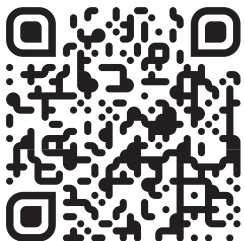
- > The GuardOne® Workstation is delivered only partly assembled to reduce transport volume by almost 50%.
- > The final assembly can be performed directly at the chosen installation site.
- > The assembly will take no longer than 15 minutes and does not require any special knowledge or any kind of tools.
- > Components of the GuardOne® Workstation are heavy. It is recommended to have a second person available for assistance.

In case of doubt, contact your local Starlab service or sales contact for assistance.

### Watch our video to GuardOne®:

<https://www.starlab.click/guardone-assembling>

## 4.1 Installing the PC Back Panel

1. Place the base plate (no. 5) on the selected installation site.
2. Take the back panel (no. 3) and insert it into the groove of the base plate. Make sure that the holes in the PC back panel are aligned with the holes in the base plate.
3. Use five of the plastic rivets (EVA02) to fix the PC back panel to the base plate. Push the plastic rivets into the aligned holes. For easiest assembly, orient the rivets with the open area facing upward. The PC back panel is installed.

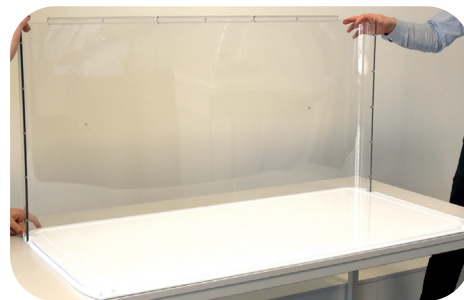
Fig. 4: Insert the back panel into the groove of the base plate

Fig. 5: Fix the back panel with plastic rivets

## 4.2 Installing the PC Side Panels and the Multifunctional Shelf

### Installing the PC left side panel

1. Take the PC left side panel (without cable duct, no. 1a) and insert it into the groove of the base plate. Make sure that the bottom holes in the PC left side panel are aligned with the holes in the base plate.
2. Make sure that the PC back panel slides fully into the groove of the aluminium profile on the PC left side panel. Make sure that the holes in the aluminium profile and the PC back panel are aligned.
3. Use five of the plastic rivets (EVA01) to fix the PC back panel to the aluminium profile of the left side panel. Push the rivets into the aligned holes.
4. Use three of the plastic rivets (EVA02) to fix the PC left side panel to the base plate. Push the rivets into the aligned holes. The left side panel is installed.

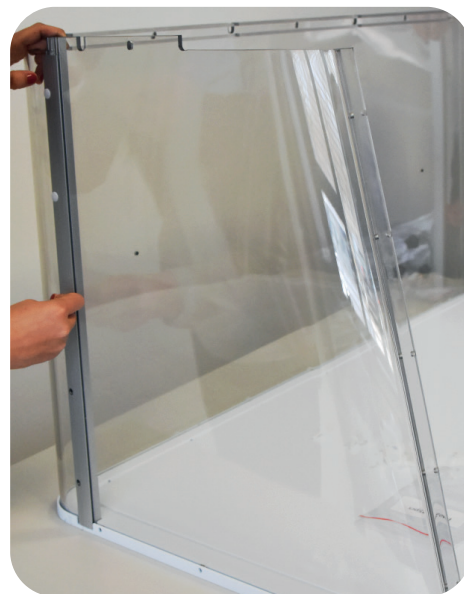
Fig. 6: Fix the PC back panel with plastic rivets

Installing the PC right sight panel (with cable duct)

Fig. 7: Fix the right side panel to the base plate

1. Take the PC right side panel (with cable duct, no. 1b) and insert it into the groove of the base plate. Make sure that the bottom holes in the PC right side panel are aligned with the holes in the base plate.
2. Make sure that the PC back panel slides fully into the groove of the aluminium profile on the PC right side panel. Make sure that the holes in the aluminium profile and the PC back panel are aligned.
3. Use five of the plastic rivets (EVA01) to fix the PC back panel to the aluminium profile of the right side panel. Push the rivets into the aligned holes.
4. Use three of the plastic rivets (EVA02) to fix the PC right side panel to the base plate. Push the rivets into the aligned holes. The right side panel (with cable duct) is installed.

Installing the PC shelf to the back panel

Fig. 8: Fix the PC shelf to the back panel

1. Fix the PC shelf (no. 4) to the PC back panel using the two black screw assemblies provided.
2. Position the PC shelf at the back panel on the inside of the PC glass chamber of the GuardOne® Workstation.

3. 

!

**NOTICE!**  
Property damage due to too high rotational torques!

Turn the two black screws into the designated holes on the PC shelf and the back panel by hand.

i

If necessary, you can provide additional support to the back panel with one hand from the outside.

The PC shelf is installed at the back panel on the inside of the glass chamber of the GuardOne® Workstation.

### 4.3 Installing the PC Front Door

**Precondition:**

> For safety reasons, have a second person available to help.

1. Take the pre-assembled front door element (no. 2) and insert it from above between the PC left and right side panels. Make sure that the entire panels slide into the groove of the aluminium profiles. Make sure that the holes in the aluminium profile and the PC side panels are aligned on both sides.

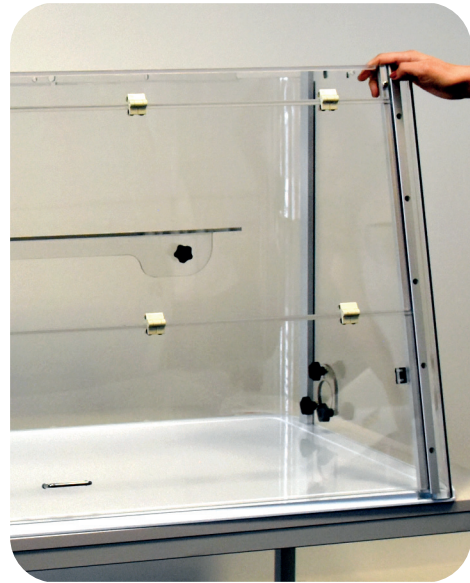
Fig. 9: Insert the front door between the side panels

2. Use five of the plastic rivets (EVA01) to fix the PC left side panel to the aluminium profile of the front door assembly. Push the rivets into the aligned holes.  
To avoid any deformation, you can provide additional support to the aluminium profile with one hand from the inside.
3. Use another five of the plastic rivets (EVA01) to fix the PC right side panel to the aluminium profile of the front door assembly. Proceed exactly as in the assembly of the PC left side panel to the front door detailed in step 2. The front door is installed between the side panels.

Fig. 10: Fix the front door with plastic rivets

## 4.4 Installing the Top Housing

Fig. 11: Assembled PC glass chamber

Fig. 12: Lift the top housing onto the PC glass chamber

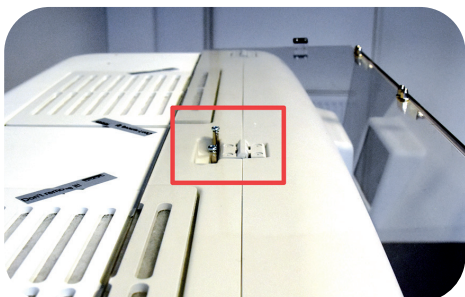
Fig. 13: Open position of the front door

Fig. 14: Install the top housing

### Preconditions:

- › The back and side panels are installed at the base plate.
- › The front door is installed between the side panels.
- › The top housing is the heaviest part of the GuardOne® Workstation. For safety reasons, have a second person available for assistance.

### NOTICE!

#### Property damage due to placement on a hard base!

The bulb holder for the UV lamp protrudes a few centimeters from the top housing unit. If you place the unit flat on a hard base, the bulb holder may be pressed in and can be damaged.

- Do not store the top housing unit unpacked on a hard base.
- After unpacking the top housing unit, lift it immediately with the help of a second person onto the preinstalled PC glass chamber.

### Lifting and fixing the top housing onto the PC glass chamber

#### 1. CAUTION!

#### Crushing hazard due to heavy weight!

Due to the heavy weight of the top housing, there is a danger of crushing your fingers.

- › Get assistance from a second person to place the top housing on the assembled PC glass chamber.
  - › Watch out for your fingers.
  - › Work carefully.
2. Take the pre-assembled top housing and lift it onto the assembled PC glass chamber.
  3. Make sure that the top housing and the PC glass panels are properly aligned. If aligned correctly, the top housing rests level on top of the aluminium profiles.
  4. Make sure that the round corners of the PC back panel are equally aligned with the top housing.
  5. For an easy and safe access to the inside of the GuardOne® Workstation, open the front door completely and secure it in the open position using the recess on the roof of the top housing.
  6. Use five plastic rivets (EVA02) to fix the top housing on the assembled PC glass chamber:
    - › Insert three plastic rivets from the back.
    - › Insert one plastic rivet from the left and right side each.
- i** If necessary, you can provide additional support to the PC glass panels with one hand from the inside.
7. Once all screws have been inserted, lower the front door again.

8. Open the flap of the top housing by reaching under the front cover with both hands. Slide the flap upwards until it clicks into place.
9. Use three of the plastic rivets (EVA02) to fix the front side of the top housing to the front door. Push the rivets into the aligned holes. The top housing is installed. The flap of the top housing must remain open.

Fig. 15: Fix the front side of the top housing

### Connecting the magnetic safety switch

1. Take the connecting cable adaptor at the right side of the opened flap of the top housing between right wall and HEPA filter.
2. Insert it from the top into the corresponding socket in the right aluminium profile of the front door assembly.
3. Plug the connector into the port. Make sure that the connector is securely plugged in.

**i** In case the magnetic safety switch has not been plugged in correctly, UV decontamination will not work!

4. Close the housing front door. To do so, first unlock the guiding rails by lifting the front door on both sides slightly upwards.

The magnetic safety switch between top housing and PC glass chamber is connected.

Fig. 16: Connecting the magnetic safety switch

## 4.5 Installing the UV Lamp

### Preconditions:

- › Remove the packaging of the UV lamp. On the packaging there is a sticker with the serial number of the UV lamp. You will need this serial number during the function test in Chapter 4.6 "Complete Installation".
1. Open the front door fully, using the recess on the top housing to safely secure it in position.
  2. Insert the new UV lamp from below into the lamp sockets. Lock the UV lamp by rotating it by approx. 90° in anticlockwise direction. As soon as the bulb has been locked, you will hear a clicking sound.
  3. In the 48" version you install two UV lamps side-by-side repeating the action detailed in step 2.
  4. Close the front door carefully.

Fig. 17: Rotate the UV lamp

4.6 Complete Installation

Fig. 18: Plug the power cord in the appropriate socket

Fig. 19: Standby mode

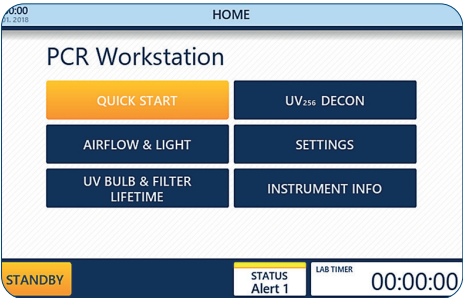
Fig. 20: Home menu

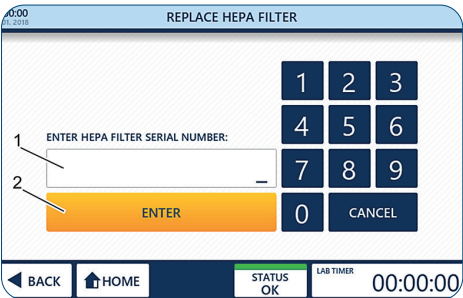
Fig. 21: Enter the serial number of the HEPA filter

Completing the assembly

1. Insert the power cord (IEC 60320 C13/C14-type with earth contact) into the socket located at the back of the top housing.
2. **NOTICE!**  
**Property damage due to increased voltage!**  
Make sure that the local strength of current conforms to the power rating of the GuardOne® Workstation.
3. Connect the power cord to the power supply.
4. To switch the GuardOne® Workstation ON (I) or OFF (O), use the main power switch located above the socket. Turn the main power switch to ON.

After switching the GuardOne® Workstation to ON, the touchpad with the graphical user interface at the front of the top housing will show the Standby screen.  
The GuardOne® Workstation is now ready for performing a function test.

Performing a function test

The function test shows if UV lamp, LED lights and HEPA filter are working.

Precondition:

- › The GuardOne® Workstation is fully assembled.
- › The main power switch is turned to ON.
- › After switching the GuardOne® Workstation to ON, the graphical user interface displays the Standby screen.

Press Tap to unlock to unlock the Standby screen.  
The Standby screen is unlocked.

You are now on the Home menu.

 When putting into service, the Home screen displays status alerts 1 and 3 when the GuardOne® Workstation is first switched on. This is to ensure that after initial installation, monitoring of the lifespan of the HEPA filter and UV lamp is activated.

Activating the HEPA filter

1. Press Settings.
2. Press Replace HEPA Filter.
3. Enter the 10-digit serial number (1) shown on the front of the filter labels.
4. Press Enter (2) to finish the installation procedure.
5. Press Home to return to the Home menu.

## Testing the Airflow

1. In the Home menu, press the Airflow & Light button.  
The page for adjusting blower speed or LED light opens.
2. To adjust the airflow, press onto the percentage value (1) in the Air Flow field.
3. Select the desired airflow rate using the + and - buttons  (2).

**i** Selecting an airflow rate of 0% will switch off the blower. To switch it on again, select an airflow rate above 0%. The available steps are 0, 50, 80 and 100%.

4. Apply the selection using Apply Settings (3).

**i** If the front door is closed, the airflow will not work. The airflow is working if the lowest segment of the front door or the whole front door is open. Even if the magnetic switch is installed improperly, LED light and blower function will be functional.

**i** When you turn on the airflow, the motor needs about ten seconds until the airflow starts to blow. When you turn off the airflow, it stops as soon as the front door is closed.

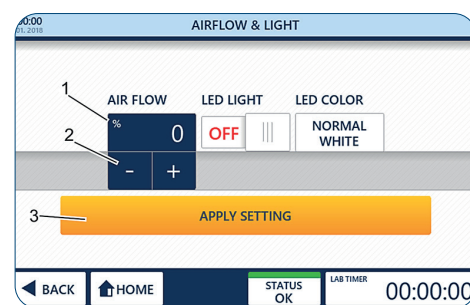
Fig. 22: Testing the Airflow

## Testing the LED light

- > You are in the Airflow & Light menu.
1. To switch on the LED light, use the LED light switch **ON** (1).  
Your selection is applied immediately.
  2. To select an LED light colour, press LED colour (2). Select the desired colour, normal white or red, by pressing the + and - button  (3).  
Your selection is applied immediately.

**i** If the LED lights are not working, please contact your local Starlab sales representative.

3. After you have made sure that the LED lights are working, switch off the LED light using the LED light switch **OFF** (1).
4. To return to the Home menu, either use Back or Home in the navigation bar.

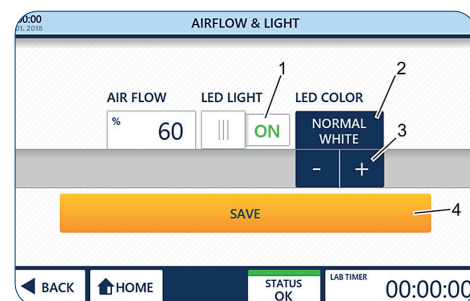
Fig. 23: LED light

## Activating the UV lamp

### Precondition:

- › The UV lamp is installed.
- › You are on the Home menu of the graphical user interface.
- › The packaging of the UV lamp including the sticker with the serial number, is available.

### Enter the serial number

1. Close the front door.

**i** As soon as the front door is closed, the airflow stops automatically.

2. Press Settings.
3. Press Replace UV Light.
4. Enter the 10-digit serial number (1) provided with the UV bulb.
5. Press Enter (2) to finish the installation procedure.
6. Press Home to get back to the Home menu.
7. Press UV254 Decon on the Home menu to open the parameter entry page.

### 8. CAUTION!

#### Health damage due to UV irradiation!

Long-term exposure to UV light may result in damage to health. For safety reasons, the UV lamp cannot be switched on while the front door is open. For safety reasons, the UV lamp will be switched off automatically as soon as the front door is being opened. Closing the door again will not restart the decontamination procedure.

- › To start the UV decontamination, switch the UV lamp On using the UV Status switch **ON** (1).
- › To stop the UV decontamination, switch the UV lamp Off using the UV Status switch **OFF** again.

**i** With correct setting and installation, the UV lamp emits a light blue light.

**i** If the UV lamp is not working, make sure that the magnetic switch connector is securely plugged into the port. If the UV light is still not working, contact your local Starlab sales representative.

**i** LED light switches off automatically after UV light is switched on.

9. After you have made sure that the UV lamp is working, switch the UV lamp off by using the UV Status switch **OFF**.
10. Press Home to return to the Home menu. After a successful function test, you can now change the pre-set parameter.

Fig. 24: Activate the automatic life time surveillance for UV light

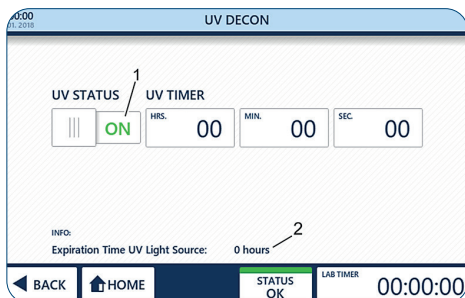
Fig. 24.1: UV 254 Decon

## 5 Putting into service

### Preconditions:

- > The GuardOne® Workstation is fully assembled.
- > The main power switch is turned to ON.
- > The serial numbers of HEPA filter and UV Lamp have been entered into the graphical user interface (Chapter 4.6 “Complete Installation”).
- > The function test was successful.

### Adjusting date and time

1. Press the date and time symbol (1) in the upper left corner of the information bar of any page to open the Date/Time page.
2. Change any value by pressing the value (2) and adjust it with the + and - button (3).
3. To save your changes, use Save Settings (4).
4. To exit the page, use Back or Home.

**i** Exiting the page without using Save Settings will discard all changes you have made!

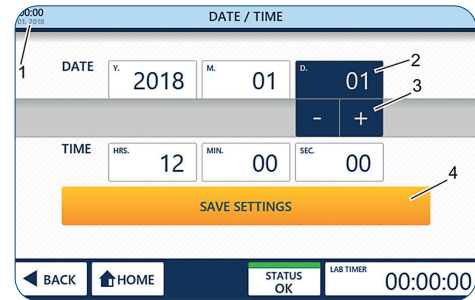
Fig. 25: Adjusting date and time

### Adjusting the default operation parameters

1. Press Settings on the Home menu.

**i** The Quick Start button on the Home menu will start the GuardOne® Workstation with default values for blower speed and LED light. The factory settings for the Quick Start are 50 % blower speed and LED lights off.

2. Press Quick Start Settings to open the setup page for changing the factory settings.
3. Change any value by pressing onto the corresponding buttons (1) and adjust the settings using the + and - button (2).
4. To save your changes, press Save (3).
5. To exit the page, use Back or Home.

**i** Exiting the page without using Save will discard all changes you have made!

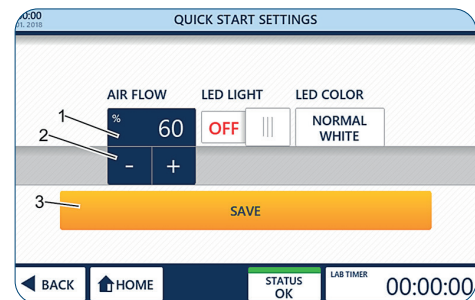
Fig. 26: Quick Start

## 6 Operation

### WARNING!

#### Incomplete decontamination due to permanent draught!

Due to the permanent draught in the glass chamber, a decontamination of 100 % by irradiation of the UV lamp is not possible. It is only possible to decontaminate the working environment by 99.995 %.

- Do not work with pathogenic laboratory samples.
- Follow the recommendations in this operating manual.
- For further information about the decontamination of the GuardOne® Workstation, refer to our homepage.

### CAUTION!

#### Health damage due to UV irradiation!

Long-term exposure to UV light results in damage to health.

- For safety reasons, the UV lamp cannot be switched on while the front door is open.
- For safety reasons, the UV lamp will be switched off automatically as soon as the front door is being opened. Closing the door again will not restart the decontamination procedure.

### NOTICE!

#### Property damage due to incorrect power supply!

An incorrect power supply will lead to property damage.

- Only connect the unit to power sources that match the electrical specifications on the nameplate of the unit.
- Only use sockets with proper electrical grounding.

## 6.1 Get to Know the Graphical User Interface

The GuardOne® Workstation is equipped with a graphical user interface which is operated via a touchscreen.

### Precondition:

- > The main power switch is turned to ON.

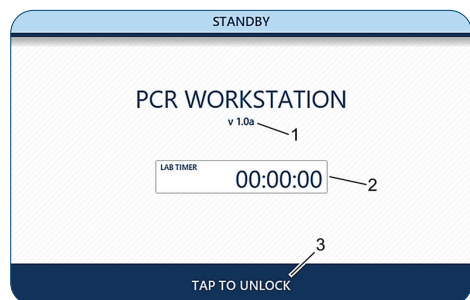
Fig. 27: Standby page

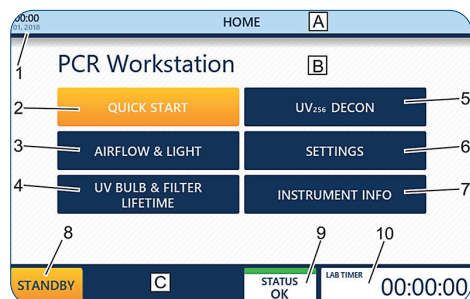
Fig. 28: Home page

### Get to know the Standby menu

#### The Standby page is shown:

- > immediately after switching on the GuardOne® Workstation.
- > after exiting the Home page using the Standby button in the navigation bar.

The Standby page informs you about the currently installed software version (1). In addition, if in use, the lab timer (2) will display the remaining time till the alarm will sound.

#### Exit the Standby page:

Press Tap to unlock (3) in the navigation bar. You are now on the Home page.

### Get to know the Home menu

The Home page offers access to the main functions of the GuardOne® Workstation.

#### Reaching the Home page from any other page:

Press Home in the navigation bar. You are now on the Home page.

#### Reaching the Home page from Standby page:

Press Tap to unlock in the navigation bar. You are now on the Home page.

<b>A</b> Information bar	<b>5</b> UV 254 Decon
<b>B</b> Display area	<b>6</b> Settings
<b>C</b> Navigation bar	<b>7</b> Instrument Info
<b>1</b> Date and time	<b>8</b> Standby
<b>2</b> Quick Start	<b>9</b> Status of the GuardOne® Workstation
<b>3</b> Airflow & Light	<b>10</b> Lab Timer
<b>4</b> UV bulb & Filter Lifetime	

## Adjusting the Lab timer

- i** The lab timer is working independent of any other function of the GuardOne® Workstation. You can switch the GuardOne® Workstation into standby mode without affecting the lab timer.

- 1 To use the lab timer, press the Lab timer symbol in the right corner of the navigation bar of any page. The Lab timer page opens.
- 2 To set the lab timer, press the buttons for hour, minute and second (1).
- 3 Enter the desired time values using the + and - button - + (2).
- 4 Press Start/Pause (3) to start or pause the timer.

- i** After the set time has elapsed, an alarm will be sounded. The alarm will stop automatically after approx. 5 seconds.

- 5 Press End + Reset (4) to stop and reset the timer to 00:00:00.
- 6 To exit the page, use Back or Home.

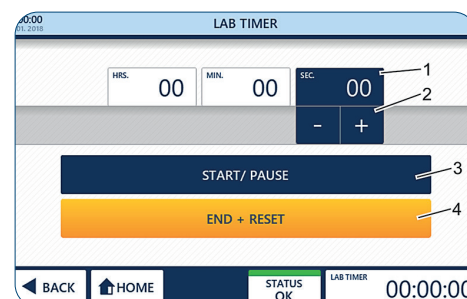
Fig. 29: Adjusting the lab timer

## 6.2 Operating the Display Area

### Quick Start and Quick Stop

When Quick Start is pressed, the GuardOne® Workstation will start working using pre-set parameters for blower speed and LED lights (factory settings). The factory settings are:

- > Blower speed is 50 %, Laminar Flow Model required.
- > LED light is turned off.

You can change the factory settings at any time during operation by pressing Settings followed by Quick Start Settings.

### Activating Quick Start

- 1 Open the front door.

- i** In case the front door is closed, the built-in magnetic switch will prevent the blower from working. The function of the LED light is independent of the front door.

- i** If the GuardOne® Workstation is in UV decontamination mode, the UV lamp will be switched off too.

- 2 Press Quick Start (1).  
The GuardOne® Workstation starts working.

### Activating Quick Stop

After Quick Start has been pressed, the button will change to Quick Stop.

Press Quick Stop (1) to switch off the blower and LED light. The GuardOne® Workstation stops working.

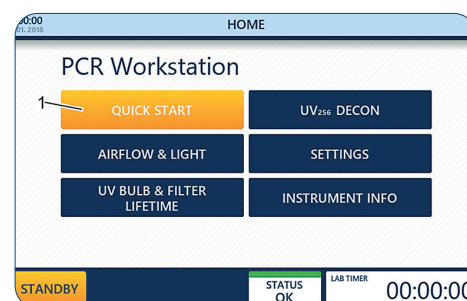
Fig. 30: Quick Start

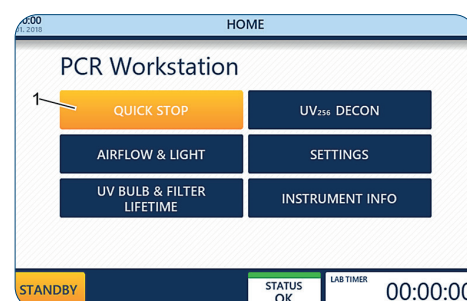
Fig. 31: Quick Stop

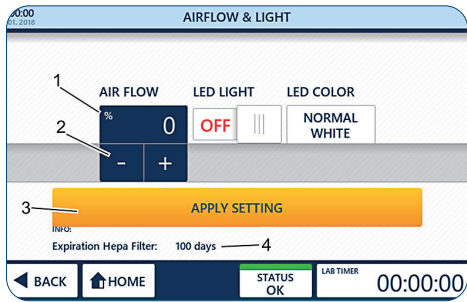
Fig. 32: Adjust the Airflow

## Airflow & Light

By pressing Airflow & Light, you can adjust the blower speed or the LED light at any time during operation.

### Precondition:

- > You are on the Home page.

1. Press Airflow & Light.

### Adjusting the Airflow

2. Press the percentage value (1).
3. Select the desired airflow rate (2) using the + and - buttons .

**i** You can choose between 0, 50, 80 and 100 % airflow rate.

**i** Selecting an airflow rate of 0 % will switch off the blower. To switch it on again, either select an airflow rate above 0 % or use the Quick Start function on the Home page ("Quick Start and Quick Stop").

**i** The Expiration HEPA Filter field (4) provides information about the remaining life time of the HEPA filter.

4. Apply the selection using Apply Settings (3).  
Your selection is applied immediately.

5. To exit the page, use Back or Home.

**i** Exiting the page without using Apply Settings will discard all changes you have made!

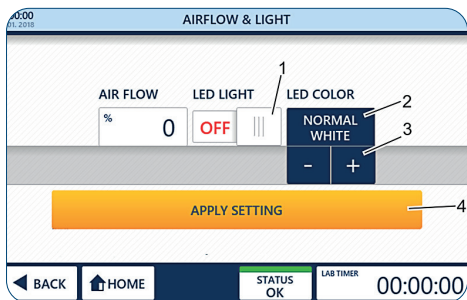
Fig. 33: Operate LED light and colour

### Operating the LED light

1. To switch on/off the LED light, press the LED Light switch (1).  
Your selection is applied immediately.
2. To select an LED light colour, press LED Colour (2).
3. Select the desired colour. You can choose between normal white (default) and red (3) by pressing the + and - button . Your selection is applied immediately.
4. To exit the page, use Back or Home.

UV Bulb & Filter Lifetime

By pressing UV bulb & Filter lifetime, you can view the information on the remaining lifetime of the UV bulb and the HEPA filter.

The life time of the UV bulb is 9000 hours.  
The life time of the HEPA filter is 600 days.

- 1. Press UV bulb & Filter lifetime.  
The remaining life times of the HEPA filter (1) and the UV bulb (2) are displayed.
- 2. To exit the page, use Back or Home.

Instrument Info

By pressing Instrument Info, you can view information about the current software version and the blower hours.

- 1. Press Instrument Info.  
The current software version (1) and the blower hours (2) are displayed.
- 2. To exit the page, use Back or Home.

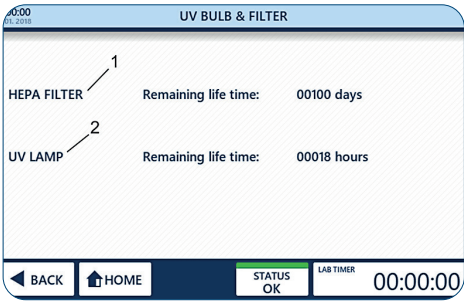
Fig. 34: Information about UV bulb and Filter lifetime

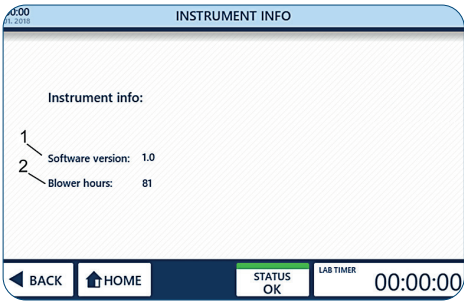
Fig. 35: Instrument Info

## Bacterial destruction through UV irradiation

UV light helps removing (or reducing the number of) airborne microorganisms from a closed environment. It is proven that UV light destroys almost all DNA-based microorganism given a sufficient UV dosage and irradiation time. But many variables like airflow, humidity, distance of microorganisms to UV light or UV dosage are present in a real-world environment that make actual calculation of the irradiation time very difficult. Before performing the steps to adjust the UV 254 Decon, refer to the calculation example for the minimum irradiation time, including the formula as well as the table of recommended times for decontaminating minimum from common organisms.

### The germicidal destruction rate is calculated as follows:

$\text{Microwatt seconds/cm}^2 = \text{microwatts/cm}^2 \times \text{seconds of exposure}$

The bacterial destruction chart below indicates the amount of shortwave (254 nm) UV energy required for complete destruction of various organisms.

**i** Please note that optimal working conditions are required for the recommended duration of irradiation. Changing only one factor results in a changed required irradiation time.

Bacteria	Required Microwatt ( $\mu\text{W}$ ) seconds/cm <sup>2</sup>	Recommended minimum exposure time in minutes in 1 m distance by a UV-C power of 48 $\mu\text{W/cm}^2$
Bacillus anthracis	8700	3
Bacillus megatherium sp. (veg.)	2500	1
Bacillus megatherium sp. (spores)	5200	2
Bacillus paratyphosus	6100	2
Bacillus subtilis (veg.)	11000	4
Bacillus subtilis (spores)	22000	8
Clostridium tetani	22000	8
Corynebacterium diphtheriae	6500	2
Eberthella typhosa	4100	1
Escherichia coli	6600	2
Micrococcus candidus	12300	4
Micrococcus sphaeroides	15400	5
Mycobacterium tuberculosis	1000	0
Neisseria catarrhalis	8500	3
Phytomonas tumefaciens	8500	3
Proteus vulgaris	6600	2
Pseudomonas aeruginosa	10500	4
Pseudomonas fluorescens	6600	2
Salmonella enteritidis	7600	3
Salmonella sp.	10000	3
Salmonella typhimurium	15200	5
Sarcina lutea	26400	9
Serratia marcescens	6160	2
Shigella paradysenteriae	3200	1
Spirillum rubrum	6160	2
Staphylococcus albus	5720	2
Staphylococcus aureus	6600	2
Streptococcus hemolyticus	5500	2
Streptococcus lactis	8800	3
Streptococcus viridans	3800	1

Yeasts	Required Microwatt ( $\mu W$ ) seconds/ $cm^2$	Recommended minimum exposure time in minutes in 1 m distance by a UV-C power of 48 $\mu W/cm^2$
Brewer's yeast	13200	5
Baker's yeast	13200	5
Saccharomyces cerevisiae	13200	5
Saccharomyces ellipsoideus	13200	5
Saccharomyces sp.	17600	6

Mould spores	Required Microwatt ( $\mu W$ ) seconds/ $cm^2$	Recommended minimum exposure time in minutes in 1 m distance by a UV-C power of 48 $\mu W/cm^2$
Aspergillus flavus	99000	34
Aspergillus glaucus	88000	31
Aspergillus niger	330000	115
Geotrichum candidum (Oospora lactis)	11000	4
Mucor racemosus	35200	12
Penicillium digitatum	88000	31
Penicillium expansum	22000	8
Penicillium roqueforti	26400	9
Rhizopus nigricans (R. stolonifer)	220000	76

Viruses	Required Microwatt ( $\mu W$ ) seconds/ $cm^2$	Recommended minimum exposure time in minutes in 1 m distance by a UV-C power of 48 $\mu W/cm^2$
Bacteriophage (E. coli)	6600	2
Influenza	6600	2
Tobacco mosaic	44000	15

Protoza	Required Microwatt ( $\mu W$ ) seconds/ $cm^2$	Recommended minimum exposure time in minutes in 1 m distance by a UV-C power of 48 $\mu W/cm^2$
Chlorella vulgaris (algae)	22000	1
Nematode eggs	9200	3
Paramecium	200000	69

UV 254 Decon

By pressing UV 254 Decon, you can activate the UV decontamination mode. To automatically stop the UV decontamination after a pre-set time, you can use the timer function.

!

**NOTICE!**

**Property damage to plastic materials from prolonged irradiation!**

Some types of plastic, such as pipette tips, can become brittle when exposed to prolonged irradiation.

- Do not place plastic containers that are not UV stable into the GuardOne® Workstation.

Operating the UV lamp

1.

To start the UV decontamination, switch the UV status (1) to **ON**.

To stop the UV decontamination, switch the UV status (1) to **OFF**.
- i

The UV lamp cannot be switched on while the front door is open.
- i

The UV lamp is switched off automatically as soon as the front door is being opened. Closing the front door again will not restart the decontamination procedure.
- i

The Info Expiration Time UV Light Source (2) gives you information about the remaining life time of the UV lamp.
2.

To exit the page, use Back or Home.

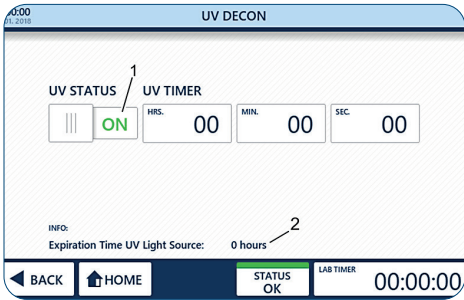
Fig. 36: UV 254 Decon

Operating the UV timer

> The timer can only be used before the UV lamp has been switched on.

**i** The UV timer is not mandatory for using UV irradiation.

- 1. Switch the UV Status to **OFF** (1).
- 2. To set the timer, press the hour, minute or second buttons (2).
- 3. Select the desired values (3) using the + and - button **- +**. The desired decontamination time has been set.
- 4. To start the UV decontamination, switch the UV status (1) to **ON**.

**i** The UV lamp is switched off automatically after the preset time has elapsed. If you do not set the timer, UV decontamination will run indefinitely.

**i** The UV lamp and timer will be switched off automatically as soon as the front door is being opened. Closing the front door again will not restart the decontamination procedure.

An acoustic signal will sound as soon as the pre-set time has elapsed.

- 5. To exit the page, use Back or Home.

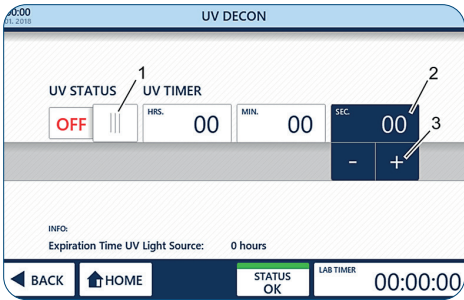
Fig. 37: UV Timer

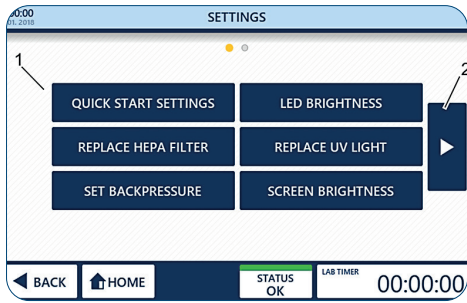
Fig. 38: Settings

## Settings

The Settings page allows you to adjust various parameters (1) of the GuardOne® Workstation. To navigate between the screens, use the arrow buttons (2) to the left and right of the main area.

### Quick Start Settings

By pressing Quick Start Settings, you can replace the factory settings for the blower and the LED lights with your preferred settings at any time during operation.

#### The factory settings are:

- > Blower speed is 50 %.
- > LED light is turned off.

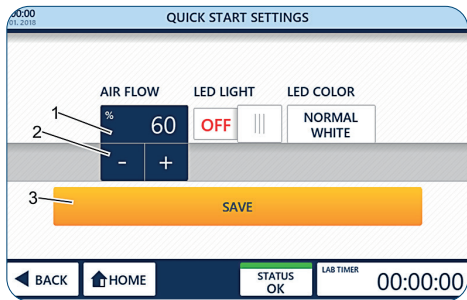
Fig. 39: Adjust the Airflow preference

### Adjusting the Airflow preference

1. Press Quick Start Settings.
2. Press the percentage value (1).
3. Select the desired airflow rate (2) using the + and - button .

**i** You can choose between 0, 50, 80 and 100 % airflow rate.

**i** Selecting an airflow rate of 0 % will switch off the blower. To switch it on again, either select an airflow rate above 0% or use the Quick Start function on the Home page Chapter 6.2 “Operating the Display Area”.

4. Apply the selection using Save (3).
5. To exit the page, use Back or Home.

**i** Exiting the page without using Save will discard all changes you have made!

## Setting the LED light preference

1. To switch on/off the LED light, press the LED Light switch (1) to ON/OFF
2. To select an LED light colour, press LED Colour (2).
3. Select the desired colour by pressing the + and - button  (3). You can choose between normal white (default) and red.
4. Apply the selection using Save (4).
5. To exit the page, use Back or Home.

**i** Exiting the page without using Save will discard all changes you have made!

## LED Brightness

By pressing LED Brightness, you can adjust the brightness of the LED light within the GuardOne® Workstation.

1. Press LED Brightness.
2. Press onto the percentage value (1).
3. Select the desired brightness rate (2) using the + and - button .

**i** You can choose from 10 % until 100 % LED brightness in 5 % increments.

**i** The selected LED brightness will be applied to all available LED colours.

4. Apply the selection using Save (3).
5. To exit the page, use Back or Home.

**i** Exiting the page without using Save will discard all changes you have made!

## Screen Brightness

By pressing Screen Brightness, you can adjust the brightness of the screen of the graphical user interface.

1. Press Screen Brightness.
2. Press the percentage value (1).
3. Select the desired brightness rate (2) using the + and - button .

**i** You can choose from 10 % to 100% Screen brightness.

4. Apply the selection using Save (3).
5. To exit the page, use Back or Home.

**i** Exiting the page without using Save will discard all changes you have made!

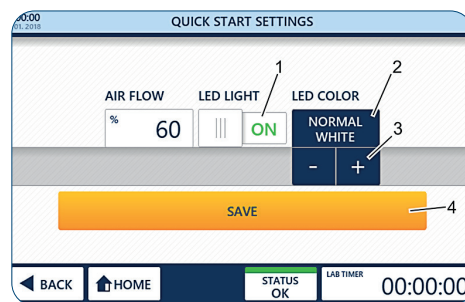
Fig. 40: Setting LED light and colour preference

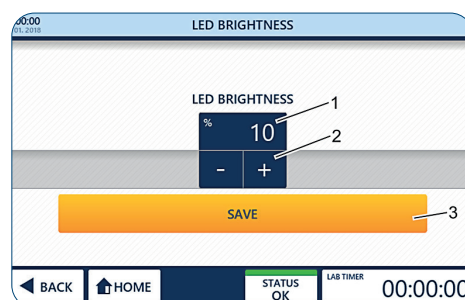
Fig. 41: Adjust the LED Brightness

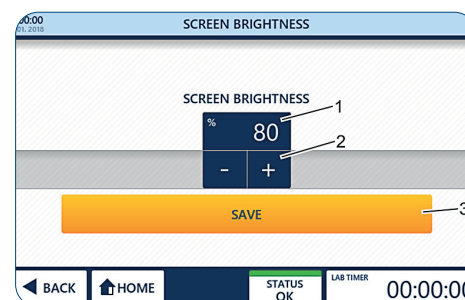
Fig. 42: Adjust the Screen Brightness

Replace HEPA Filter

**NOTICE!**  
**Restricted function due to the installation of a non-original filter!**

When using any other filter than the original Starlab HEPA filter, the proper function of the air filtration cannot be guaranteed.

- Only use an original Starlab HEPA filter designed for the GuardOne® Workstation.
- In case of questions, contact your local Starlab service or sales contact.

For ordering information, visit the website or use the contact information “Starlab Service Hotline” or the back page.

**i** The HEPA filter life time counter must only be reset after replacing the current old HEPA filter with a new HEPA filter.

For information on replacing the HEPA filter and activating the automatic life time surveillance for the HEPA filter, refer to Chapter 7.2 “Maintenance” .

Replace UV Light

**NOTICE!**  
**Restricted function due to the installation of a non-original UV lamp!**

When using any other UV lamp than the original Starlab UV lamp, the proper function of the UV decontamination cannot be guaranteed.

- Only use an original Starlab UV lamp designed for the GuardOne® Workstation.
- In case of any questions, contact your local Starlab service or sales representative.

For ordering information, visit the website or use the contact information “Starlab Service Hotline” or the back page.

**i** The UV light life time counter must only be reset after the replacement of the current old UV light with a new UV light.

For information on replacing the UV light and activating the automatic life time surveillance for the UV light, refer to Chapter 7.2 “Maintenance”.

Set Backpressure

**NOTICE!**  
**Limited effectiveness of the air filtration due to adjusting the HEPA filter airflow alarm threshold!**

Adjusting the threshold for the HEPA filter blockage alarm will change the effectiveness of the air filtration. Changing the factory settings may affect optimal performance of the GuardOne® Workstation.

- If you have the feeling that the alarm settings should be changed, contact a Starlab service technician or your local Starlab sales representative.

Reset

When Reset is pressed, the GuardOne® Workstation is reset to the factory settings.

Any changes made to screen brightness, LED brightness, quick start settings, time and date adjustment and backpressure adjustment are lost. After restoring the factory settings you will also have to calibrate the touchscreen.

1. Press Reset.
2. Confirm by pressing Reset (1).  
The GuardOne® Workstation turns off.
3. To calibrate the touchscreen, press onto the crosses that will subsequently appear in the four corners of the touchscreen with a pencil.  
The workstation has been reset to factory settings.  
To change the factory settings see Chapter 5 “Putting into service”.

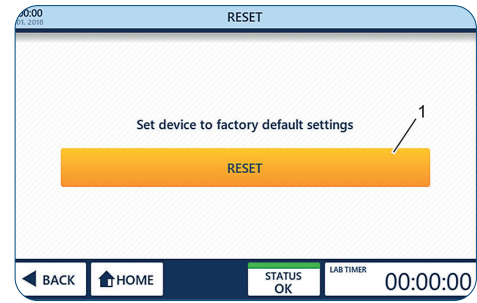
Fig. 43: Reset the GuardOne® Workstation

## 7 Cleaning and Maintenance

### 7.1 Safety Indications

#### Contamination indications

##### NOTICE!

##### **Contamination of the pre-filter due to contamination of the air!**

Due to the contamination of the air, the pre-filter can become dirty. A sterile working environment inside the GuardOne® Workstation is no longer guaranteed.

- To ensure optimal function of the pre-filter, it is recommended to visually inspect the pre-filter in regular intervals.
- It must be changed as soon as it visibly dirty, latest after a period of 6 months.

##### NOTICE!

##### **Contamination of the inside area of the GuardOne® Workstation due to bacterial contamination!**

Due to contamination of the work area, a repeated operation without cleaning the work area can lead to contamination or destruction of the sample material and falsification of the results.

- It is necessary to clean the inside area and the work surface after each use.

#### Property damage

##### NOTICE!

##### **Top housing damage by placing objects on it!**

If laboratory materials, samples or liquids are placed on the surface of the top housing, damage may occur if something gets into the prefilter and the HEPA filter.

- The surface of the top housing is not suitable as a storage area.

## 7.2 Maintenance

### Replacing the pre-filter

The pre-filter of the GuardOne® Workstation is located at the top of the device above the blower motor. Its function is to reduce the particle load for the main HEPA filter. To ensure the optimal function of the pre-filter, it is recommended to visually inspect the pre-filter in regular intervals. It must be exchanged as soon as it visibly dirty, latest after a period of 6 months.

##### NOTICE!

##### **Property damage due to non-simultaneous replacement of pre-filter and HEPA filter!**

Only replacing the HEPA filter but not the pre-filter can result in damage to the new HEPA filter.

- Always replace the pre-filter along with the HEPA filter.

Fig. 44: Lift the protective cover

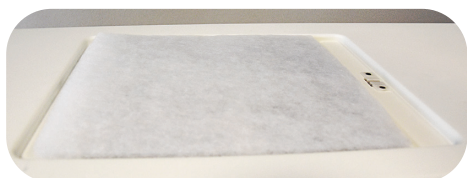
Fig. 45: Insert the new pre-filter

1. Lift the protective cover off from the top housing.
2. Remove the used pre-filter.
3. Insert the new pre-filter. In the 48" version, you replace two pre-filters side by side.
4. Re-insert the protective cover to fix the new pre-filter in place.

## Replacing the HEPA filter

The GuardOne® Workstation of size 32" is equipped with a single HEPA filter, while the 48" model has two HEPA filters. Both models are designed with the intention to make the exchange of the filter as easy and fast as possible.

### NOTICE!

#### Loss of filtration efficiency due to the incorrect version of HEPA filter!

Installing the wrong type of HEPA filter (e.g. different width, height or brand) may result in a partial or complete loss of filtration efficiency.

- Make sure, you have the correct version of replacement filter available.

### NOTICE!

#### Restricted function due to the installation of a non-original filter!

When using any other filter than the original Starlab HEPA filter, the proper function of the air filtration cannot be guaranteed.

- Only use original Starlab HEPA filters designed for the GuardOne® Workstation.
- In case of any questions, contact your local Starlab service or sales contact.

For ordering information, visit the website or use the contact information "Starlab Service Hotline" or the back page.

## Removing the used HEPA filter

1. Switch off the GuardOne® Workstation by using the main power switch on the back of the housing.
2. Open the front door of the top housing until the guiding rails on its left and right side lock and keep the door open.
3. Unlock the HEPA filter housing which is fixed with two black screws on its left and right side between the HEPA filter housing and the walls of the top housing.

**i** It is sufficient to open the screws by a few turns. It is not necessary to remove the screws.

4. Lift up the HEPA filter housing.
  - > To do so, simply pull the two lifting helps on the left and right side of the housing towards you until they lock into the groove in the filter housing bottom.
5. Remove the old HEPA filter by pulling it on its guiding rails towards you.

Fig. 46: Open the front door of the top housing

Fig. 47: Open the screws by two or three turns

Fig. 48: Lift the HEPA filter housing up

Fig. 49: Remove the HEPA filter

Inserting the new HEPA filter

1. Place the new HEPA filter on top of the guiding rails with the label containing the serial number facing towards you. Make sure, the sides of the HEPA filter and the guiding rails are aligned.
2. Slide the new HEPA filter back into the housing.
3. In the 48" version you replace two HEPA filters side by side.
4. Make a note of the serial number at the front of the HEPA filter. You will need it later for the life time surveillance activation.
5. Push back the lifting helps of the HEPA filter housing into their original position.  
The housing now rests directly on top of the isolation foam of the filter.
6. Secure the housing by retightening the two black screws on the left and right side between HEPA filter housing and walls of the top housing.
7. Make sure that the HEPA filter is placed flush inside the housing and cannot be moved.  
If secured tightly, the isolation foam on top and bottom of the filter will now prevent any airflow leakage.
8. Close the housing front door. To do so, first unlock the two lifting helps on the right and the left side by slightly lifting the front door on both sides upwards.
9. Switch on the power.

**i** The front door is equipped with a safety switch that automatically interrupts the power. If the touchscreen does not work after switching on the power, ensure that the front door is fully closed.

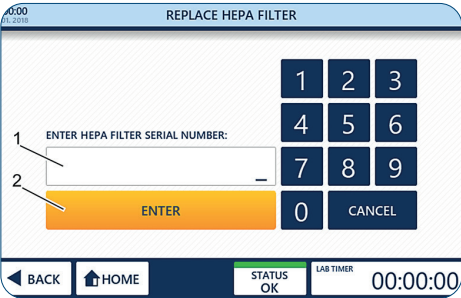
Fig. 50: Activate the automatic life time surveillance

Activating the automatic life time surveillance

1. To get to the Home page, press Tap to unlock.
2. Press Settings.
3. Press Replace HEPA Filter.
4. Enter the 10-digit serial number (1) shown on the front of the filter labels.
5. Press Enter (2) to finish the installation procedure.  
The GuardOne® Workstation is now ready to be used.  
The remaining life time of the HEPA filter is 600 days.

## Replacing the UV lamp

The 32" GuardOne® Workstation is equipped with a single UV lamp while the 48" model has two UV lamps installed. Both models are designed with the intention to make the exchange of the lamp as easy and fast as possible.

### Precondition:

- > Remove the packaging of the UV lamp. On the packaging, there is a sticker with the serial number of the UV lamp. You will need this serial number for the life time surveillance activation.

#### NOTICE!

##### **Risk of malfunctioning when exchanging only one UV lamp!**

The 48" model has two UV lamps installed. If you replace only one of them, there is a risk of malfunctioning.

- Always replace both UV lamps at the same time.

#### NOTICE!

##### **Loss of UV decontamination efficiency due to the incorrect version of UV lamp!**

Installing the wrong type of UV lamp (e.g. different power or brand) results in a partial or complete loss of UV decontamination efficiency.

- Make sure that you have the correct version of replacement UV lamp available.

#### NOTICE!

##### **Restricted function due to the installation of non-original UV lamp!**

When using any other UV lamp than the original Starlab UV lamp, the proper function of the UV decontamination cannot be guaranteed.

- Only use original Starlab UV lamp designed for the GuardOne® Workstation.
- In case of any questions, contact your local Starlab service or sales contact.

For ordering information, visit the website or use the contact information "Starlab Service Hotline" or the back page.

## Removing the used UV lamp

1. Switch off the GuardOne® Workstation by using the main power switch on the back of the housing.
2. Open the front door fully, using the recess on the roof of the top housing to safely secure it in this position.
- 3.

#### CAUTION!

##### **Risk of burns due to the heat of the UV lamp during decontamination!**

UV lamps can heat up during operation.

Touching the UV lamp immediately after use can cause minor burns and even blistering of the skin.

Wait at least 10 minutes before removing the cooled lamp from the holder.

4. Rotate the UV lamp approx. 90° in clockwise direction to release it from its socket. As soon as the lamp has been released, you will hear a clicking sound.
5. Withdraw the UV lamp downwards.

Fig. 51: Rotate the UV lamp

## Inserting the new UV lamp

1. Insert the new UV lamp from below into the lamp sockets. Lock the UV lamp by rotating it by approx. 90° in anticlockwise direction. In the 48" version you replace two UV lamps side by side according to the same principle. As soon as the lamp has been locked, you will hear a clicking sound.
2. Close the front door carefully.
3. Switch on the power.

Fig. 52: Activate the automatic life time surveillance

### Activating the automatic life time surveillance

1. To navigate to the Home page, press Tap to unlock.
2. Press Settings.
3. Press Replace UV Light.
4. Enter the 10-digit serial number shown on the sticker on the packaging of the UV lamp.
5. Press Enter to finish the installation procedure.  
The GuardOne® Workstation is now ready to be used again.  
The remaining life time of the UV bulb is 9000 hours.

### Exchanging the battery

The GuardOne® is equipped with a battery permitting device data storage when the device is switched off. Depending on the service life of the battery, this should be exchanged every 3–5 years.

#### WARNING

##### Data loss when removing the battery

Software version V2.9 and lower: Contact your local Starlab sales partners or Starlab service technician, as a software update to version V3.0 must first be performed before removing the battery to prevent the loss of the remaining service life of the UV lamp and the HEPA filter.

Information on the currently installed software version is displayed on the „Standby“ screen (see figure 27). Additionally, this information is also displayed under „Instrument Info“.

### Removing the battery

Abb. 53: Opening the lid of the upper housing part

Abb. 54: Rear of the LCD touch screen with battery opening

1. Switch off the GuardOne® workstation via the mains/power switch on the rear of the housing.
2. Open the lid of the upper housing part until the guiding rails to the left and right engage and keep the lid open.
3. The battery is located on the rear of the LCD touch screen on the lid of the upper housing part.

If your GuardOne® has no opening on the rear of the touch screen, contact your local Starlab sales partner or Starlab service technician. Contact information can be found on the status screen by pressing „Starlab Service & Order Information“.

4. Pressing the spring blade to the left releases the battery for removal.

## Inserting the battery

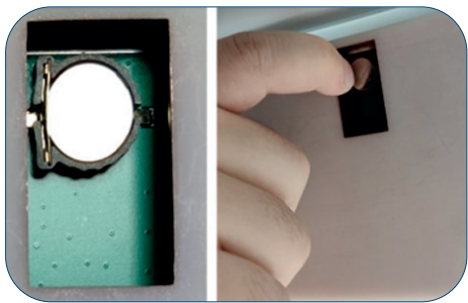
Abb. 55: Releasing the battery by pressing the spring blade on the left side of the battery to the left

1. Insert a new battery (type CR1220, button cell Ø 12.5 mm, 3 V / 35 mAh).
2. Close the lid of the upper housing part. To do this, first unlatch the two lifting aids on the left and right housing sides by slightly lifting the lid on both sides.
3. Switch on the mains/power supply.

**i** The lid is equipped with a safety switch that automatically disconnects the mains/power supply. If the touch screen does not function after the mains/power supply has been switched on, check that the lid has been closed correctly.

4. Enter the date and the time (see 5., „Commissioning – Adjusting the date and the time“)

7.3    Cleaning and Decontamination

Cleaning agents and tissues

To clean the GuardOne® Workstation, use the following cleaning agents and cleaning equipment:

- > lint-free cleaning tissue
- > alcohol-based disinfectant

 For cleaning, a wipe disinfection with alcohol based disinfectants is recommended. Further information can be found on the Starlab homepage.

Cleaning interval

The following intervals must be adhered to:

- > After each use of the GuardOne® Workstation, carry out the cleaning according to use. Clean the workstation with a recommended cleaning agent. Depending on the intended use of the GuardOne® Workstation, carry out the decontamination of the PC glass chamber by the UV lamp.
- > In case of a visible contamination, carry out appropriate cleaning.
- > To ensure the optimal function of the pre-filter, it is recommended to visually inspect the pre-filter in regular intervals. Exchange it as soon as it is visibly dirty, latest after a period of 6 months.

Cleaning the outside area of the GuardOne® Workstation

Preconditions:

- > The recommended cleaning intervals have been observed.
- > The GuardOne® Workstation is switched off or disconnected from the power supply.

1. Use a lint-free cleaning tissue to dust off the outer glass chamber of the GuardOne® Workstation.
2. If necessary, gently wipe the outside of the glass chamber with a mild detergent.

Cleaning the touchscreen

Preconditions:

- > The recommended cleaning intervals have been observed.
- > The GuardOne® Workstation is switched off or disconnected from the power supply.

1. Use a lint-free, slightly damp cleaning tissue to gently wipe the touchscreen.
2. In case of heavy soiling, apply a little alcohol-based detergent to the cleaning tissue. Gently wipe the touchscreen with the cleaning tissue.

## Cleaning the inside area and glass chamber of the GuardOne® Workstation

### CAUTION!

#### Health damage due to UV irradiation!

Long-term exposure to UV light results in damage to health.

- For safety reasons, the UV lamp cannot be switched on while the front door is open.
- For safety reasons, the UV lamp is switched off automatically as soon as the front door is being opened. Closing the door again does not restart the decontamination procedure.

### NOTICE!

#### Complete decontamination impossible due to permanent draught!

Due to the permanent draught in the glass chamber, a decontamination of 100 % by irradiation by the UV lamps is not possible.

- Follow the recommendations in this operating manual.
- Further information regarding the decontamination of the GuardOne® Workstations can be found on the Starlab homepage.

## Wet cleaning

### Preconditions:

- > The recommended cleaning intervals have been observed.
- > The GuardOne® Workstation is switched off.

1. Use a lint-free, damp cleaning tissue to wipe the interior of the glass chamber with a mild detergent and water.
2. Use a lint-free cleaning tissue to wipe dry.

## Cleaning by UV decontamination

1. Switch on the power.
2. Press Tap to unlock the Standby screen. The Standby screen is unlocked. You are now in the Home menu.
3. On the Home menu, press UV 254 Decon. The page for activating the UV decontamination opens.
4. To start the UV decontamination, switch on the UV lamp using the UV Status switch **ON**.  
To stop the UV decontamination, switch off the UV lamp using the UV Status switch **OFF** again.
5. To automatically stop the UV decontamination after a pre-set time, you can use the timer function:

**i** The timer can only be used before the UV lamp has been switched on.

**i** For a general decontamination of the internal space of the glass chamber, 20 minutes of irradiation is recommended. More information about decontamination after working with specific organisms can be found in "Bacterial destruction through UV irradiation".

- > To set the timer, press the hour, minute or second buttons.
  - > Select the desired values using the + and - button .
  - > After setting the desired decontamination time, switch on the UV lamp using the UV Status switch.
  - > The UV lamp will be switched off automatically after the pre-set time has elapsed.
  - > Recommended decontamination times can be found in "Bacterial destruction through UV irradiation".
6. After cleaning and decontamination, you can switch off the GuardOne® Workstation.

8    Troubleshooting

The GuardOne® Workstation was built with the intention to maximize user comfort and to reduce the amount of maintenance and service to the absolute necessity. To this end, the GuardOne® Workstation is constantly monitoring its vital parameters. The current device status is shown in the navigation bar.

**CAUTION!**  
**Health damage due to insufficient decontamination when returning device!**

An insufficient decontamination when returning the device to Starlab can constitute a health risk to any person handling the GuardOne® Workstation.

- If a device is returned to Starlab, a decontamination certificate must be sent along with it.
- Everything that is sent back must be cleaned first.
- The return must be registered with Starlab. Contact your local Starlab sales representative for further information. Contact details can be found by pressing Starlab Service & Order Information on the Status page.

**NOTICE!**  
**Property damage due to unauthorized repair of the GuardOne® Workstation!**

Independent repairs of the GuardOne® Workstation with an error status without the assistance of a Starlab service technician can lead to property damage.

- In case of an error status, contact your local Starlab sales representative. Contact details can be found by pressing Starlab Service & Order Information on the Status page.

**i** A fault is visually indicated by the colouring of the corresponding field.

- Green colour for an OK status.
- Orange colour for an Alert status.
- Red colour for an Error status.

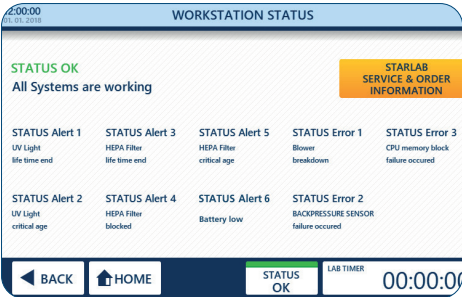
Fig. 56: Workstation status

**Workstation status**

1. Press Status (1) in the navigation bar.
  - › A green Status with OK indicates that all vital parameters are within their specifications and the GuardOne® Workstation is operating normally.
  - › An orange Status with Alert indicates that at least one parameter is outside the specifications. Nevertheless, the GuardOne® Workstation is still fully operational.
  - › A red Status with Error indicates that at least one parameter is outside the specifications and the GuardOne® Workstation cannot be used anymore.
2. For further information and contact details press Starlab Service & Order Information (2).

## 8.1 Alerts

- i** In case of an incoming alert, you can contact your local Starlab sales representative or Starlab service technician for further information and assistance at any time.

Status alerts reflect information about the functionality and life time of the HEPA filter and the UV lamp.

- > **UV light life time end** (Alert 1)
  - UV light life time end means that the UV lamp has reached the end time of 9000 hours and must be replaced to ensure the functionality of the decontamination process.
- > **UV light critical age** (Alert 2)
  - UV light critical age means that the UV lamp is close to the end of its life time of 9000 hours but has not reached it yet. 10% of the life time remains.
  - The UV decontamination process can still be used but it is recommended to have a spare UV lamp ready for future replacement.
  - To order a new UV lamp, contact your local Starlab sales representative. Contact details may be found by pressing Starlab Service & Order Information on the Status page.
- > **HEPA filter life time end** (Alert 3)
  - HEPA filter life time end means that the HEPA filter has reached its final life time of 600 days and must be replaced to ensure the functionality of the GuardOne® Workstation.
- > **HEPA filter blocked** (Alert 4)
  - HEPA filter blocked means that the GuardOne® Workstation has registered an impaired airflow that is usually caused by a blocked HEPA filter.
  - The HEPA filter must be replaced, irrespective of any remaining filter life time.
  - It is necessary to replace the HEPA filter immediately. Continuing the operation is not recommended.
  - Contact your Starlab service technician for further assistance.
- > **HEPA filter critical age** (Alert 5)
  - HEPA filter critical age means that the HEPA filter is close to the end of its life time of 600 days but has not reached it yet. 10 % of the life time remains.
  - The functionality of the GuardOne® Workstation is not impaired but it is recommended to have a spare HEPA filter ready for future replacement.
  - To order a new HEPA filter, contact your local Starlab sales representative. Contact details may be found by pressing Starlab Service & Order Information on the Status page.
- > **Battery low** (Alert 6)
  - „Battery Low“ means that the battery is nearing, but has not yet reached the end of its service life.
  - This does not affect the function of the GuardOne® workstation, but exchanging the battery soon is recommended.

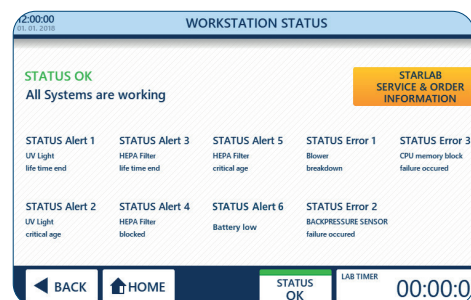
Fig. 57: Alerts

## 8.2 Errors

### NOTICE!

#### Property damage due to ignoring error messages!

In the event of an error message, any further use of the GuardOne® Workstation may result in material damage to the device and the sample material.

- Contact your Starlab service technician for further assistance.
- Contact details may be found by pressing Starlab Service & Order Information on the status page.

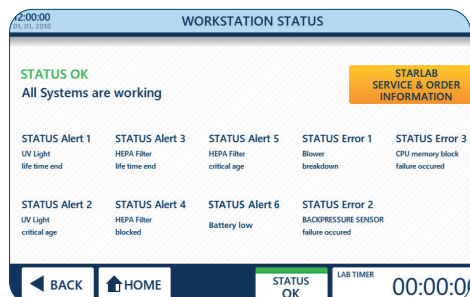
Fig. 58: Errors

Status errors inform about property damage that prevents the continued operation of the GuardOne® Workstation until repair by a Starlab service technician.

- › **Blower breakdown** (Error 1)
  - Blower breakdown means that the GuardOne® Workstation has detected a malfunction of the blower motor.
- › **Backpressure sensor failure occurred** (Error 2)
  - Backpressure Sensor failure occurred means that the GuardOne® Workstation has detected a malfunction of the airflow control.
- › **CPU memory block failure occurred** (Error 3)
  - CPU memory block failure occurred means that the GuardOne® Workstation has detected a malfunction of the device memory.

#### Behaviour in case of a status error

1. Contact your Starlab service technician for further assistance.
2. For contact details, press Starlab Service & Order Information on the status page.

## 8.3 Identifying non-automatically monitored malfunctions

### UV light does not work

In the case of a malfunctioning UV lamp, check the correct fitting of the magnetic switch cable.

1. Open the front door of the top housing.
2. Check the connection of the magnetic switch cable.
3. Make sure that the magnetic switch cable is connected.
4. Close the front door carefully.

**i** Make sure that the magnetic switch cable does not get caught between the front and the side flap of the top housing.

5. Enter the settings for the UV lamp again.

In case the UV lamp is still not working, contact your local Starlab sales representative or Starlab service technician. For contact details, press Starlab Service & Order Information on the status page.

### Incorrect date and time display

If the time and date entered is not saved or is no longer correct after switching the device off and on again, the battery must be exchanged.

**i** An incorrect date or time does not affect the function of the GuardOne®.

The exchange of the battery is described under 7.2, „Maintenance – battery exchange“. If problems with the date and time persist, please contact your local Starlab sales partner or Starlab service technician. Contact information can be found on the status screen by pressing „Starlab Service & Order Information“.

### Behaviour in case of a malfunctioning engine

If you hear abnormal engine running noise over a long period of time, there may be a malfunction of the engine.

1. Turn the main power switch to OFF.
2. Wait a few minutes before turning the GuardOne® Workstation on again.
3. Turn the main power switch to ON.

**i** If the noise of the engine persists, contact your local Starlab sales representative or Starlab service technician. For contact details, press Starlab Service & Order Information on the status page.

### Behaviour in case of a malfunctioning door switch

If the door switch is defective, the UV light can no longer be switched on. In addition, the airflow works even when the front door is closed. Contact your local Starlab sales representative or Starlab service technician for further assistance. For contact details, press Starlab Service & Order Information on the status page.

### Behaviour in case of a power failure

1. Turn the main power switch to ON.
2. Continue with your application.

**i** You do not have to reset previous settings. Previous settings are retained in the event of a power failure.

### Behaviour in case of a malfunctioning LED light

In case of a malfunctioning LED light bar, faults will not be displayed on the control display.

1. Turn the main power switch to OFF.
2. Wait a few minutes before turning on the GuardOne® Workstation again.
3. Turn the main power switch to ON.
4. Enter the settings for the LED light again.

**i** If the LED light still does not work, contact your local Starlab sales representative or Starlab service technician. For contact details, press Starlab Service & Order Information on the status page.

### Behaviour in case of a malfunctioning display

In case of a malfunctioning display, the touchscreen with the access to the graphical user interface does not respond to touches.

1. Turn the main power switch to OFF.
2. Wait a few minutes before turning on the GuardOne® Workstation again.
3. Turn the main power switch to ON.

**i** If the touchscreen still does not work, contact your local Starlab sales representative or Starlab service technician. Contact details can be found "STARLAB Service Hotline" or the back page.

### Behaviour in case of breakage due to transport

Contact your local Starlab sales representative or Starlab service technician. Contact details can be found "Starlab Service Hotline" or the back page.

## 9 Disposal

There is a warranty of two years on the GuardOne® Workstation. Disassembly and disposal is the responsibility of the operator.

**i** The life expectancy of the GuardOne® Workstation, excluding consumables, is at least five years, depending on the frequency of use and environmental conditions.

Unless there is a return or disposal agreement, lead in disassembled components to recycling:

- > Scrap metals.
- > Give plastic elements for recycling.
- > Dispose of other components sorted according to material quality.

### Disposal of the UV lamps

Defective UV lamps are hazardous waste. They contain quicksilver and cannot be recycled.

**i** Contact your in-house or local waste management system representative for assistance or further information.

In the event of disposing of the product, please observe the applicable legal regulations.

#### Information on the disposal of electrical and electronic devices in the European Community:

The disposal of electrical devices is regulated within the European Community by national regulations based on EU Directive 2012/19/EU pertaining to waste electrical and electronic equipment (WEEE).

In accordance with this, any devices delivered after 13/08/2005 on a business-to-business basis, which includes this product, may no longer be disposed of in household waste. To document this they have been marked with the icon on the left.

Because disposal regulations may differ from country to country within the EU, please contact your local distributor for information.

### Disposal of batteries

The adjacent symbol means that batteries / accumulators (rechargeable batteries) and electronic devices must be disposed of separately from household waste (unsorted municipal waste) at the end of their service life.

Electronic devices must be disposed of properly in accordance with Directive 2002/96/EC of the European Parliament and of the Council of January 27, 2003 on waste electrical and electronic equipment and in compliance with national disposal regulations.

Batteries and accumulators (rechargeable batteries) contain substances that can be harmful to the environment and to human health. Therefore, they must be disposed of properly in accordance with Directive 2006/66/EC of the European Parliament and of the Council of September 6, 2006 on batteries and accumulators and in compliance with national disposal regulations. Only dispose of fully discharged batteries and accumulators (rechargeable batteries).

**Warning!** Do not short-circuit batteries and accumulators (rechargeable batteries) to discharge them!

## 10 Specifications

### 10.1 Operating conditions

Operating conditions	Specifications	Value	Unit
	Permissible ambient temperature	+15 - +40	°C
	Permissible humidity	40 - 65	%
	Minimum life time	> 5	years
	Max. installation altitude above sea level	2000	m
Storage conditions	Specifications	Value	Unit
	Permissible storage temperature	-5 - +55	°C
	Permissible humidity	10 - 95	%

## 10.2 Measurements and weight

### External dimensions and weight package 32"

Specifications	Value	Unit
Length	920	mm
Width	680	mm
Height	510	mm
Weight		
Laminar Flow	40	kg
Dead Air	36	kg

### External dimensions and weight 32"

Specifications	Value	Unit
Length	813	mm
Width	610	mm
Height	877	mm
Weight		
Laminar Flow	36	kg
Dead Air	32	kg

### Internal dimensions 32"

Specifications	Value	Unit
Length	800	mm
Width	600	mm
Height	600	mm

### External dimensions and weight package 48"

Specifications	Value	Unit
Length	1330	mm
Width	680	mm
Height	520	mm
Weight		
Laminar Flow	47	kg
Dead Air	40	kg

### External dimensions and weight 48"

Specifications	Value	Unit
Length	1220	mm
Width	610	mm
Height	877	mm
Weight		
Laminar Flow	42	kg
Dead Air	35	kg

### Internal dimensions 48"

Specifications	Value	Unit
Length	1200	mm
Width	600	mm
Height	600	mm

### 10.3 Power requirements

#### Power requirements 32"

Specifications	Value	Unit
Maximum power consumption (Laminar Flow)	90	W
Maximum power consumption (Dead Air)	38	W
Voltage	200 - 240	V
Frequency	50/60	Hz
Protection class	5 (ISO 14644-1)	
Light intensity	1400	Lux
Airflow speed	0.2 - 0.4	m/s
HEPA filtration efficiency	>99.995	%
Sound level	55	dBA
UV-C Light	254-256	nm

#### Power requirements 48"

Specifications	Value	Unit
Maximum power consumption (Laminar Flow)	160	W
Maximum power consumption (Dead Air)	38	W
Voltage	200 - 240	V
Frequency	50/60	Hz
Protection class	5 (ISO 14644-1)	
Light intensity	1400	Lux
Airflow speed*	0.2 - 0.4	m/s
HEPA filtration efficiency	>99.995	%
Sound level	55	dBA
UV-C Light	254-256	nm

### 10.4 Product model label

 starlab

Order no.: SXXXX-0000

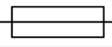
9206

Serial no.: XXXXXXXXXXXXXXX

200-240V~

50-60 Hz

180W

Fuse  T5A/250V

LABORATORY EQUIPMENT

Barcode area

CE UK CA

! (Warning symbol)

Made in China

Fig. 59: Product model label

The product model label contains the following information:

- > Order number
- > Serial number
- > Voltage
- > Frequency
- > Power output
- > Fuse
- > CE marking
- > WEEE label: Do not dispose with household waste.

## 11 Index

### A

Adjust	
Airflow .....	25, 30
Date and time .....	27
Default operation parameters .....	27
Lab timer .....	29
LED light .....	25, 30
UV lamp .....	26, 34
Airflow & Light	
Airflow .....	25, 30
LED light .....	25, 30
Alert	
Battery low .....	49
HEPA filter blocked .....	49
HEPA filter critical age .....	49
HEPA filter life time end .....	49
UV light critical age .....	49
UV light life time end .....	49
Assembly .....	18

### B

Back panel .....	6, 8, 19
Bacterial destruction .....	32
Base plate .....	6, 8, 19
Behaviour in case of emergency	
Accidents .....	15
Fire outbreak .....	15
Power failure .....	15

### C

Cleaning .....	40
Cleaning agents and tissues .....	44
Cleaning interval .....	44
Cleaning liquids .....	44
Cleaning the inside area and glass chamber of the GuardOne® Workstation .....	45
Cleaning the outside area of the GuardOne® Workstation .....	44
Cleaning the touchscreen .....	44
Cleaning and maintenance	
Safety indications .....	40
Coarse filter .....	8, 40

### D

Dead Air .....	6
Decontamination .....	32, 44
Display .....	6, 9, 28
Disposal .....	17, 52

### E

Error	
Backpressure sensor failure occurred .....	50
Behaviour in case of a malfunctioning display .....	50
Behaviour in case of a malfunctioning door switch .....	50
Behaviour in case of a malfunctioning engine .....	50
Behaviour in case of a malfunctioning LED lights .....	50
Behaviour in case of breakage due to transport .....	50
Behaviour in case of power failure .....	50
Blower breakdown .....	50
CPU memory block failure occurred .....	50
Date and Time are not displayed correctly .....	50
UV light does not work .....	50

### F

Front door .....	6, 8, 21
Function of the GuardOne® Workstation .....	8
Function test .....	24

### G

Graphical user interface .....	6, 9, 28
--------------------------------	----------

### H

Hazards .....	13
Cleaning liquids .....	15, 44
Dangers to persons .....	13
Dangers to the environment .....	14
Property damage .....	14
UV light .....	14, 52
HEPA filter .....	9, 24, 38
Home menu .....	24, 28

### I

In-house transport .....	17
Inspection	
Delivery .....	16
Transport .....	16
Installation .....	18
Installation site .....	18
Multifunctional shelf .....	20
PC back panel .....	19
PC front door .....	21
PC side panel .....	19
Top housing .....	22
UV lamp .....	23
Instrument Info .....	31
Intended use .....	11

### L

Laminar Flow .....	6
LED light .....	9, 25, 30, 51

### M

Main power connection .....	9, 24
Main power switch .....	9, 11, 24
Maintenance	
Replace the Battery .....	45
Replace the HEPA filter .....	41
Replace the pre-filter .....	40
Replace the UV lamp .....	43
Measurements and weight .....	43
Multifunctional shelf .....	8, 20

### O

Operating conditions .....	52
Operation	
Safety indications .....	28
Operator	
Obligations .....	11
Overview .....	6

### P

Packaging material .....	17
Packing .....	16
PC glass chamber .....	6, 8, 22
Personal protective equipment .....	12
Personnel	
Requirements .....	12
Power requirements .....	54
Pre-filter .....	9
Product model label .....	54
Putting into service .....	27

### Q

Quick Start .....	29
Quick Stop .....	29

### R

Recess .....	9, 43
--------------	-------

### S

Safety	
Danger to persons .....	13
Installations .....	11
Instructions .....	10
Procedures .....	10
Safety labelling	
GuardOne® Workstation .....	10
Scope of delivery .....	8
Settings .....	36
LED Brightness .....	37
Quick Start Settings .....	36
Replace HEPA filter .....	38
Replace UV light .....	38
Reset .....	39
Screen Brightness .....	37
Set Backpressure .....	38
Setup of the GuardOne® Workstation .....	6, 24
Side panel .....	6, 8, 19
Spare parts .....	9, 41, 43
Standby menu .....	24, 28
Storage .....	17
Conditions .....	52
Location .....	17
Requirements .....	17
Symbols	
At the GuardOne® Workstation .....	10
In the manual .....	10

### T

Top housing .....	6, 9, 22
Transport	
Safety .....	16
Troubleshooting .....	48
Control device status .....	48

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**Starlab GmbH**
**Starlab S.r.l.**