



PIPETTES

ErgoOne[®] E Electronic Pipettes

Reliable and ergonomic

ErgoOne^E

Single and multi-channel electronic pipettes with a lightweight and ergonomic design.



Reliable and Ergonomic

ErgoOne® E

ErgoOne® E single and multi-channel pipettes provide reliable pipetting every day. ErgoOne® E are microprocessor controlled, battery-operated, piston-stroke pipettes, which have been optimised for ergonomic operation and ease of use.

The lightweight, ergonomic design of ErgoOne® E pipettes are perfect for comfortable pipetting during long periods of use.

Five operating modes

1. Pipetting
2. Reverse Pipetting
3. Mixing
4. Multi Dispense
5. Electrophoresis
for loading electrophoresis gels. The gel is aspirated at high, adjustable speed and then slowly discharged

ErgoOne® E



Ergonomic, easy to reach pipetting button

Colour-coded ejector button
provides significantly reduced ejection forces

Adjustable finger rest
ErgoOne® E pipettes rest easily and ergonomically in your hand. The finger rest is height adjustable to allow the fine-tuning of the hand position for effortless operation of the function buttons

Comfortable, ergonomic hand grip

Autoclavable pipette shaft

Significantly reduced tip loading forces

Universal tip fit



Charging stand for single-channels



Single stand for single-channels



Single stand for multi-channels

ErgoOne® E Electronic Pipettes



Easy to read digital display

Large, digital volume display and easy to master intuitive menu structure makes the ErgoOne® E the researcher's choice for easy, everyday pipetting.



Long lasting battery

~4000 pipetting cycles on one charge! Special 'Battery Refresh Function' extends the service life and optimises battery performance. Auto shut-off after 10 minutes of inactivity helps save battery charge.



Can be recharged during use

You will never be without the use of your pipette – the ErgoOne® E can be recharged during use!

ErgoOne^E



Ordering Information

ErgoOne^E



ErgoOne^E Electronic Pipettes

Cat. No.	Volume Range	Increments	Volume μ l	Systematic Error (Inaccuracy)* $\leq \pm \%$	Random Error (Imprecision)* $\pm \%$	List Price
Single-Channel Pipettes						
G9001-0010	0.5–10 μ l (red)	0.01 μ l	1.0 / 5.0 / 10	5.0 / 1.5 / 1.0	2.0 / 0.8 / 0.4	£401.70
G9001-0020	2–20 μ l (red)	0.02 μ l	2.0 / 10 / 20	5.0 / 1.5 / 1.0	2.5 / 0.8 / 0.4	£401.70
G9001-0200	20–200 μ l (yellow)	0.2 μ l	20 / 100 / 200	4.0 / 1.2 / 0.8	0.6 / 0.3 / 0.2	£401.70
G9001-1000	100–1000 μ l (blue)	1.0 μ l	100 / 500 / 1000	3.0 / 1.0 / 0.6	0.6 / 0.3 / 0.2	£401.70
G9001-5000	500–5000 μ l (violet)	5.0 μ l	500 / 2500 / 5000	3.0 / 1.0 / 0.6	0.6 / 0.3 / 0.2	£427.45
8-Channel Pipettes						
G9008-0010	0.5–10 μ l (red)	0.01 μ l	1.0 / 5.0 / 10	8.0 / 2.0 / 1.2	4.0 / 1.5 / 0.80	£669.50
G9008-0020	1–20 μ l (red)	0.02 μ l	2.0 / 10 / 20	8.0 / 2.0 / 1.0	3.0 / 1.0 / 0.50	£669.50
G9008-0100	5–100 μ l (yellow)	0.1 μ l	10 / 50 / 100	4.0 / 1.6 / 0.8	1.5 / 0.4 / 0.25	£669.50
G9008-0300	15–300 μ l (green)	0.5 μ l	30 / 150 / 300	3.0 / 1.2 / 0.6	1.2 / 0.4 / 0.25	£669.50
12-Channel Pipettes						
G9012-0010	0.5–10 μ l (red)	0.01 μ l	1.0 / 5.0 / 10	8.0 / 2.0 / 1.2	4.0 / 1.5 / 0.80	£772.50
G9012-0020	1–20 μ l (red)	0.02 μ l	2.0 / 10 / 20	8.0 / 2.0 / 1.0	3.0 / 1.0 / 0.50	£772.50
G9012-0100	5–100 μ l (yellow)	0.1 μ l	10 / 50 / 100	4.0 / 1.6 / 0.8	1.5 / 0.4 / 0.25	£772.50
G9012-0300	15–300 μ l (green)	0.5 μ l	30 / 150 / 300	3.0 / 1.2 / 0.6	1.2 / 0.4 / 0.25	£772.50

Single-channel pipettes supplied with battery, AC Adapter and silicone oil. Multi-channel pipettes supplied with battery, AC Adapter, individual stand, a rack of TipOne[®] tips and silicone oil. Tested in accordance with ISO 8655. Pipettes supplied with one-year warranty. All list prices valid until 31 December 2016.

Accessories

Cat. No.	Description	Pack Size	List Price
G9999-1000	AC Adapter for Europe/UK (230V/50 Hz)	1	£32.96
G9999-1001	Charging Stand with AC Adapter for Europe/UK (230V/50 Hz) for Three Single-Channel Pipettes up to 1000 μ l	1	£194.67
G9999-1002	Individual Stand (for Single-Channel Pipettes, 10 μ l to 1000 μ l)	1	£41.20
G9999-1003	Individual Stand (for 5000 μ l Single-Channel Pipette)	1	£41.20
G9999-1004	Individual Stand (for all Multi-Channel Pipettes)	1	£26.78
G9999-1005	Replacement Battery	1	£22.66
G9999-1006	Silicone Oil for Piston (for Single-Channel Pipettes, 10 μ l to 1000 μ l)	1	£7.21
G9999-1007	Silicone Oil for Piston (for 5000 μ l Single-Channel and all Multi-Channel Pipettes)	1	£11.69
G9999-1008	Fluorstatic Grease for Guiding Axis (for Multi-Channel Pipettes)	1	£2.63



TipOne[®] Pipette Tips

Use STARLAB's TipOne[®] tips with ErgoOne[®] E pipettes for the perfect pipetting partnership. For ordering information, see our catalogue or www.starlab.co.uk