Magnetic Stirrer

Instruction Manual (v.1)

N2400-2010

Magnetic Stirrer

CE

www.starlabgroup.com

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STARLAB certifies that

N2400-2010 Magnetic Stirrer

has been manufactured and inspected in accordance

with the following European standards:

2006/95/EC Low Voltage Directive

2004/108/EC Electromagnetic Compatibility Directive

Harmonized Standards:

EN 61010-1:2010

EN 61326-1:2006

As an ISO 9001 and ISO 13485 certified company,
our working procedures and processes are checked and audited regularly
and the quality of our products is continuously monitored.

STARLAB International GmbH

MAGNETIC STIRRER

Cat. No. N2400-2010

1. Safety instructions

Before operation

- Read the Instruction Manual carefully before use.
- Ensure that only trained staff use the instrument. The instrument must only be opened by a qualified and trained technician.
- Make sure the socket is earthed before use.
- Set the instrument on a stable, clean, dry, non-slip and fireproof surface.
- Do not operate in hazardous areas.
- Check instrument for damage before every use. Do not use damaged components.
- The operation voltage noted on the label must match.
- Always disconnect the plug before fitting accessories.
- When working, wear personal protection to avoid the risk from splashing of liquids and possible glass breakage which can result from mechanical vibrations.

During operation

- Set the "speed-control knob" to the lowest position first and then increase the speed.
- Reduce the speed if the medium splashes out of the vessel or the instrument is unstable on the bench.

2. Product description

The instrument is designed for stirring. Normal use: The instrument is designed for use in professional laboratories. Do not use it in residential areas.

3. Inspection

Unpack the equipment carefully and check for any damages. If the instrument is faulty you should contact STARLAB or your STARLAB distributor promptly.

The packing includes:

- One Magnetic Stirrer N2400-2010
- One EU-power supply and UK-power supply
- One Instruction manual

Please check all items are present.

4. Ambient conditions

- Power supply: 100 – 240 V 50/60Hz
- Ambient temperature: 4 – 40 °C
- Relative humidity: < 80 %

Place the instrument on a safe and stable surface.

5. Operation

5.1 Control elements

5.2 Operating

Switch on the instrument at the power switch. The function "stirring" is switched on or off by rotating the speed control knob. The speed range is 100 to 1500 rpm.

6. Maintenance and cleaning

- Proper maintenance and cleaning will keep the instrument in good working order and will lengthen its life time.
- Be careful not spray cleaning fluid into the instrument. Unplug the power cord when cleaning.
- Please use only Isopropyl alcohol and water containing tenside. Before using other methods or cleaning agents the user must ascertain with the manufacturer that this will not destroy the instrument.
- Wear protective gloves during cleaning of the instrument.
- When not in use for long periods, switch off the instrument and store it in a dry, clean and stable place at room temperature.

Warranty

This instrument is under warranty to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty claim expires if the instrument is opened by untrained or unqualified staff. Our "General terms and conditions" apply.

Service

If help is needed, please contact STARLAB or your STARLAB distributor.

See www.starlabgroup.com for your local contact.

Please provide the customer care representative with the following information:

- Serial Number (on the rear panel)
- Description of problem (i.e., hardware or software)
- Methods and procedures adopted to resolve the problems
- Your contact information

7. Technical specifications

Parameter	Data
Voltage	100 – 240 V
Frequency	50/60 Hz
Power	10 W
Stirring points	1
Max. stirring quantity	3 l
Max. length, magnetic bar	50 mm
Motor type	DC motor
Speed	100 – 1500 rpm
Protection class	IP42
Permissible ambient temperature	4 – 40 °C
Permissible relative humidity	< 80 %
Weight	0.9 kg
Dimensions (WxDxH)	160 x 220 x 80 mm

8. Ordering information

Cat. No.	Description
N2400-2010	Magnetic Stirrer

Please visit our website www.starlabgroup.com, for details on STARLAB's other products for stirring and shaking.

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STARLAB GmbH
STARLAB (UK), Ltd
STARLAB S.r.l.