# *Microplate Mixer*

Instruction Manual (v.1)

**N2400-8040**

Microplate Mixer

CE

[www.starlabgroup.com](http://www.starlabgroup.com)

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*STARLAB certifies that*

### **N2400-8040 Microplate Mixer**

*has been manufactured and inspected in accordance*

*with the following European standards:*

**2006/95/EC** Low Voltage Directive

**2004/108/EC** Electromagnetic Compatibility Directive

*Harmonized Standards:*

**EN 61010-1:2010**

**EN 61326-1:2006**

As an ISO 9001 and ISO 13485 certified company,  
our working procedures and processes are checked and audited regularly  
and the quality of our products is continuously monitored.

**STARLAB International GmbH**

# MICROPLATE MIXER

Cat. No. N2400-8040

## 1. Safety instructions

### Before operation

- Read the Instruction Manual carefully before use.
- Ensure that only trained staff use the instrument. The instrument must only be opened by a qualified and trained technician.
- Make sure the socket is earthed before use.
- Set the instrument on a stable, clean, dry, non-slip and fireproof surface.
- Do not operate in hazardous areas.
- Safe operation is only guaranteed with the original accessories from STARLAB. You will find the complete list in chapter 8.
- Accessories must be securely attached.
- Before each use, check both the instrument and the accessories for damage. Do not use damaged components.
- The operation voltage noted on the label must match.
- Always disconnect the plug before fitting accessories.
- When working, wear personal protection to avoid the risk from splashing of liquids and possible glass breakage which can result from mechanical vibrations.

### During operation

- Set the "speed-control knob" to the lowest position first and then increase the speed.
- Reduce the speed if medium splashes out of the vessel or the instrument is unstable on the bench.
- Keep a minimum distance of 100 mm clear space around the instrument and away from walls and other instruments.

## 2. Product description

The instrument is designed for mixing liquids.

Normal use:

The instrument is designed for use in professional laboratories. Do not use it in residential areas.

### 3. Inspection

Unpack the equipment carefully and check for any damages. If the instrument is faulty you should contact STARLAB or your STARLAB distributor promptly.

The packing includes:

- One Microplate Mixer N2400-8040
- One EU-power supply and one UK-power supply
- One Instruction Manual

Please check all items are present.

### 4. Ambient conditions

Power supply: 100 – 240 V 50/60Hz

Ambient temperature: 4 – 40 °C

Relative humidity: < 80 %

Place the instrument on a safe and stable surface.

### 5. Operation

#### 5.1 Control elements

#### 5.2 Control

- Make sure that the required operating voltage and power supply voltage match.
- Plug in the power adapter.
- Switch on the device at the power switch.

- Set the rated mix speed and the device begins to work.
- Switch off the device at the power switch.

### **5.3 Speed range**

- Microplate Clamp: 0 – 1500 rpm
- Double Microplate Clamp: 0 – 1000 rpm

## **6. Maintenance and cleaning**

- Proper maintenance and cleaning will keep the instrument in good working order and will lengthen its life time.
- Be careful not spray cleaning fluid into the instrument. Unplug the power cord when cleaning.
- Please use only Isopropyl alcohol and water containing tenside. Before using other methods or cleaning agents the user must ascertain with the manufacturer that this will not destroy the instrument.
- Wear protective gloves during cleaning of the instrument.
- When not in use for long periods, switch off the instrument and store it in a dry, clean and stable place at room temperature.

### **Warranty**

This instrument is under warranty to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty claim expires if the instrument is opened by untrained or unqualified staff. Our “General terms and conditions” apply.

### **Service**

If help is needed, please contact STARLAB or your STARLAB distributor.

See **[www.starlabgroup.com](http://www.starlabgroup.com)** for your local contact.

Please provide the customer care representative with the following information:

- Serial Number (on the rear panel)
- Description of problem (i.e., hardware or software)
- Methods and procedures adopted to resolve the problems
- Your contact information

## 7. Technical specifications

Parameter	Data
Voltage	100 – 240 V
Frequency	50/60 Hz
Power	20 W
Mixing movement	Orbital
Orbital diameter	4.5 mm
Max. mixing weight (with attachments)	0.5 kg
Motor type	Brushless motor
Motor rating input	18 W
Motor rating output	10 W
Permissible ON time	100 %
Speed range	0 – 1500 rpm
Speed display	Scale
Dimensions (WxDxH)	150 x 260 x 80 mm
Weight	3 kg
Permitted ambient temperature	4 – 40 °C
Permitted relative humidity	< 80 %
Protection class acc. to DIN 60529	IP21

## 8. Ordering information

Cat. No.	Description
N2400-8040	Microplate Mixer

### 8.1 Accessories

Cat. No.	Description
N2400-8041	Replacement attachment for one microtitre plates
N2400-8042	Attachment for two microtitre plates
N2400-8043	Microplate Mixer Adapter for use with N2400-6012 & N2400-6013

Please visit our website [www.starlabgroup.com](http://www.starlabgroup.com), for details on our other instruments and our comprehensive range of consumables for the laboratory, including PlateOne deepwell and microplates.

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### **STARLAB GmbH**
### **STARLAB (UK), Ltd**
### **STARLAB S.r.l.**