

Figure 1.



Figure 2.

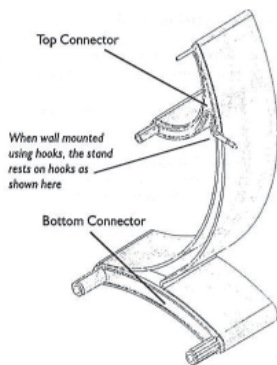


Figure 3.

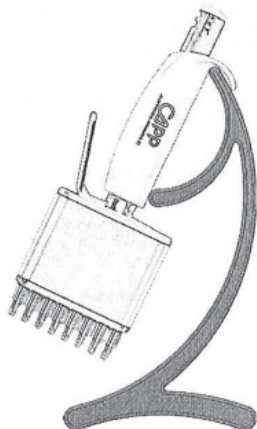


Figure 4.

Congratulations with your new CappAero CombiStand. The modular stand adds a pleasing design to your workspace as well as offering an elegant way to store your pipettes upright, which is very important in order to ensure the continuing functionality of your pipettes.

Contents

- 2 stand modules
- 2 connector sets (shown in fig. 1)
- One hook for wall mounting (requires three or more modules)
- Two strips of double adhesive tape for wall mounting without hooks
- User guide (this document)

Connecting modules

1. Push a top and a bottom connector into one module as shown in fig. 3.
2. Align the second module next to the first and push them firmly together.
3. Subsequent modules can be added using the second connector set.

Any number of stand modules can be connected this way to generate a pipette stand of a size to hold the desired number of CappAero pipettes.

Placing pipettes on the CombiStand

Pipettes are simply placed by hooking the curved finger rest of the pipette handle over the top of the stand and allowing the pipette to rest against the front edge of the stand (see fig. 4)

CappAero Multi pipettes are most optimally placed as shown in fig. 4 – in order to take up the smallest amount of space on the stand.

Wall mounting using tape

Wall mounting can be done using the double sided adhesive tape delivered with the CombiStand.

1. Please ensure that the wall surface is smooth, clean and non-porous.
2. Peel off the protective paper backing on one side of the tape and affix the tape to the CombiStand as shown in fig. 2.
3. Peel off the other paper backing and press the stand firmly onto the wall in the position desired.

Please take care to verify that the stand is positioned horizontally and in the position desired before attaching it to the wall as the tape does not allow repositioning.

Wall mounting using hooks

A stand made by connecting three or more modules can be wall mounted using the hooks supplied.

The stand rests on the hooks as shown in fig. 3.

1. Put the desired stand together using the appropriate number of modules (minimum 3).
2. Hold the stand to the wall in the desired position, checking that it is placed horizontally.
3. Mark where the hooks will be placed in the wall.
4. Setting the stand aside, proceed to mount the hooks into the wall (tools, plugs and other materials needed are not supplied).
5. Place the stand on the hooks.

Please note that you will need at least two hooks for the stand to hang stably and horizontally. For longer stands, more hooks may be needed to make the stand hang securely.