

STARTAG LARGE LASER SHEETS 90TV

LABEL SIZE: 32.5MM X 12.5MM

The following details will facilitate accurate configuration of the above laser labels with **Microsoft WORD 2007** for windows.

1. Open Word
2. Click on 'Mailings and Labels'
3. Select 'Labels' then 'Options'
4. Select 'Page Printers' and Tray - hand feeding is recommended – then 'Label Information' select 'Other / Custom'
5. Click on Details or New Label to customize your label
Top margin: 1.1 cm
Side margin: 1.5 cm
Vertical pitch: 1.55 cm
Horizontal pitch: 3.75 cm
Label height: 1.25 cm
Label width: 3.25 cm
Number across: 5
Number down: 18

The above settings can be used for previous versions of Word, however, the terminology may vary.

7. Set page size to A4 (21 x 29.7cm)
8. Select a name for the label; Large Lab Tags

Please Note

We recommend that prior to printing the labels, a print test is carried out on a plain A4 sheet, then held against label sheet to ensure that all print is included within the label area. Rigorous testing has shown that a slight variation may occur with different makes and models of Laser Printers.

STORAGE – PRIOR TO ATTACHMENT SHEETS (PRINTED & UNPRINTED) SHOULD BE STORED IN THE GRIP-SEAL POLYBAG AND CARDBOARD WALLET PROVIDED AT BETWEEN 10°C/25°C AND 35%/65% RH

THESE LABELS ARE NOT DESIGNED FOR USE WITH INKJET PRINTERS

BATCH NO: XXXXXX