Rocking Shaker and 3D Rocking Shaker

Instruction Manual (v.1)

N2400-8010 Rocking Shaker

N2400-8020 3D Rocking Shaker

CE

www.starlabgroup.com

ENGLISH3
DEUTSCH 9
FRANCAIS 14
ITALIANO 19

EC DECLARATION OF CONFORMITY

STARLAB certifies that

N2400-8010 Rocking Shaker

N2400-8020 3D Rocking shaker

has been manufactured and inspected in accordance

with the following European standards:

2006/95/EC Low Voltage Directive

2004/108/EC Electromagnetic Compatibility Directive

Harmonized Standards:

EN 61010-1:2010

EN 61326-1:2006

As an ISO 9001 and ISO 13485 certified company,
our working procedures and processes are checked and audited regularly
and the quality of our products is continuously monitored.

STARLAB International GmbH

ROCKING SHAKER Cat. No. N2400-8010**3D ROCKING SHAKER** Cat. No. N2400-8020**1. Safety instructions****Before operation**

- Read the Instruction Manual carefully before use.
- Ensure that only trained staff use the instrument. The instrument may only be opened by a qualified and trained technician.
- Make sure the socket is earthed before use.
- Set the instrument on a stable, clean, dry, non-slip and fireproof surface.
- Do not operate in hazardous areas.
- Safe operation is only guaranteed with the original accessories from STARLAB. You will find the complete list in chapter 8.
- Accessories must be securely attached.
- Check instrument and accessories beforehand for damage each time you use it. Do not use damaged components.
- The operation voltage mentioned on the label must match.
- Always disconnect the plug before fitting accessories.
- When working, wear personal protection to avoid the risk from splashing of liquids and possible glass breakage which can result from mechanical vibrations.

During operation

- Do not touch any running parts (fingers can get caught!)
- Do not attempt to move the instrument during operation.
- Set the "speed-control knob" to the lowest position first and then gradually increase the speed.
- Reduce the speed if the instrument does not run smoothly.
- Firmly secure the accessories and vessels in place to avoid damage or risk.
- Preparation of certain samples may lead to dangerous flammability. Only process samples that will not react dangerously.
- To ensure safety, follow the instructions for use for accessories. Please switch off the power before assembly of accessories and ensure both the instrument and the accessories are intact before switching on again.

2. Product description

The instrument is designed for mixing liquids. Normal use: The instrument is designed for use in professional laboratories. Do not use it in residential areas.

3. Inspection

Unpack the equipment carefully and check for any damages. If the instrument is faulty you should contact STARLAB or your STARLAB distributor promptly.

The packing includes:

- One Rocking Shaker or one 3D Rocking Shaker
- One EU-power supply and one UK-power supply
- One Shaking Table Attachment (N2400-8021)
- One Instruction Manual

Please check all items are present.

4. Ambient conditions

Power supply: 100 – 240 V 50/60Hz

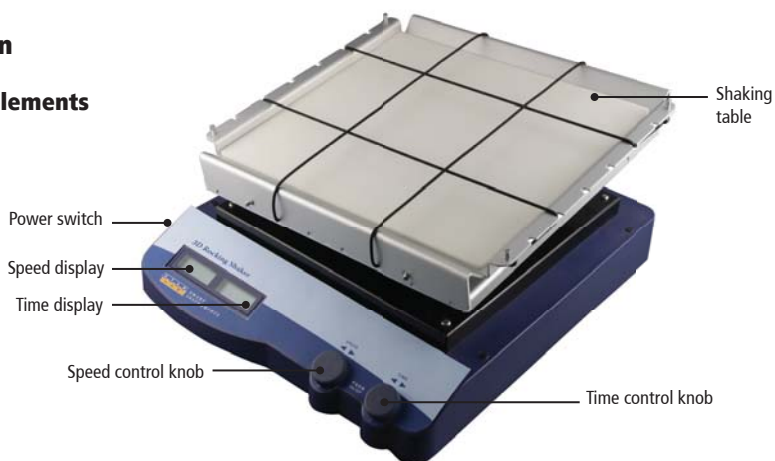
Ambient temperature: 4 – 40° C

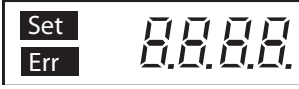
Relative humidity: < 80 %

Place the instrument in safe and stable surface.

5. Operation

5.1 Control elements

Item	Description
Speed control knob:	Sets target speed. Push on/off to start/stop the shaking function.
Timer control knob:	Sets target time. Push on/off to start/stop the shaking function.
Speed display:	Displays the set speed and actual speed values.
Time display:	Displays the set and remaining time values.
Power switch:	Switches the instrument ON/OFF.
Shaking table:	Place your sample containers here.
Set 	Displays when setting target value.
Err 	Displays when there is an error.
Display area 	When Set display, this area shows the value being set. When Set disappears, this area shows the running value.

5.2 Operating

- Place the instrument on a safe and stable surface. Ensure the required operating voltage and power supply voltage are the same, and that the socket is earthed.
- Connect the power cable.
- Switch ON instrument.
- The instrument begins self-checking.
- When initialization is over, the LCD displays the last running speed.
- Rotate the speed control knob to set the target speed.
- Press the speed control knob to start the rocking function. Press the speed control knob again to stop the rocking function.
- Rotate the timer control knob to set the target time. Press the timer control button to start the rocking function at the set speed and time. During operation, motion can be stopped at any time by pressing the speed control knob or timer control knob. If the timer control knob is pressed again, the

rocking function will start again and the timer will restart the countdown. When the timer reaches zero, the unit automatically stops and a beep will sound.

- If the rocking is smooth and regular, the instrument is working normally. If not, the instrument may be damaged, in which case contact STARLAB or your STARLAB distributor.

Troubleshooting

- The instrument does not work when switched ON:
 - check that the power cable is connected.
 - check that the fuse is not faulty.
- The rocking function suddenly stops:
 - if LCD displays "Er 03", this indicates the instrument is overloaded. Press the speed control knob or timer control knob first to deactivate the rocking function, and then restart with a lighter load. If the rocking function does not start, then repeat the process.

6. Maintenance and cleaning

- Proper maintenance and cleaning can keep instruments in good working order and will lengthen its lifetime.
- Be careful not to spray cleaning fluid into the instrument. Unplug the power cord when cleaning.
- Please use only Isopropyl alcohol and water containing tenside. Before using other methods or cleaning agents, the user must ascertain with the manufacturer that this will not destroy the instrument.
- Wear protective gloves during cleaning of the instrument.
- When not in use for long periods, switch off the instrument and store it in a dry, clean and stable place at room temperature.

Warranty

This instrument is under warranty to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty claim expires if the instrument is opened by untrained or unqualified staff. Our "General terms and conditions" apply.

Service

If help is needed, please contact STARLAB or your STARLAB distributor. See www.starlabgroup.com for your local contact. Please provide the customer care representative with the following information:

- Serial Number (on the rear panel)

- Description of problem (i.e., hardware or software)
- Methods and procedures adopted to resolve the problems
- Your contact information

7. Technical specifications

Parameter	Data for N2400-8010	Data for N2400-8020
Voltage	100 – 240 V	100 – 240 V
Frequency	50/60 Hz	50/60 Hz
Power	40 W	50 W
Angle	9°	9°
Max. Load	10 kg	5 kg
Motor Type	DC Motor	DC Motor
Speed range	10 – 70 rpm	10 – 70 rpm
Speed display	LCD	LCD
Time setting range	1 – 1199 min	1 – 1199 min
Time display	LCD	LCD
Operating modes	Two: timed & continuous	Two: timed & continuous
Dimensions (WxDxH)	370 x 420 x 106 mm	370 x 420 x 200 mm
Weight	9 kg	9 kg
Permissible ambient temperature	4 – 40 °C	4 – 40 °C
Permissible relative humidity	< 80 %	< 80 %
Safety class DIN EN60529	IP21	IP21

8. Ordering information

Cat. No.	Description
N2400-8010	Rocking Shaker incl. N2400-8021
N2400-8020	3D Rocking Shaker incl. N2400-8021

8.1 Accessories

Cat. No.	Description
N2400-8021	Universal/ Replacement attachment with bungee cords
N2400-8011	Universal attachment with clamping rolls

Please visit our website www.starlabgroup.com, for details on our range of other instruments and consumables for the laboratory.

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STARLAB GmbH
STARLAB (UK), Ltd
STARLAB S.r.l.